

Instruction Manual


## **ZEISS Axiolab 5, Axiolab 5 materials**

Upright Microscope for Routine and Entry-Level Research



## ZEISS Axiolab 5, Axiolab 5 materials

### Original Manual

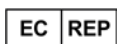

Carl Zeiss Microscopy GmbH  
Carl-Zeiss-Promenade 10  
07745 Jena  
Germany  
[info.microscopy.de@zeiss.com](mailto:info.microscopy.de@zeiss.com)  
[www.zeiss.com/microscopy](http://www.zeiss.com/microscopy)



Carl Zeiss AG  
Feldbachstr. 81  
8714 Feldbach  
Switzerland

### UK Responsible Person

Carl Zeiss Ltd  
1030 Cambourne Business Park, Cambourne  
CB23 6DW Cambridge  
United Kingdom



Carl Zeiss Suzhou Co., Ltd.  
Modern Industrial Square 3-B, No.333 XingPu Road SIP  
215126 Suzhou  
China

Document Name: Instruction Manual ZEISS Axiolab 5, Axiolab 5 materials

Order Number: 430037-7444-001

Revision: 11

Language: en-US

Effective from: 05/2024



© 2024 Without the prior written consent of ZEISS, this document or any part of it must neither be translated nor reproduced or transmitted in any form or by any means - including electronic or mechanic methods, by photocopying, recording or by any information or filing system. The right to make backup-copies for archiving purposes shall remain unaffected thereby. Any violations may be prosecuted as copyright infringements.

The use of general descriptive names, registered names, trademarks, etc. in this document does not imply that such names are exempt from the relevant intellectual property laws and regulations and therefore free for general use. This shall also apply if this is not specifically referred to. Software programs shall entirely remain the property of ZEISS. No program or subsequent upgrade thereof may be disclosed to any third party, copied or reproduced in any other form without the prior written consent of ZEISS, even if these copies or reproductions are destined for internal use at the customer's only, the only exception being one single back-up copy for archiving purposes.

# Table of Contents

<b>1</b>	<b>About this Instruction Manual.....</b>	<b>7</b>
1.1	Text Conventions and Link Types.....	7
1.2	Explanation of Warning Messages and Additional Information .....	8
1.3	Explanation of Symbols .....	9
1.4	Further Applicable Documents.....	10
1.5	Contact .....	10
<b>2</b>	<b>Safety .....</b>	<b>11</b>
2.1	Intended Purpose .....	11
2.2	General Safety Information.....	12
2.2.1	Requirements for Operators .....	12
2.2.2	Safe Operating Condition .....	12
2.2.3	Order and Use of Spare Parts .....	13
2.2.4	EMC Information.....	13
2.2.5	Optical Risk Grouping.....	14
2.2.6	Lifetime.....	14
2.3	Prevention of Hazards .....	14
2.3.1	Mechanical Hazards .....	14
2.3.2	Electrical Hazards .....	15
2.3.3	Hazards Generated with the Operating Environment.....	15
2.3.4	Ergonomic Hazards .....	15
2.3.5	Hazards Generated by Materials and Substances .....	16
2.3.6	Hazards Generated by Radiation.....	17
2.3.7	Thermal Hazards .....	17
2.4	Labels and Lights.....	17
2.4.1	Labels on the Axiolab 5 .....	18
2.5	Safety Devices and Interlocks.....	19
<b>3</b>	<b>Product and Functional Description .....</b>	<b>20</b>
3.1	Main Components.....	21
3.1.1	Axiolab 5 Bio-TL .....	21
3.1.2	Axiolab 5 Bio-TL/FL.....	22
3.1.3	Axiolab 5 Pol-TL .....	22
3.1.4	Axiolab 5 Pol-TL/conoscopy.....	23
3.1.5	Axiolab 5 Pol-TL/RL .....	23
3.1.6	Axiolab 5 Mat-TL/RL .....	24
3.2	Controls and Functional Elements on Components .....	25
3.2.1	Axiolab 5 Bio-TL Stand .....	25
3.2.2	Axiolab 5 Bio-TL/FL Stand.....	26
3.2.3	Axiolab 5 Pol-TL Stand .....	27
3.2.4	Axiolab 5 Pol-TL/conoscopy Stand.....	28
3.2.5	Axiolab 5 Pol-TL/RL Stand.....	29
3.2.6	Axiolab 5 Mat-TL/RL Stand .....	31
3.2.7	Functions of Stand Keys and Display Elements.....	32
3.2.8	Binocular Tube .....	34
3.2.9	Eyepieces .....	38
3.2.10	Eyepieces with Eyepiece Reticles.....	38
3.2.11	Nosepiece with Objectives.....	39
3.2.12	Condenser Carrier .....	39
3.2.13	Condensers .....	40

3.2.14	Stages .....	42
3.2.15	Reflector Turret .....	43
3.2.16	Storage Compartment .....	44
3.3	Light Manager Function .....	44
3.4	Microscopy and Contrast Techniques .....	45
3.4.1	Transmitted Light Brightfield Microscopy Using the KÖHLER Method ....	45
3.4.2	Transmitted Light Darkfield Microscopy Using the KÖHLER Method ....	45
3.4.3	Transmitted Light Phase Contrast Microscopy .....	45
3.4.4	Transmitted Light Polarization .....	46
3.4.5	Transmitted Light Polarization for Coniscope Observation.....	50
3.4.6	Reflected Light Brightfield Microscopy Using the KÖHLER Method .....	50
3.4.7	Reflected Light Darkfield Microscopy Using the KÖHLER Method .....	51
3.4.8	Reflected Light Polarization Microscopy.....	51
3.4.9	Reflected Light Fluorescence Microscopy.....	51
<b>4</b>	<b>Installation .....</b>	<b>52</b>
4.1	Unpacking and Setting up the Microscope .....	52
4.2	Assembling the Binocular Tube.....	53
4.3	Assembling the Binocular Tube with an Intermediate Plate .....	54
4.4	Assembling Components into the Binocular Tube .....	55
4.5	Assembling the Objectives .....	56
4.6	Assembling Condensers.....	57
4.6.1	Assembling the Condenser 0.9/1.25 BF, DF, Ph1, Ph2, Ph3 with Modu- lator Disk.....	57
4.6.2	Assembling the Dry Darkfield Condenser .....	58
4.6.3	Assembling the Condenser, Chromatic-aplanatic 0.9 BF DF PhC DIC.....	59
4.6.4	Assembling the Modulator Disk in the Condenser 0.9 BF Pol .....	60
4.7	Loading the Reflector Turret.....	61
4.7.1	Assembling Reflector Modules .....	61
4.7.2	Removing Reflector Modules.....	62
4.8	Connecting the Microscope to the Mains .....	62
<b>5</b>	<b>Operation .....</b>	<b>63</b>
5.1	Prerequisites for Commissioning and Operation.....	63
5.2	Switching On the Microscope.....	63
5.3	Adjusting .....	64
5.3.1	Adjusting the Position of the Eyepieces .....	64
5.3.2	Adjusting for Ametropia when Using Eyepiece Reticles .....	64
5.3.3	Adjusting the Height Stop on the Condenser Carrier .....	65
5.3.4	Using the Light Manager Function.....	66
5.3.5	Setting the ECO/Permanent Mode .....	68
5.4	Setting Up for Transmitted Light Techniques .....	68
5.4.1	Setting Up for Transmitted Light Brightfield Microscopy .....	68
5.4.2	Setting Up for Transmitted Light Darkfield Microscopy .....	71
5.4.3	Setting Up for Transmitted Light Phase Contrast Microscopy.....	74
5.4.4	Setting Up for Transmitted Light Polarization .....	74
5.5	Setting Up for Reflected Light Techniques .....	83
5.5.1	Setting Up for Reflected Light Brightfield Microscopy .....	83
5.5.2	Setting Up for Reflected Light Darkfield Microscopy .....	86
5.5.3	Setting Up for Reflected Light Polarization Microscopy .....	87
5.5.4	Setting Up for Reflected Light Fluorescence Microscopy .....	89
5.6	Switching Off the Microscope .....	92

<b>6</b>	<b>Care and Maintenance.....</b>	<b>93</b>
6.1	Safety during Cleaning and Maintenance .....	93
6.2	Maintenance Schedule .....	94
6.3	Maintenance Work.....	94
6.3.1	Cleaning an Optical Surface.....	94
6.3.2	Removing Water-Soluble Contamination .....	95
6.3.3	Exchanging the Fuses T 15 A/H 250 V in the Stand.....	95
6.3.4	Mechanical Stages .....	96
<b>7</b>	<b>Troubleshooting .....</b>	<b>98</b>
7.1	Resetting the Microscope to Factory Settings .....	101
<b>8</b>	<b>Decommissioning and Disposal .....</b>	<b>102</b>
8.1	Decommissioning .....	102
8.2	Transport and Storage.....	102
8.3	Disposal .....	103
8.4	Decontamination.....	103
<b>9</b>	<b>Technical Data and Conformity .....</b>	<b>104</b>
9.1	Performance Data and Specifications.....	104
9.2	Applicable Standards and Regulations .....	106
<b>10</b>	<b>Accessories and System Expansions .....</b>	<b>108</b>
10.1	Assembling the Reversible Eyecups .....	110
10.2	Binocular Tube .....	111
10.2.1	Binocular Tube 30°/23.....	111
10.2.2	Binocular Photo Tube Pol 20°/23 (100:0/0:100).....	111
10.2.3	Binocular Ergo Photo Tube 20°/23 (100:0/0:100) .....	112
10.2.4	Binocular Photo Tube 30°/23 (98:2), Reversed Image .....	113
10.2.5	Binocular Photo Tube 30°/23 (100:0/30:70/0:100), Reversed Image .....	114
10.3	Stages .....	114
10.3.1	Assembling a Mechanical Stage and Sample Holder .....	115
10.3.2	Removing a Mechanical Stage and Sample Holder.....	115
10.3.3	Rackless Mechanical Stage, 75x30 R.....	116
10.3.4	Mechanical Stage, 75x30 R .....	117
10.3.5	Rotary Stage Pol 360° with Sample Guide .....	118
10.4	Analyzer Sliders .....	123
10.4.1	Analyzer Slider TL/RL, Fixed .....	123
10.4.2	Analyzer Slider TL/RL, with Lambda Plate, 360° Rotatable .....	123
10.4.3	Analyzer Slider TL/RL with Lambda Plate, each Rotatable +/- 10° .....	123
10.5	Polarizers.....	124
10.5.1	Polarizer D, Fixed, Removable.....	124
10.5.2	Polarizer D, 90°, Rotatable, Removable .....	124
10.5.3	Polarizer, Fixed, with Lambda Plate, Rotatable.....	125
10.5.4	Polarizer, Rotatable, with Color Filter Carrier .....	125
10.5.5	Circular Polarizer D .....	126
10.5.6	Low-power System for Objectives 2.5x/4x .....	126
10.5.7	Color Filter Carrier 3x for Filter d=32mm .....	127
10.5.8	Assembling Polarizer or Color Filter Carrier onto the Condenser Carrier ..	128
10.6	DIC Slider C 6x20 .....	129

10.7	Condenser, achromatic-aplanatic 0.9 BF DF PhC DIC .....	129
10.7.1	Assembling the Condenser, achromatic-aplanatic 0.9 BF DF PhC DIC ....	130
10.7.2	Centering the Darkfield Diaphragm of the Condenser.....	131
10.7.3	Centering the Annular Phase Diaphragm of the Condenser .....	131
10.8	Light Source for Transmitted Light.....	132
10.8.1	Removing the Cover (Transmitted Light Source) .....	132
10.8.2	Exchanging the Halogen Lamp .....	133
10.8.3	Exchanging the LED Light Source .....	134
10.9	Light Source for Reflected Light.....	136
10.9.1	Removing the Cover (Reflected Light Source).....	136
10.9.2	Exchanging the Halogen Lamp .....	137
10.9.3	Exchanging the LED Light Source .....	137
10.9.4	Exchanging the LED Modules of the Fluorescence Light Source .....	140
10.10	Axiocam 202 Mono/208 Color .....	142
10.10.1	Assembling the Axiocam 202 mono or Axiocam 208 color .....	142
10.10.2	Operating Modes Using the Axiocam 202 mono/208 color .....	143
10.11	Assembling the Filter 32x4 mm on the Luminous-field Diaphragm Operating Ring .....	148
10.12	Dual Observation.....	149
10.12.1	Preparing the Binocular Tubes .....	150
10.12.2	Assembling the Dual Observation .....	152
10.12.3	Operating the Dual Observation .....	153
<b>Revision History.....</b>		<b>154</b>
<b>Glossary.....</b>		<b>155</b>
<b>Index .....</b>		<b>156</b>

# 1 About this Instruction Manual

This Instruction Manual (further called "document") is considered to be part of the Axiolab 5, Axiolab 5 materials, herein after referred to as "microscope".

This document contains basic steps and safety information that must be observed during operation and maintenance. Therefore, the document must be read by the operator prior to commissioning and must always be available at the place of use of the microscope.

This document is an essential part of the microscope and, if the microscope is resold, the document must remain with the microscope or be handed over to the new owner.

Axiolab 5 microscopes include:

- Axiolab 5 Bio-TL (430037-9011-000, 430037-9110-000, 430037-9060-000)
- Axiolab 5 Bio-TL/FL (430037-9021-000, 430037-9120-000, 430037-9070-000)
- Axiolab 5 Pol-TL (430037-9130-000)

Axiolab 5 materials microscopes include:

- Axiolab 5 Pol-TL/conoscopy (430037-9042-000)
- Axiolab 5 Pol-TL/RL (430037-9032-000)
- Axiolab 5 Mat-TL/RL (430037-9052-000)

## 1.1 Text Conventions and Link Types

Explanation	Example
Software controls and GUI elements.	Click <b>Start</b> .
Hardware controls and elements.	Press the <b>Standby</b> button.
Key on the keyboard.	Press <b>Enter</b> on the keyboard.
Press several keys on the keyboard simultaneously.	Press <b>Ctrl + Alt + Del</b> .
Follow a path in the software.	Select <b>Tools &gt; Goto Control Panel &gt; Airlock</b> .
Text to be entered by the user.	Enter <i>example.pdf</i> in this field.
Anything typed in literally during programming, for example macro codes and keywords.	Enter <code>Integer</code> in the console.
Link to further information within this document.	See: <i>Text Conventions and Link Types</i> [▶ 7].
Link to a website.	<a href="https://www.zeiss.com">https://www.zeiss.com</a>

## 1.2 Explanation of Warning Messages and Additional Information

DANGER, WARNING, CAUTION, and NOTICE are standard signal words used to determine the levels of hazards and risks of personal injury and property damage.

Always observe the safety and warning messages in **all** chapters of this document. Failure to comply with these instructions and warnings may result in personal injury, property damage, and the loss of any claims for damages.

The following warning messages indicating dangerous situations and hazards are used in this document.

### **DANGER**

#### **Type and source of danger**

DANGER indicates an imminently hazardous situation which, if not avoided, will result in death or serious injury.

### **WARNING**

#### **Type and source of danger**

WARNING indicates a potentially hazardous situation which, if not avoided, may result in death or serious injury.

### **CAUTION**

#### **Type and source of danger**

CAUTION indicates a potentially hazardous situation which, if not avoided, may result in minor or moderate injury.

### **NOTICE**

#### **Type and source of danger**













NOTICE indicates a potentially harmful situation which, if not avoided, may result in property damage.

### **Info**

Provides additional information or explanations to help the operator better understand the contents of this document.



### 1.3 Explanation of Symbols

	CE marking (Conformité Européene)
	CSA label: product tested by CSA to meet U.S. and Canadian standards. CSA approval master number optionally given adjacent to this symbol
	UKCA marking (UK conformity assessed)
	Manufacturer
	Country of manufacture. "CC" is the country code, e.g. "DE" for Germany, "CN" for China. Date of manufacture optionally given adjacent to this symbol
	Importer
	Authorized representative in the European Community
	Swiss authorized representative
	In vitro diagnostic medical device
	Serial number
	Catalogue number
	WEEE label: Do not discard as unsorted waste. Send to separate collection facilities for recovery and recycling

## 1.4 Further Applicable Documents

Please take also note of the following documents:

<b>Brochures and Certificates</b>	For brochures, certificates (e.g. ISO, CSA, SEMI), and declarations of conformity (e.g. EU, UK) ask your ZEISS Sales & Service Partner.
<b>Local and National Health and Safety Regulations</b>	Observe local and national health and safety regulations for the location of installation and during the use of the microscope. Consult with your ZEISS Sales & Service Partner if these regulations are in conflict with the installation requirements of the microscope.
<b>Software</b>	For detailed information on how to use Microscopy Apps, refer to its manual (e.g. Online Help, Software Manual) or ask your ZEISS Sales & Service Partner.
<b>System and Third-Party Components, Accessories</b>	Information about the individual components, enhancements, and accessories can be obtained from your ZEISS Sales & Service Partner. Also refer to the documentation of third-party manufacturers. <ul style="list-style-type: none"> <li>▪ Axiocam 202 mono</li> <li>▪ Axiocam 208 color</li> <li>▪ Multidiscussion Units</li> </ul>

## 1.5 Contact

If you have any questions or problems, contact your local ZEISS Sales & Service Partner or one of the following addresses:

### Headquarters

Phone:	+49 1803 33 63 34
Fax:	+49 3641 64 3439
Email:	info.microscopy.de@zeiss.com

### Microscopy Courses, Training, and Education

For information on microscopy courses, training, and education visit the ZEISS Academy Microscopy (<https://www.zeiss.com/microscopy/en/service-support/training-education/academy-microscopy.html>).

### ZEISS Portal

The ZEISS Portal (<https://portal.zeiss.com/>) offers various services that simplify the daily work with your ZEISS systems (machines and software).

### Service Germany

Phone:	+49 7364 20 3800
Fax:	+49 7364 20 3226
Email:	service.microscopy.de@zeiss.com

## 2 Safety

This chapter contains general requirements for safe working practices. Any person using the microscope or commissioned with installation or maintenance must read and observe these general safety instructions. Knowledge of basic safety instructions and requirements is a precondition for safe and fault-free operation. Operational safety of the supplied microscope is only ensured if it is operated according to its intended use.

If any work is associated with residual risks, this is mentioned in the relevant parts of this document in a specific note. When components must be handled with special caution, they are marked with a warning label. These warnings must always be observed.

Improper use of the microscope and its components can easily lead to impairment of their function or even damage them. Damage caused by incorrect operation, negligence, or unauthorized intervention, in particular by removing, modifying, or replacing parts of the microscope or its components, cannot be held liable by the device manufacturer. Third-party devices or components that are not expressly approved by ZEISS may not be used.

Any serious incident that has occurred in relation to the microscope and its components shall be reported to these institutions:

- the competent authority of the Member State in which the user is established
- ZEISS
  - for users within the EU:  
Carl Zeiss Microscopy GmbH, Jena, Germany
  - for users outside the EU:  
Carl Zeiss Suzhou Co., Ltd., Suzhou, China

### 2.1 Intended Purpose

The Axiolab 5 microscopes are instruments for general microscopic imaging for the in vitro examination of various biological samples including samples collected from humans or animals. This imaging provides information to further assess physiological and pathological conditions.

The microscopes are intended to be used by trained professionals only.

Axiolab 5 microscopes include:

- Axiolab 5 Bio-TL (430037-9011-000, 430037-9110-000, 430037-9060-000)
- Axiolab 5 Bio-TL/FL (430037-9021-000, 430037-9120-000, 430037-9070-000)
- Axiolab 5 Pol-TL (430037-9130-000)

The Axiolab 5 materials microscopes are universally usable microscopes for applications such as materials analyses. It is not intended to either directly or indirectly generate medical diagnostic results.

Axiolab 5 materials microscopes include:

- Axiolab 5 Pol-TL/conoscopy (430037-9042-000)
- Axiolab 5 Pol-TL/RL (430037-9032-000)
- Axiolab 5 Mat-TL/RL (430037-9052-000)

#### Info

The catalogue number can be found on the type plate, see *Labels and Lights* [▶ 17].

## 2.2 General Safety Information

This document must be read before commissioning in order to ensure safe and uninterrupted operation. Pay particular attention to all listed safety notes. Make sure, that

- the operating personnel has read and understood this manual, associated documents and particularly all safety regulations and instructions, and applies them.
- the local and national safety and accident prevention regulations must be observed, as well as the applicable laws and regulations in your country.
- this document is always available at the place of use of the microscope.
- the microscope is always in perfect condition.
- in case of defect or damage, the affected parts and the microscope are taken out of operation immediately and are secured against unintentional use.
- maintenance and repair work, retrofitting, removal or replacement of components, as well as any other intervention in the microscope not described in this document, may only be carried out by the manufacturer ZEISS or persons expressly authorized by ZEISS to do so.

### 2.2.1 Requirements for Operators

The microscope, components, and accessories may only be operated and maintained by authorized and trained personnel. The microscope may only be used in accordance with this document. If the microscope is not used as described, the safety of the user may be impaired and/or the microscope may be damaged.

Any unauthorized intervention or use other than within the scope of the intended use shall void all rights to warranty claims. The regional regulations on health protection and accident prevention must be observed at all times and during all work on and with the microscope.

**Training** Authorized ZEISS personnel will provide basic training in operating the microscope, as well as information on equipment safety and maintenance work that can be conducted by the operator. The training will be documented by ZEISS and its completion is to be confirmed by the operator. Special application training is offered for a fee. Current training dates, additional information and the registration form can be found at [www.zeiss.com](http://www.zeiss.com) or at the [ZEISS Portal](#).

### 2.2.2 Safe Operating Condition

If circumstances occur which impair safety and cause changes in operating behavior, the microscope and its components must be shut down immediately and a ZEISS service representative should be informed.

The microscope may only be operated if the operating conditions are adhered to.

- Do not operate the microscope and its components until you have completely read and understood the entire documentation.
- Make sure that all protective cover panels are installed and all warning labels are available and legible.
- Ensure conditions and take measures to prevent the build up of electrostatic charge on the workplace.

### 2.2.3 Order and Use of Spare Parts

Using spare parts that are not provided by ZEISS can be hazardous or can lead to property damage.

- Unless authorized by ZEISS, all spare parts should be installed by a ZEISS service representative.
- Contact your ZEISS service representative for information on spare parts order.
- Only genuine parts supplied by ZEISS are to be used in servicing the microscope and its components.

### 2.2.4 EMC Information

#### 2.2.4.1 Axiolab 5

The following EMC information is valid for the Axiolab 5 microscopes. For a definition of the included microscopes, see *Intended Purpose* [▶ 11].

The microscope is intended to be used in an industrial electromagnetic environment for non-clinical applications or in a home healthcare environment for clinical applications.

Use of this microscope in a dry environment, especially if synthetic materials are present (synthetic clothing, carpets, etc.), may cause electrostatic discharges that may cause erroneous results.

Do not use the microscope in proximity to sources of strong electromagnetic radiation, as these can interfere with proper operation.

If it is suspected that performance is affected by electromagnetic interference, correct operation may be restored by increasing the distance between the equipment and the source of the interference.

The microscope complies with the emission and immunity requirements as a CISPR 11 / EN 55011 / class B group 1 system according to IEC 61326-1 and IEC 61326-2-6. Emissions, which exceed the levels required by CISPR 11 / EN 55011, can occur when the microscope is connected to other devices.

The electromagnetic environment should be evaluated prior to operation of the microscope.

The following EMC user notice is for Korea only:

기종별	사용자안내문
B급기기 (가정용 방송통신기자재)	이 기기는 가정용(B급) 전자파적합기기로서 주로 가정에서 사용하는 것을 목적으로 하며, 모든 지역에서 사용할 수 있습니다.

#### 2.2.4.2 Axiolab 5 Materials

The following EMC information is valid for the Axiolab 5 materials microscopes. For a definition of the included microscopes, see *Intended Purpose* [▶ 11].

The microscope is intended to be used in an industrial electromagnetic environment.

The microscope complies with the emission and immunity requirements as a CISPR 11 / EN 55011 / class B group 1 system according to IEC 61326-1. Emissions, which exceed the levels required by CISPR 11 / EN 55011, can occur when the microscope is connected to other devices.

The following EMC user notice is for Korea only:

기종별	사용자안내문
B급기기 (가정용 방송통신기자재)	이 기기는 가정용(B급) 전자파적합기기로서 주로 가정에서 사용하는 것을 목적으로 하며, 모든 지역에서 사용할 수 있습니다.

### 2.2.5 Optical Risk Grouping

According to IEC 62471 sources of optical radiation are classified into risk groups subject to their potential photobiological hazard. Sources are classified into the following four groups according to hazard, based on the emission limit as well as permissible exposure time before hazard exceeded.

Risk group	Description
Exempt	No photobiological hazard.
1	No hazard due to normal behavioural limitations on exposure.
2	No hazard due to the aversion response to very bright light sources or thermal discomfort.
3	Hazardous even for momentary exposure.

The following table lists the risk grouping of the available light sources/illumination units according to the mentioned standard:

Light source/Illumination unit	Risk group
LED module 385 nm	3 (high risk)
LED modules 470 nm, 505 nm, 565 nm, 625 nm	2 (moderate risk)

### 2.2.6 Lifetime

A microscope is an opto-electronic device. Its availability for use is significantly determined by the performed maintenance. ZEISS guarantees the ability for maintenance and repair within eight years after initial operation. This is ensured by a corresponding service and spare parts concept, thus enabling the intended purpose within this duration.

## 2.3 Prevention of Hazards

This section summarizes potential hazards and recommended safety precautions. Failure to follow the safety instructions and instructions may result in personal injury and property damage.

### 2.3.1 Mechanical Hazards

#### Property Damage due to Transport

There is a risk of injury and property damage if the microscope is improperly handled and transported.

- Only use the handle, if applicable, for transport of the microscope. Otherwise hold the microscope with one hand and the base plate with the other hand.

### 2.3.2 Electrical Hazards

- Voltage Hazards** Risk of electric shock in case of contact with live parts.
- Detachable mains supply cords must not be replaced with inadequately rated cords.
  - Disconnect all power cords before cleaning.
  - Only connect electrical systems that are authorized by ZEISS to the supplied power supply cord.
  - Set up and operate the microscope so that the connectors are easily accessible.
  - Position the microscope in a way so that you can easily unplug the power cord at any time.
  - The microscope must be plugged into a properly installed power socket with protective earth contact using the supplied mains cords. The protective earth connection must not be impaired by the use of extension cables.
  - Always use the power cords supplied by ZEISS. When an unsuitable power cord is used, ZEISS can no longer guarantee the electrical safety and functionality of the microscope.
  - Shut down the microscope when not using the microscope.
  - Safe disconnection from the power supply is only ensured by pulling out the power cord. The switch on the microscope only switches to standby mode.

### 2.3.3 Hazards Generated with the Operating Environment

- Explosive Hazard** Fire hazard due to explosive or flammable environment.  
Do not operate the microscope and its components in a potentially explosive atmosphere, in the presence of volatile anesthetics or flammable solvents such as alcohol, petrol, or similar substances.
- Dirt, Dust, and Moisture** Dirt, dust, and moisture can impair the microscope's functionality.
- Shut down the microscope whenever it is not used and cover it with a dust protection cover.
  - Always cover unused openings/ports with the corresponding system component or with blind caps.
  - Perform regular maintenance and cleaning according to the instructions in this manual.
  - Make sure that no cleaning liquid or moisture gets inside the microscope.
  - Make sure that the electrical parts never come into contact with moisture.
  - Never expose the microscope to inadmissible climate conditions (high humidity and temperature).

### 2.3.4 Ergonomic Hazards

- Prevention of Musculoskeletal Disorders** Musculoskeletal disorders (MSDs) affect the muscles, nerves, blood vessels, ligaments and tendons. Workers in many different industries and occupations can be exposed to risk factors at work, such as lifting heavy items, bending, reaching overhead, pushing and pulling heavy loads, working in awkward body postures and performing the same or similar tasks repetitively. Employers are responsible for providing a safe and healthful workplace for their workers.

### 2.3.5 Hazards Generated by Materials and Substances

- |                              |  |
|------------------------------|--|
| <b>Infection Hazards</b>     | <p>Direct contact with the eyepieces can be a potential way of passing on bacterial and viral infections.</p> <ul style="list-style-type: none"> <li>▪ The risk can be lowered by using personal eyepieces or eyecups. If eyepieces need to be disinfected frequently, ZEISS recommends to use the eyepieces without eyecups.</li> <li>▪ To avoid infections, the use of personal protective equipment (PPE), e.g. gloves, for operation, cleaning, and decontamination is highly recommended. Disposable gloves can be decontaminated with alcohol for example, if necessary, or should be changed frequently to minimize the risk of contamination.</li> </ul>   |
| <b>Biological Hazard</b>     | <p>Biological substances/agents may pose a risk to the health of humans and other living organisms.</p> <ul style="list-style-type: none"> <li>▪ Keep a logbook of the known biological substances/agents used when working with the microscope and show it to the ZEISS service representative before they perform any work on the microscope.</li> </ul>   |
| <b>Consumable Hazards</b>    | <p>Incorrect handling of consumables and cleaning agents can lead to property damage or skin and eye injuries. Consumables that are not approved by ZEISS can lead to property damage. Consult your ZEISS Sales &amp; Service Partner to learn what consumables you can order and how to handle them.</p>  |
| <b>Immersion Oil</b>         | <p>Ensure that no immersion oil enters the surface water or the sewage system. The immersion oil can cause skin irritation.</p> <ul style="list-style-type: none"> <li>▪ Avoid any contact with skin, eyes and clothes.</li> <li>▪ Read and observe the safety data sheet of the immersion fluid.</li> <li>▪ In the event of skin contact, wash the oil off with plenty of water and soap.</li> <li>▪ In the event of eye contact, flush eyes with copious amounts of water for a minimum of 5 minutes. See a medical specialist if the irritation persists.</li> </ul>  |
| <b>Contamination Hazards</b> | <p>The microscope and other components can come into contact with various samples and substances that can be hazardous to humans and the environment. The microscope is not equipped with special user protection against substances that are corrosive, potentially infectious, toxic, radioactive or otherwise hazardous to health. When handling such substances, make sure to observe all legal regulations, particularly the relevant national accident prevention regulations.</p> <ul style="list-style-type: none"> <li>▪ Make sure that the microscope was not in contact with hazardous substances (check the laboratory logbook); otherwise, the microscope must be cleaned/decontaminated/disinfected.</li> <li>▪ Check the components also. If necessary, clean the components as meticulously as possible. Label contaminated/infected components that cannot be properly cleaned.</li> <li>▪ Direct contact with the eyepieces can be a potential way of passing on bacterial and viral infections. The risk can be lowered by using personal eyepieces or eyecups. If eyepieces need to be disinfected frequently, ZEISS recommends to use the eyepieces without eyecups.</li> <li>▪ Contaminated parts shall not be returned to any ZEISS department. Decontaminated parts can be sent to ZEISS accompanied by a signed „Customer Declaration of Decontamination“.</li> <li>▪ Wear gloves.</li> <li>▪ Observe all legal requirements, in particular the relevant national accident prevention regulations.</li> </ul> |



### 2.3.6 Hazards Generated by Radiation

**Optical Radiation Hazards** Gas discharge lights, LED lights and other sources of white light emit strong optical radiation (e.g. UV, VIS, IR). Optical radiation may cause damage to the skin and eyes. The extent of the damage depends on the parameters such as wavelength, exposure time, mode of operation (continuous or pulsed), etc.

- Avoid exposure of eyes and skin to radiation.
- Do not introduce reflective objects into the beam path.
- Never remove covers or cover panels during operation.
- Do not disable any interlock system elements.
- Use suitable protective equipment / protective clothing if required.

**Electromagnetic Radiation Hazards** The microscope may cause radio interference, which may be mitigated by relocating or re-orienting the equipment. The use of non-specified accessories, cables, or other auxiliary parts from the field of information technology may lead to increased electromagnetic emissions and reduced immunity to interference. Any integration into the system may result in a degradation of the EMC performance.

### 2.3.7 Thermal Hazards

**Burning Hazards** Hot surfaces, radiation and/or aggressive chemicals can cause burns.

- Use suitable protective equipment / protective clothing if mandatory.
- Always observe the cooling time of the hot surfaces.

**Heat Accumulation** Covering the ventilation openings can lead to heat accumulation that may damage the microscope and its components and, in extreme cases, can cause a fire.

- Keep ventilation openings unobstructed at all times.
- Do not cover devices or openings emitting heat.
- Comply with minimum distance from walls.

## 2.4 Labels and Lights

This chapter shows labels and, where applicable, indicator lights.

All parts that may pose specific hazards are marked with warning labels.

Always observe **all** warning labels!

- Check all warning labels for availability and legibility.
- Immediately replace damaged or illegible warning labels.

In case a label is missing, contact your ZEISS service representative for free of charge replacement.

2.4.1 Labels on the Axiolab 5

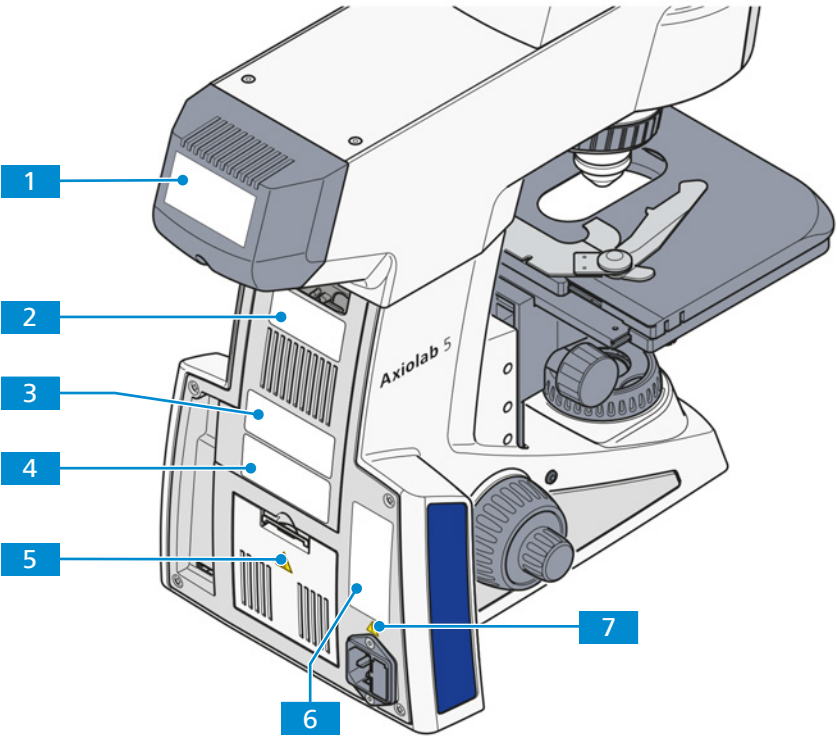







Fig. 1: Position of warning labels Axiolab 5 for transmitted light and reflected light

Pos.	Label or light	Explanation
1		Only for reflected light: CAUTION LED Radiation Do not stare at operating lamp. May be harmful to the eyes.
		Only for reflected light when the 385 nm LED is used: CAUTION UV Radiation Do not stare at operating lamp. May be harmful to the eyes.
2		UDI label
3		Microscope type label

Pos.	Label or light	Explanation
4		Microscope type label not applicable for Axiolab 5 materials microscopes
		Microscope type label only applicable for Axiolab 5 materials microscopes
5		Hot surface! Do not touch.
6		Axiolab 5, Axiolab 5 materials EU Representative Carl Zeiss Microscopy GmbH Carl-Zeiss-Promenade 10 07745 Jena, Germany
7		Observe the safety notes in the instruction manual and the supplied documents.

## 2.5 Safety Devices and Interlocks

In order to prevent injuries and/or property damage, the microscope and its components are equipped with several safety devices and interlocks. In case of defect or damage, the affected parts and the microscope must be taken out of operation immediately and must be secured against unintentional use.

To verify the safety of the microscope and its components, contact your ZEISS service representative and keep the service logs and logbooks.

## 3 Product and Functional Description

The Axiolab 5, Axiolab 5 materials is a transmitted-light microscope of compact design with a small footprint. The microscope provides high resolution, infinity-corrected objectives for transmitted-light techniques and for reflected-light techniques, depending on the microscope type.

Depending on the configuration of the microscope, the following microscopy and contrast techniques are available:

- |                               |   |
|-------------------------------|---|
| <b>Transmitted light (TL)</b> | <ul style="list-style-type: none"> <li>▪ <i>Brightfield (BF)</i> [<a href="#">▶ 45</a>]</li> <li>▪ <i>Darkfield (DF)</i> [<a href="#">▶ 45</a>]</li> <li>▪ <i>Phase contrast (Ph)</i> [<a href="#">▶ 45</a>]</li> <li>▪ <i>Polarization (Pol)</i> [<a href="#">▶ 46</a>]</li> <li>▪ <i>Polarization (Conoscopy)</i> [<a href="#">▶ 50</a>]</li> </ul> |
| <b>Reflected light (RL)</b>   | <ul style="list-style-type: none"> <li>▪ <i>Brightfield (BF)</i> [<a href="#">▶ 50</a>]</li> <li>▪ <i>Darkfield (DF)</i> [<a href="#">▶ 51</a>]</li> <li>▪ <i>Polarization (Pol)</i> [<a href="#">▶ 51</a>]</li> <li>▪ <i>Fluorescence (FL)</i> [<a href="#">▶ 51</a>]</li> </ul>   |

The following microscope types are available:

Axiolab 5 Bio-TL	Transmitted light stand for bioscience
Axiolab 5 Bio-TL/FL	Transmitted light and reflected light fluorescence stand for bioscience
Axiolab 5 Pol-TL	Transmitted light stand for polarization
Axiolab 5 Pol-TL/conoscopy	Transmitted light stand for polarization/conoscopy
Axiolab 5 Pol-TL/RL	Transmitted light and reflected light stand for polarization
Axiolab 5 Mat-TL/RL	Transmitted light and reflected light materials stand

### Typical Applications

Axiolab 5

- examination of blood and tissue samples taken from the human body, from plants, or animals
- medical examinations in laboratories, hospitals, and doctors' offices
- academic and practical education in medicine and biology
- industrial applications, e.g. in pharmacology, food technology, and wastewater examination

Axiolab 5 materials

- metallographic laboratories
- automotive industry
- microsystems engineering
- geoscientific institutes
- mineral exploration industry

### Info

Additional information about the hardware configuration and optional enhancements can be obtained from your ZEISS Sales & Service Partner.

### 3.1 Main Components

#### 3.1.1 Axiolab 5 Bio-TL

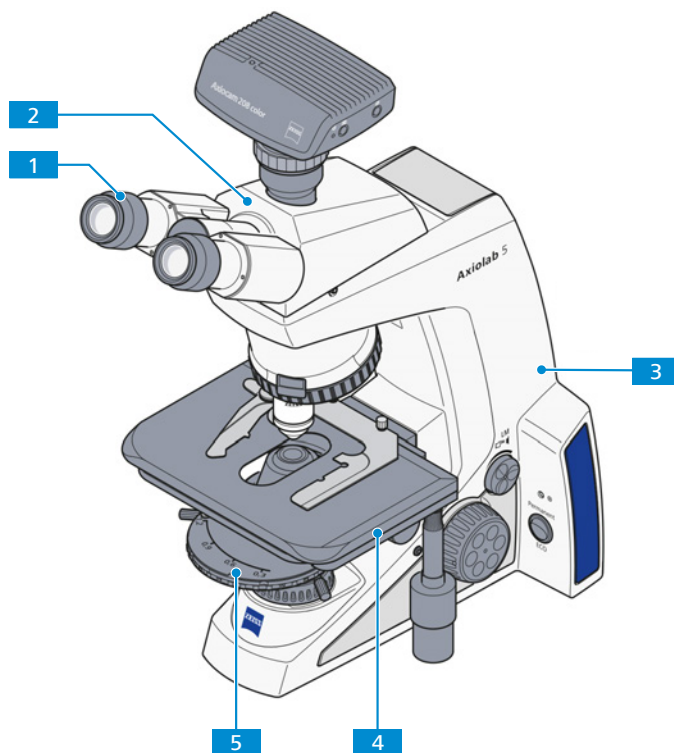


Fig. 2: Main components - Axiolab 5 Bio-TL

- |   |  |
|---|--|
| <b>1</b> Eyepieces [ <a href="#">▶ 38</a> ] | <b>2</b> Binocular tube [ <a href="#">▶ 34</a> ]   |
| <b>3</b> Stand Bio/TL                       | <b>4</b> Mechanical stage [ <a href="#">▶ 42</a> ] |
| <b>5</b> Condenser [ <a href="#">▶ 40</a> ] |  |

### 3.1.2 Axiolab 5 Bio-TL/FL

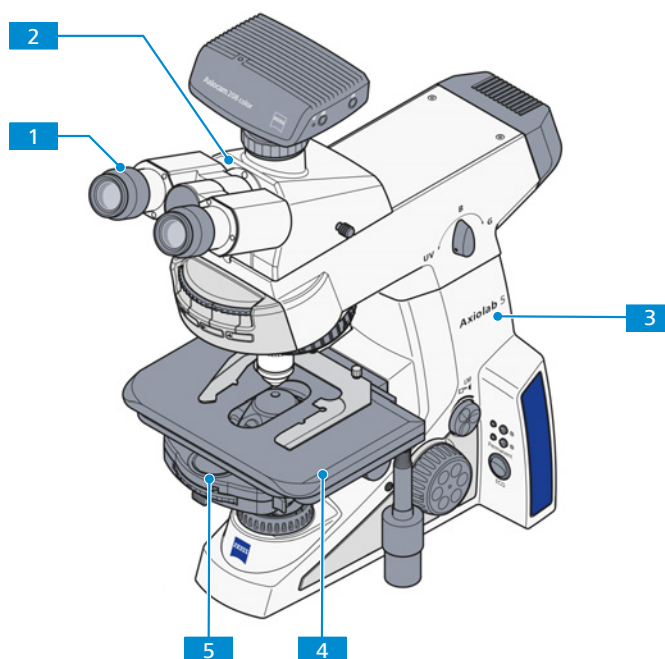


Fig. 3: Main components - Axiolab 5 Bio-TL/FL

- |                           |                                  |
|---------------------------|----------------------------------|
| <b>1</b> Eyepieces [▶ 38] | <b>2</b> Binocular tube [▶ 34]   |
| <b>3</b> Stand Bio-TL/FL  | <b>4</b> Mechanical stage [▶ 42] |
| <b>5</b> Condenser [▶ 40] |                                  |

### 3.1.3 Axiolab 5 Pol-TL

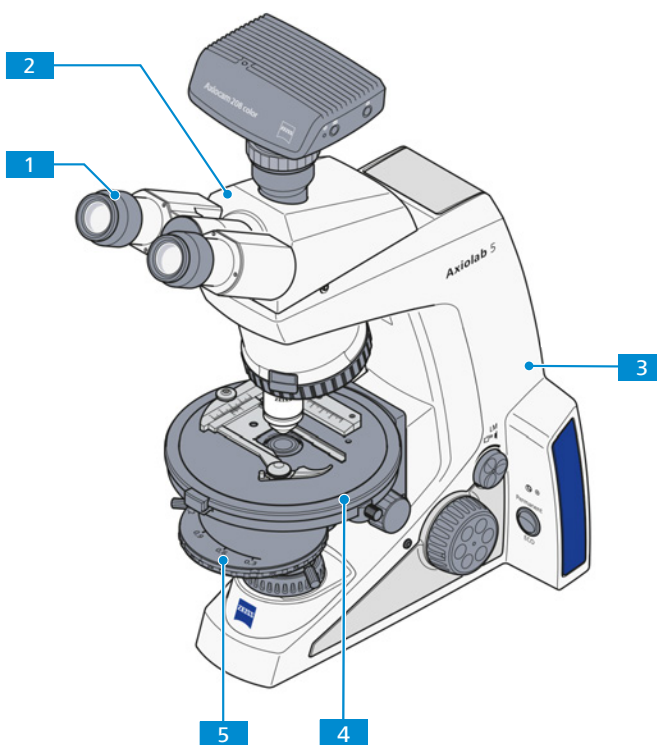


Fig. 4: Main components - Axiolab 5 Pol-TL

- |                           |                                |
|---------------------------|--------------------------------|
| <b>1</b> Eyepieces [▶ 38] | <b>2</b> Binocular tube [▶ 34] |
| <b>3</b> Stand Pol-TL     | <b>4</b> Rotary stage [▶ 42]   |
| <b>5</b> Condenser [▶ 40] |                                |

### 3.1.4 Axiolab 5 Pol-TL/conoscopy

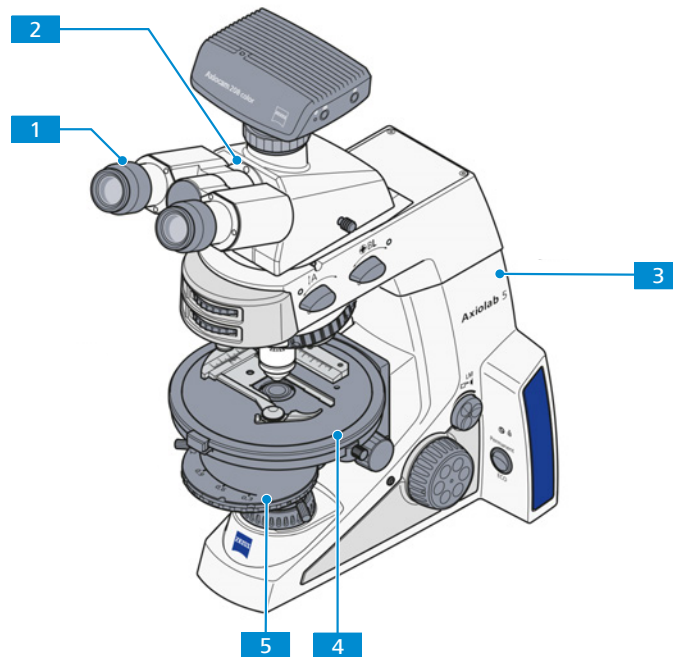


Fig. 5: Main components - Axiolab 5 Pol-TL/conoscopy

- |                                 |                                |
|---------------------------------|--------------------------------|
| <b>1</b> Eyepieces [▶ 38]       | <b>2</b> Binocular tube [▶ 34] |
| <b>3</b> Stand Pol-TL/conoscopy | <b>4</b> Rotary stage [▶ 42]   |
| <b>5</b> Condenser [▶ 40]       |                                |

### 3.1.5 Axiolab 5 Pol-TL/RL

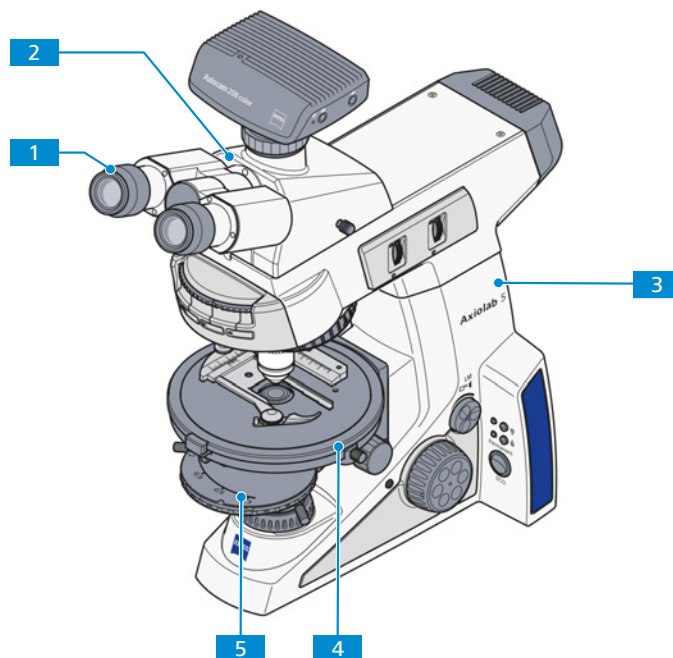


Fig. 6: Main components - Axiolab 5 Pol-TL/RL

- |                           |                                |
|---------------------------|--------------------------------|
| <b>1</b> Eyepieces [▶ 38] | <b>2</b> Binocular tube [▶ 34] |
| <b>3</b> Stand Pol-TL/RL  | <b>4</b> Rotary stage [▶ 42]   |
| <b>5</b> Condenser [▶ 40] |                                |



## 3.1.6 Axiolab 5 Mat-TL/RL

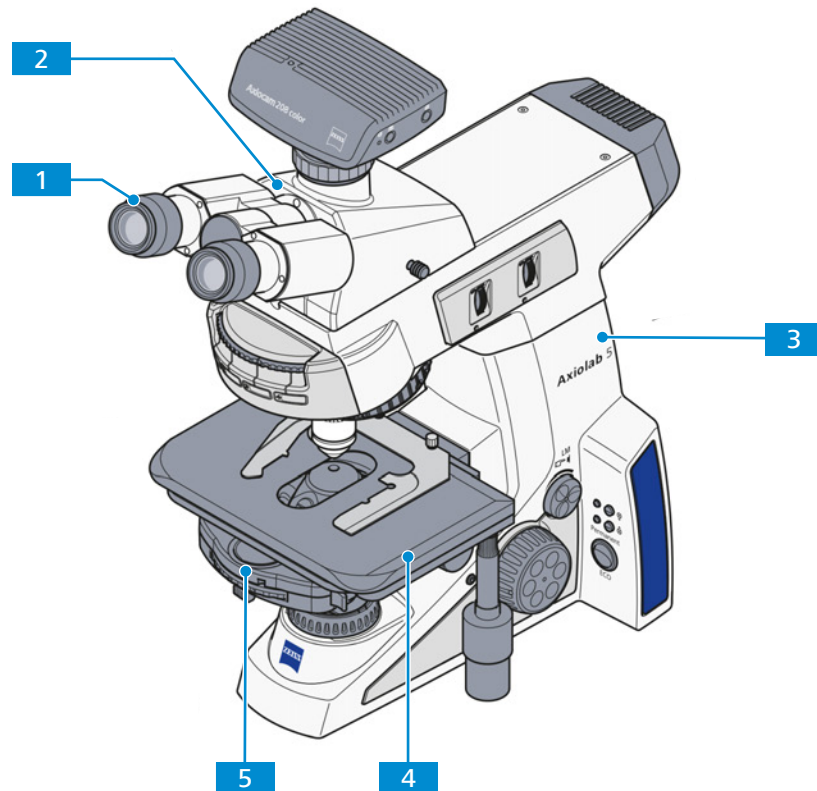


Fig. 7: Main component - Axiolab 5 Mat-TL/RL

- |                           |                                  |
|---------------------------|----------------------------------|
| <b>1</b> Eyepieces [▶ 38] | <b>2</b> Binocular tube [▶ 34]   |
| <b>3</b> Stand Mat-TL/RL  | <b>4</b> Mechanical stage [▶ 42] |
| <b>5</b> Condenser [▶ 40] |                                  |



## 3.2 Controls and Functional Elements on Components

### 3.2.1 Axiolab 5 Bio-TL Stand

This section shows the functional components and controls of stand Bio-TL.

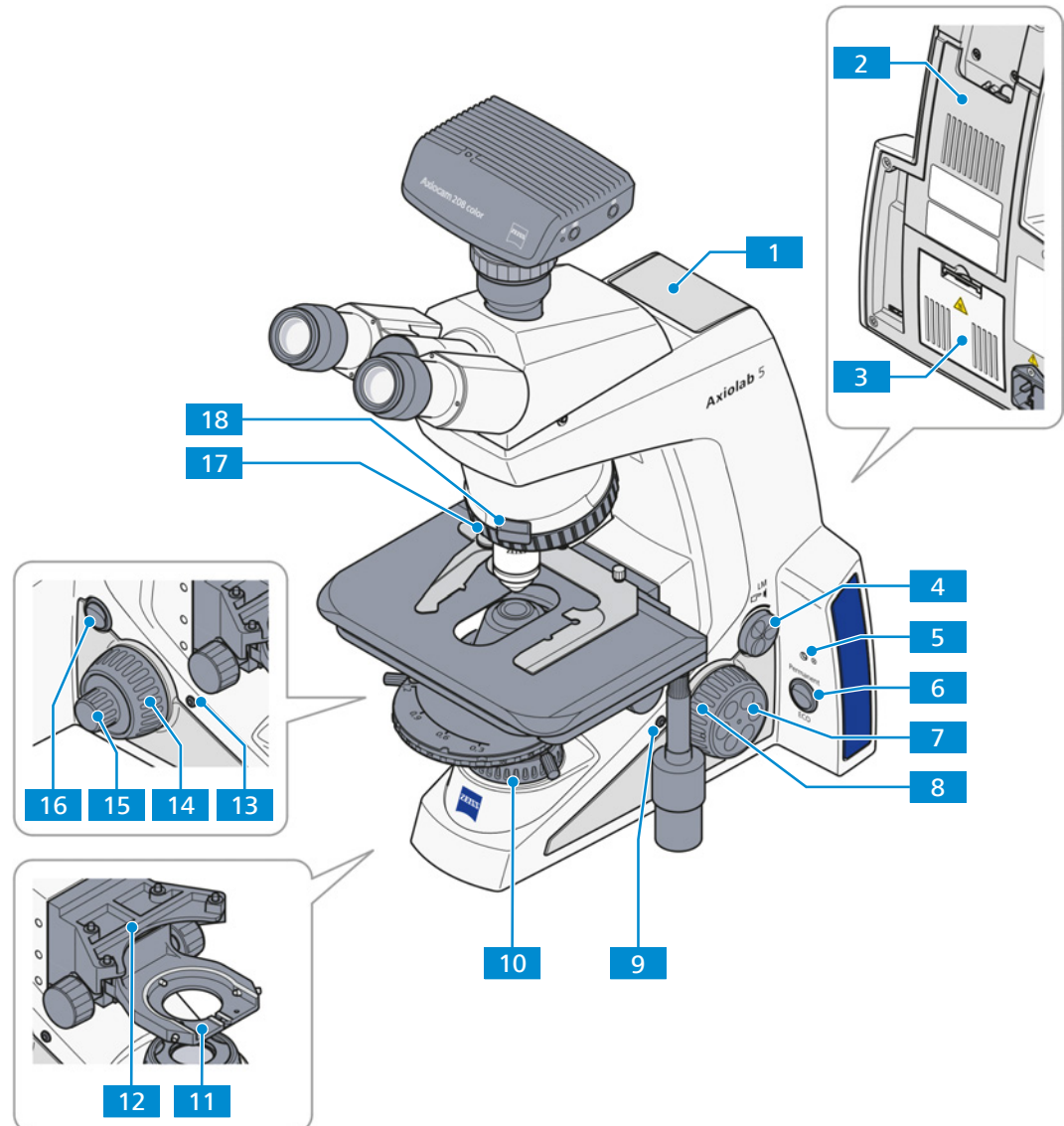


Fig. 8: Functional components and controls - stand Bio-TL

- |  |  |
|--|--|
| <b>1</b> Transport handle  | <b>2</b> Tool kit storage/cable holder                   |
| <b>3</b> Transmitted light light source (inside)                     | <b>4</b> <b>Intensity/LM</b> knob                        |
| <b>5</b> Indicator light   | <b>6</b> <b>Permanent/ECO</b> mode switch                |
| <b>7</b> Focusing drive – fine adjustment (right side, finger wheel) | <b>8</b> Focusing drive – coarse adjustment (right side) |
| <b>9</b> <b>Snap</b> button (right side)                             | <b>10</b> Luminous-field diaphragm                       |
| <b>11</b> <i>Condenser carrier</i> [▶ 39]                            | <b>12</b> Stage carrier for mechanical stages            |
| <b>13</b> <b>Snap</b> button (left side)                             | <b>14</b> Focusing drive – coarse adjustment (left side) |
| <b>15</b> Focusing drive – fine adjustment (left side)               | <b>16</b> Power switch On/Off                            |
| <b>17</b> <i>Nosepiece</i> [▶ 39] with 5 positions BF                | <b>18</b> Slot for 6x20mm slider                         |

### 3.2.2 Axiolab 5 Bio-TL/FL Stand

This section shows the functional elements and controls of stand Bio-TL/FL.

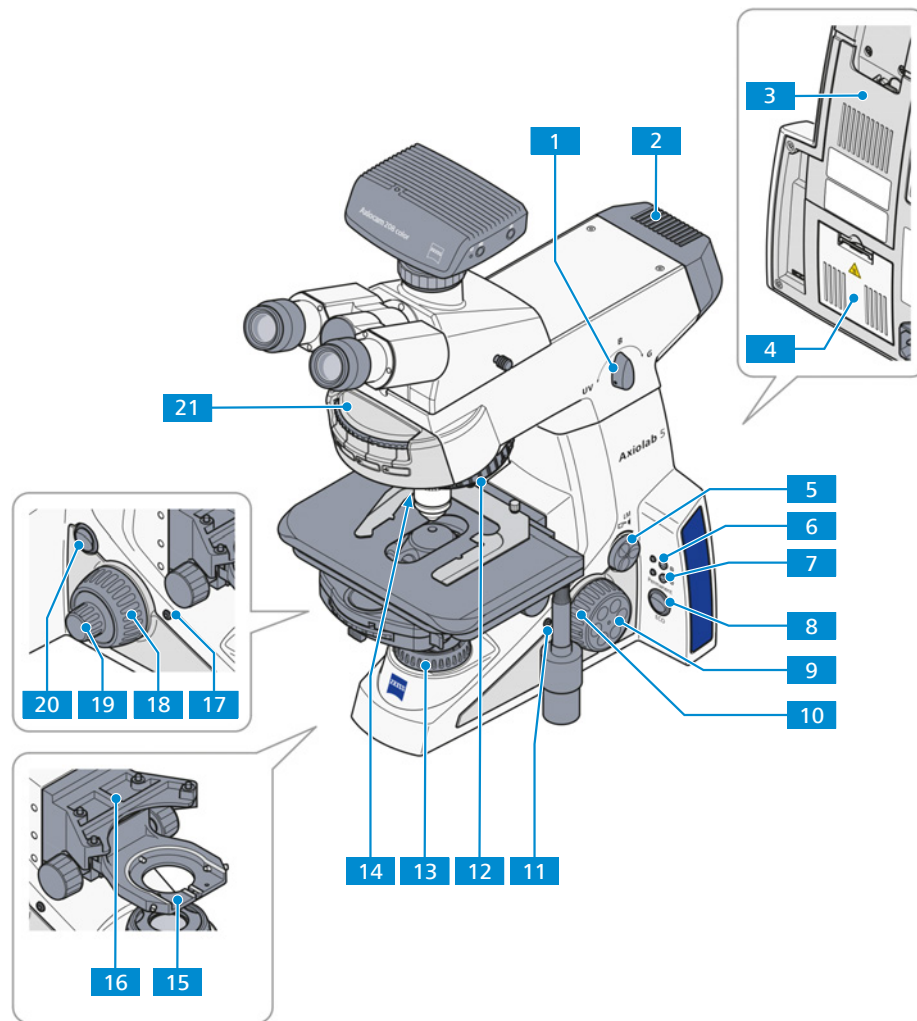


Fig. 9: Functional components and controls - stand Bio-TL/FL

- |   |   |
|---|---|
| <b>1</b> LED selection knob for 3 positions ( <b>UV</b> , <b>B</b> , <b>G</b> )           | <b>2</b> FL-LED reflected light light source  |
| <b>3</b> Tool kit storage/cable holder  | <b>4</b> Transmitted light light source (inside)                                      |
| <b>5</b> <b>Intensity/LM</b> knob   | <b>6</b> Reflected light ( <b>RL</b> ) button and indicator light for reflected light |
| <b>7</b> Transmitted light ( <b>TL</b> ) button and indicator light for transmitted light | <b>8</b> <b>Permanent/ECO</b> mode switch   |
| <b>9</b> Focusing drive – fine adjustment (right side, finger wheel)                      | <b>10</b> Focusing drive – coarse adjustment (right side)                             |
| <b>11</b> <b>Snap</b> button (right side)   | <b>12</b> Nosepiece with 5 positions BF   |
| <b>13</b> Luminous-field diaphragm  | <b>14</b> Slot for 6x20mm slider  |
| <b>15</b> Condenser carrier   | <b>16</b> Stage carrier for mechanical stages   |
| <b>17</b> <b>Snap</b> button (left side)  | <b>18</b> Focusing drive – coarse adjustment (left side)                              |
| <b>19</b> Focusing drive – fine adjustment (left side)                                    | <b>20</b> Power switch On/Off   |
| <b>21</b> 4-position reflector turret   |   |

### 3.2.3 Axiolab 5 Pol-TL Stand

This section shows the functional elements and controls of stand Pol-TL.

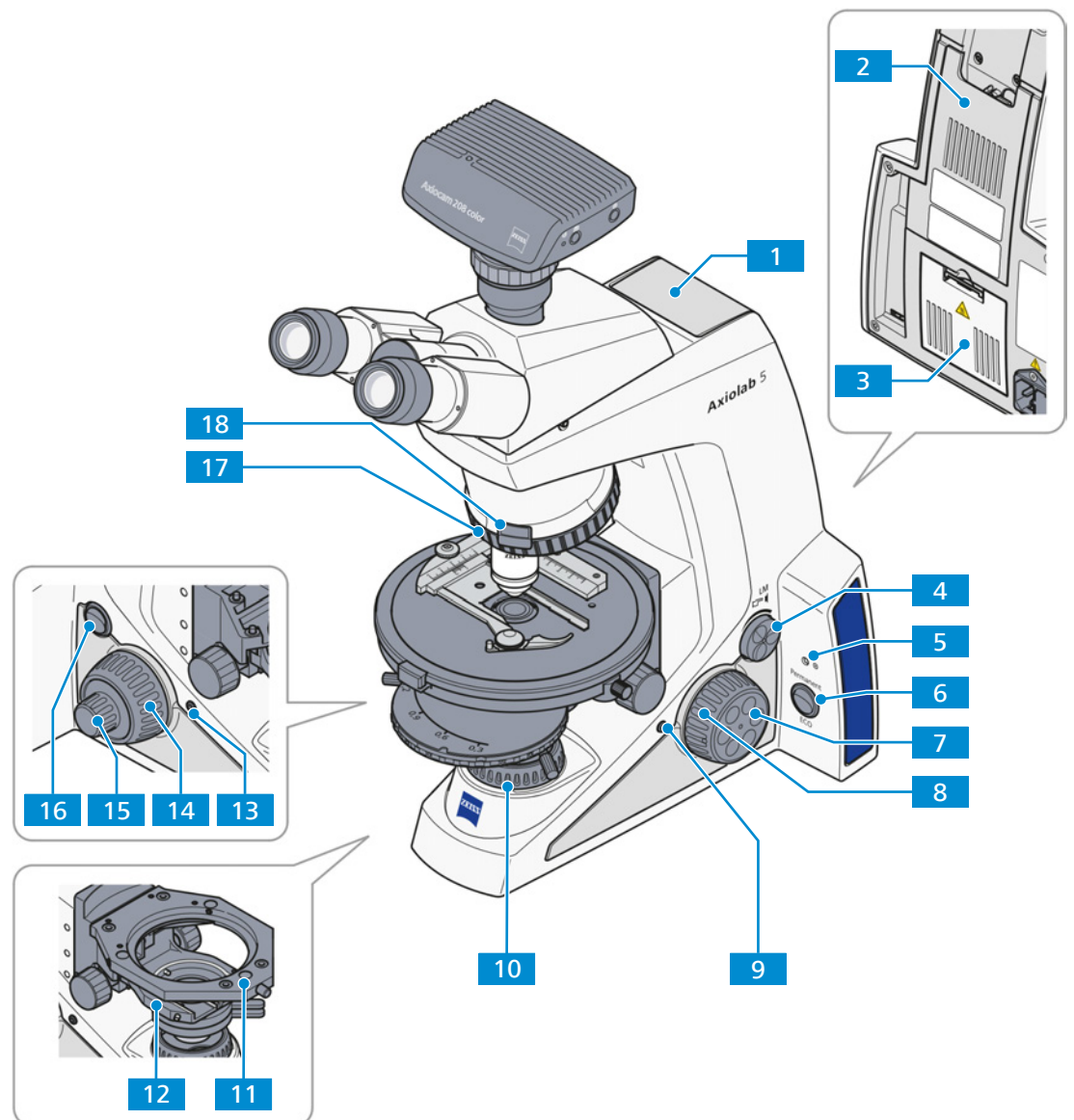


Fig. 10: Functional components and controls - stand Pol-TL

- |  |  |
|--|--|
| <b>1</b> Transport handle  | <b>2</b> Tool kit storage/cable holder                   |
| <b>3</b> Transmitted light light source (inside)                     | <b>4</b> <b>Intensity/LM</b> knob                        |
| <b>5</b> Indicator light   | <b>6</b> <b>Permanent/ECO</b> mode switch                |
| <b>7</b> Focusing drive – fine adjustment (right side, finger wheel) | <b>8</b> Focusing drive – coarse adjustment (right side) |
| <b>9</b> <b>Snap</b> button (right side)                             | <b>10</b> Luminous-field diaphragm                       |
| <b>11</b> Stage carrier for rotary stages                            | <b>12</b> Condenser carrier with mounted polarizer       |
| <b>13</b> <b>Snap</b> button (left side)                             | <b>14</b> Focusing drive – coarse adjustment (left side) |
| <b>15</b> Focusing drive – fine adjustment (left side)               | <b>16</b> Power switch On/Off                            |
| <b>17</b> Nosepiece with 5 positions BF                              | <b>18</b> Slot for 6x20mm slider                         |

### 3.2.4 Axiolab 5 Pol-TL/conoscopy Stand

This section shows the functional elements and controls of stand Pol-TL/conoscopy.

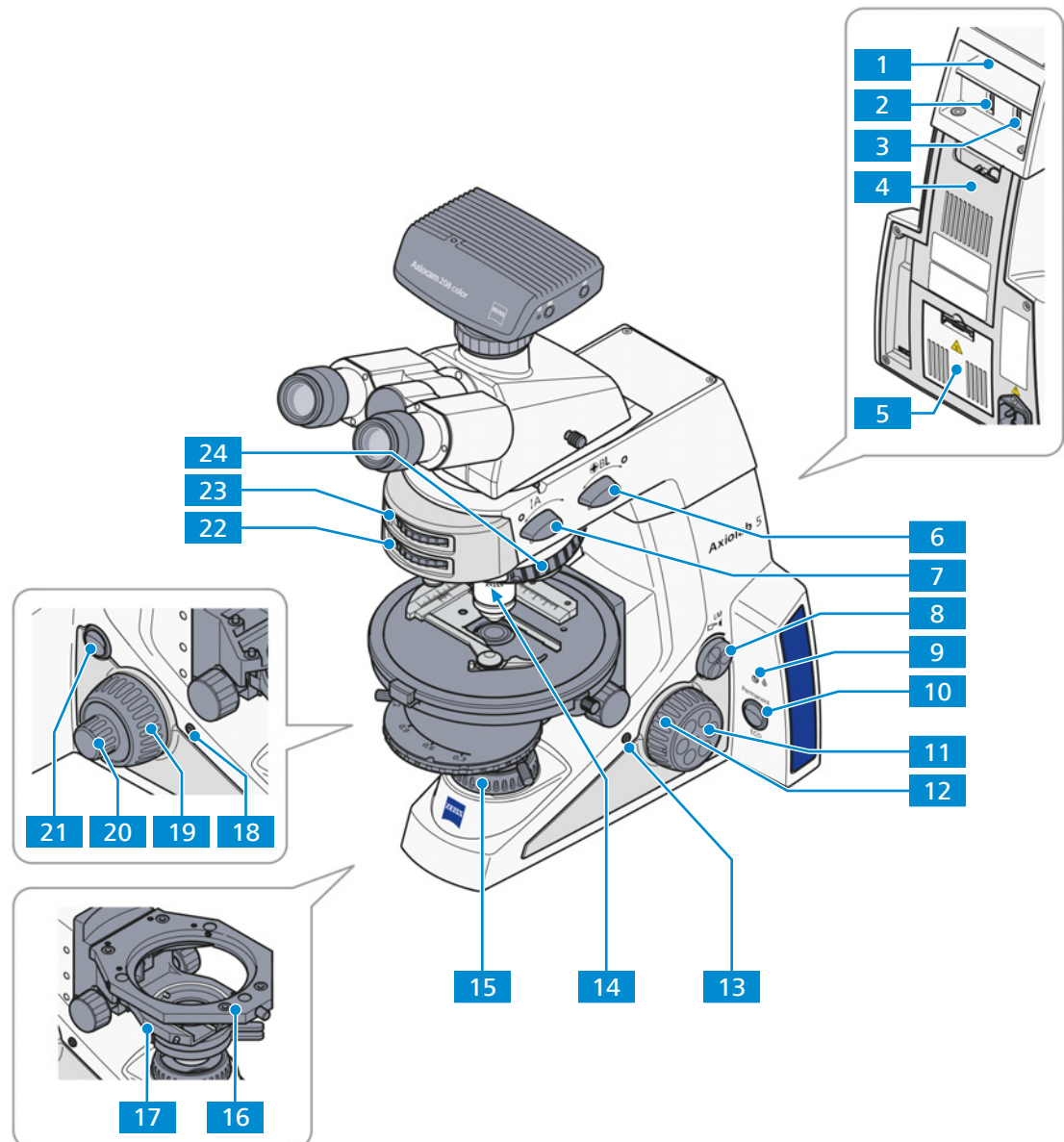


Fig. 11: Functional elements and controls - stand Pol-TL/conoscopy

- |   |   |
|---|---|
| <b>1</b> Transport handle   | <b>2</b> Storage compartment for one 6x20mm slider                              |
| <b>3</b> Storage compartment for one 6x20mm slider                    | <b>4</b> Tool kit storage/cable holder  |
| <b>5</b> Transmitted light light source (inside)                      | <b>6</b> Rotary knob <b>BL</b>  |
| <b>7</b> Rotary knob <b>A</b>   | <b>8</b> <b>Intensity/LM</b> knob   |
| <b>9</b> Indicator light  | <b>10</b> <b>Permanent/ECO</b> mode switch                                      |
| <b>11</b> Focusing drive – fine adjustment (right side, finger wheel) | <b>12</b> Focusing drive – coarse adjustment (right side)                       |
| <b>13</b> <b>Snap</b> button (right side)                             | <b>14</b> Slot for 6x20mm slider  |
| <b>15</b> Luminous-field diaphragm                                    | <b>16</b> Stage carrier for rotary stages (also suitable for mechanical stages) |



- |  |  |
|--|--|
| <b>17</b> Condenser carrier with mounted polarizer       | <b>18</b> Snap button (left side)  |
| <b>19</b> Focusing drive – coarse adjustment (left side) | <b>20</b> Focusing drive – fine adjustment (left side)                     |
| <b>21</b> Power switch On/Off                            | <b>22</b> Setting wheel for polarization direction of analyzer             |
| <b>23</b> Setting wheel for focusing of Bertrand lens    | <b>24</b> Nosepiece with 5 positions BF Pol (with 4x centerable, 1x fixed) |

### 3.2.5 Axiolab 5 Pol-TL/RL Stand

This section shows the functional elements and controls of stand Pol-TL/RL.

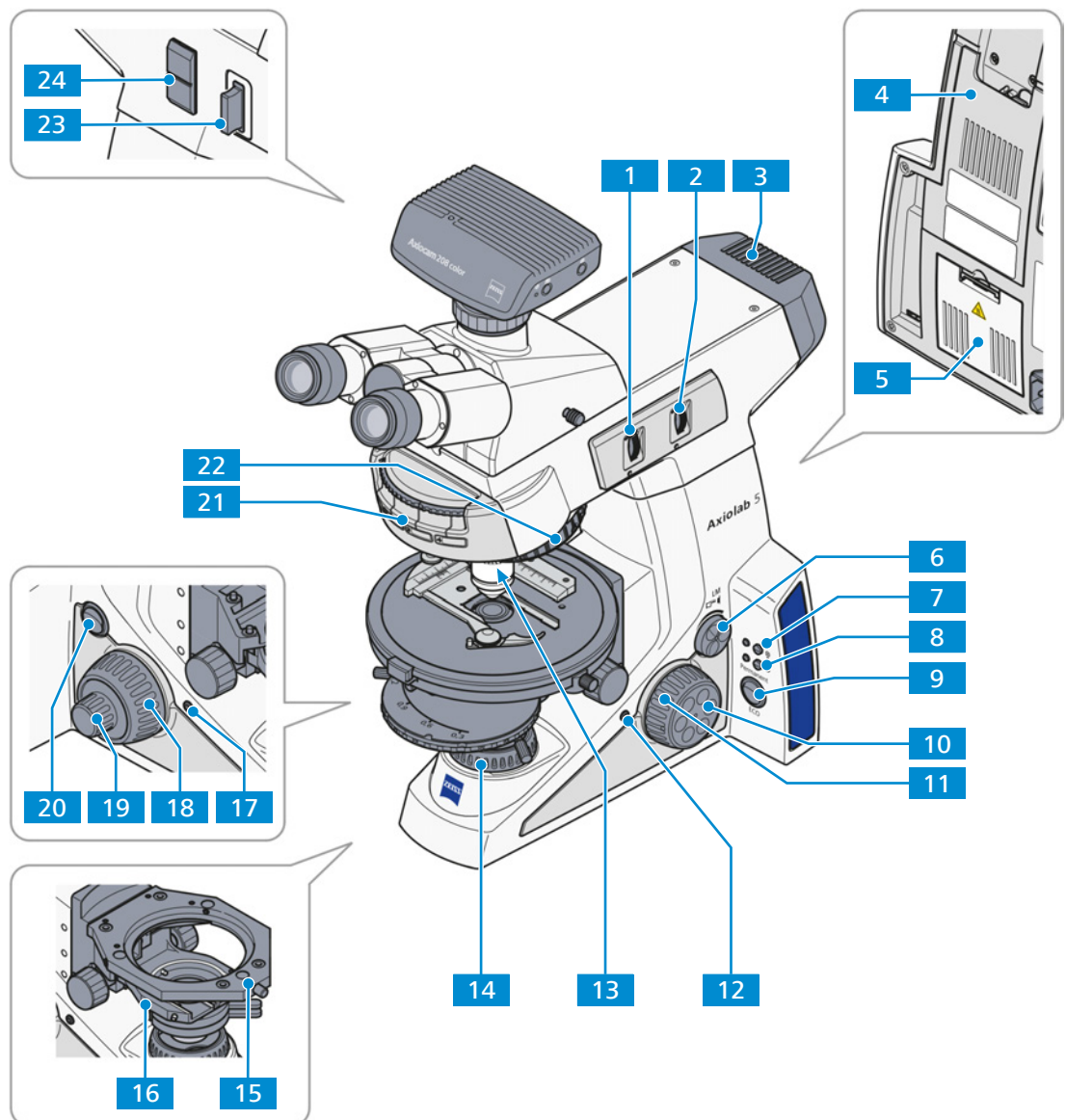


Fig. 12: Functional elements and controls - stand Pol-TL/RL

- |   |   |
|---|---|
| <b>1</b> Luminous-field diaphragm for reflected light (centered)                      | <b>2</b> Aperture diaphragm for reflected light (centered)                                |
| <b>3</b> Reflected light light source   | <b>4</b> Tool kit storage/cable holder  |
| <b>5</b> Transmitted light light source (inside)                                      | <b>6</b> <b>Intensity/LM</b> knob   |
| <b>7</b> Reflected light ( <b>RL</b> ) button and indicator light for reflected light | <b>8</b> Transmitted light ( <b>TL</b> ) button and indicator light for transmitted light |
| <b>9</b> <b>Permanent/ECO</b> mode switch   | <b>10</b> Focusing drive – fine adjustment (right side, finger wheel)                     |
| <b>11</b> Focusing drive – coarse adjustment (right side)                             | <b>12</b> <b>Snap</b> button (right side)   |
| <b>13</b> Slot for 6x20mm slider  | <b>14</b> Luminous-field diaphragm  |
| <b>15</b> Stage carrier for rotary stages (also suitable for mechanical stages)       | <b>16</b> Condenser carrier with mounted polarizer  |
| <b>17</b> <b>Snap</b> button (left side)  | <b>18</b> Focusing drive – coarse adjustment (left side)                                  |
| <b>19</b> Focusing drive – fine adjustment (left side)                                | <b>20</b> Power switch On/Off   |
| <b>21</b> 4-position reflector turret   | <b>22</b> Nosepiece, with 5 positions BF DF Pol   |
| <b>23</b> Slot for polarizer slider 6x30mm, reflected light (slider inserted)         | <b>24</b> Slot for 2-position filter slider, reflected light (covered)                    |

### 3.2.6 Axiolab 5 Mat-TL/RL Stand

This section shows the functional elements and controls of stand Mat-TL/RL.

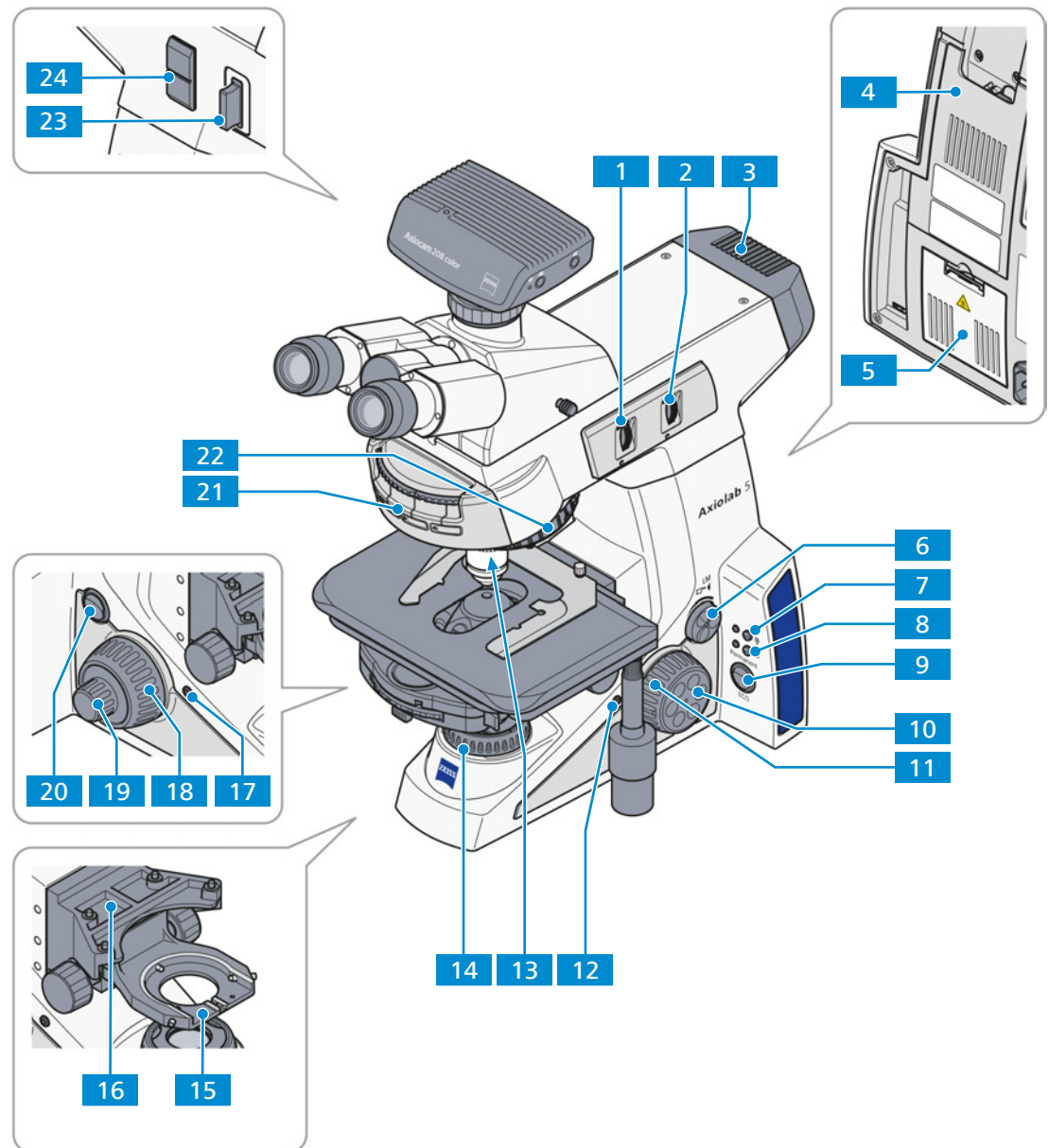


Fig. 13: Functional elements and controls - stand Mat-TL/RL

- |   |   |
|---|---|
| <b>1</b> Luminous-field diaphragm for reflected light (centered)                      | <b>2</b> Aperture diaphragm for reflected light (centered)                                |
| <b>3</b> Reflected light light source   | <b>4</b> Tool kit storage/cable holder  |
| <b>5</b> Transmitted light light source (inside)                                      | <b>6</b> <b>Intensity/LM</b> knob   |
| <b>7</b> Reflected light ( <b>RL</b> ) button and indicator light for reflected light | <b>8</b> Transmitted light ( <b>TL</b> ) button and indicator light for transmitted light |
| <b>9</b> <b>Permanent/ECO</b> mode switch   | <b>10</b> Focusing drive – fine adjustment (right side, finger wheel)                     |
| <b>11</b> Focusing drive – coarse adjustment (right side)                             | <b>12</b> <b>Snap</b> button (right side)   |
| <b>13</b> Slot for 6x20mm slider  | <b>14</b> Luminous-field diaphragm  |

<b>15</b>	Condenser carrier	<b>16</b>	Stage carrier for mechanical stages
<b>17</b>	<b>Snap</b> button (left side)	<b>18</b>	Focusing drive – coarse adjustment (left side)
<b>19</b>	Focusing drive – fine adjustment (left side)	<b>20</b>	Power switch On/Off
<b>21</b>	4-position reflector turret	<b>22</b>	Nosepiece, with 5 positions BF DF
<b>23</b>	Slot for polarizer slider 6x30mm, reflected light (slider inserted)	<b>24</b>	Slot for filter slider, reflected light (covered)

### 3.2.7 Functions of Stand Keys and Display Elements

Key	Action	Functionality/Description
Power switch On/Off	<b>I</b> = on; <b>O</b> = off	Switches the microscope on/off.
<b>Permanent/ECO</b> mode switch	Toggle	<p>Switches between Permanent (continuous) mode and ECO mode of the microscope illumination.</p> <ul style="list-style-type: none"> <li>Permanent mode active: illumination is continuously switched on.</li> <li>ECO mode active: illumination switches off after 15 minutes without action.</li> </ul> <p>Don't use ECO mode for experiments involving time-lapse or video recording.</p>
Indicator light	Blinking***	Indicates whether microscope is working in TL or RL mode.
<b>RL</b> button, <b>TL</b> button	Short press*	<p>Switches RL/TL illumination alternately on/off.</p> <p>The respective indicator light is continuously illuminated.</p>
<b>Intensity/LM</b> knob	Turn	Controls the light intensity of the active light source.
	Long press**	<p>Light Manager function:</p> <p>Saves the light intensity; after saving is done, the illumination switches off for 300 ms (darkness indicates action to user).</p>
	Long press for 20 s	<p>Activates the factory default settings (enables/disables Light Manager (LM) functionality).</p> <p>The indicator light starts blinking in RED after 3 s until 20 s is reached. After 20 s, the indicator light turns to GREEN continuously.</p>



Key	Action	Functionality/Description
Left <b>Snap</b> button or right <b>Snap</b> button (only if Ax-iocam 202 or 208 is installed)	Short press*	Snaps an image; after snap is finished, the attached monitor appears in BLACK for 50 ms.
	Long press**	Starts video recording; another short press is required to stop recording. After recording is finished, the attached monitor appears in BLACK for 300 ms.
<b>Snap</b> button + <b>Intensity/LM</b> knob	Long press** (simultaneously)	Enables/disables the Light Manager (LM) functionality: <ul style="list-style-type: none"> <li>Disabling: The indicator light blinks GREEN / ORANGE / GREEN in sequence.</li> <li>Enabling: The indicator light blinks GREEN / GREEN / GREEN in sequence.</li> </ul>
Left <b>Snap</b> button	press > 1.5 s	Disables/enables the dazzle protection function: <ul style="list-style-type: none"> <li>Disabling: The indicator light blinks ORANGE twice.</li> <li>Enabling: The indicator light blinks GREEN twice.</li> </ul> <p>By factory default, the dazzle protection function is enabled.</p>

\* Short press means: hold less than 1 second, then release.

\*\* Long press means: hold at least 1.5 seconds.

\*\*\* Blinking: the indicator light alternately goes on/off at 500 ms intervals

### 3.2.8 Binocular Tube

Various tubes with different inclination angles enable suitable eye levels to be selected for observation.

#### 3.2.8.1 Binocular Photo Tube 30°/23 (50:50)

**Purpose** Binocular photo tubes are used to visualize the microscopic image by means of the eyepieces. Using the camera port, the microscopic image can be sent to a connected camera.

**Position** The binocular photo tubes are mounted on the top of the stand.

**Function** The pupil distance and the viewing height can be adjusted with the aid of the binocular section. The following features and controls are available:

- reversed image
- camera port with fixed light graduation (50:50)
- viewing angle 30°
- field of view 23 mm

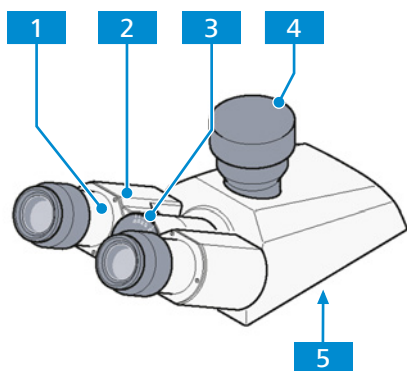


Fig. 14: Binocular Photo Tube 30°/23 (50:50)

- |                              |                                |
|------------------------------|--------------------------------|
| <b>1</b> Eyepiece socket     | <b>2</b> Binocular section     |
| <b>3</b> Angle scale         | <b>4</b> Camera port (covered) |
| <b>5</b> Dovetail ring mount |                                |

### 3.2.8.2 Binocular Photo Tube 30°/23 (100:0/0:100)

**Purpose** Binocular photo tubes are used to visualize the microscopic image by means of the eyepieces. Using the camera port, the microscopic image can be sent to a connected camera.

**Position** The binocular photo tubes are mounted on the top of the stand.

**Function** The pupil distance and the viewing height can be adjusted with the aid of the binocular section. The following features and controls are available:

- reversed image
- camera port with toggle graduation (100:0/0:100)
- viewing angle 30°
- eyepiece shutter
- field of view 23 mm

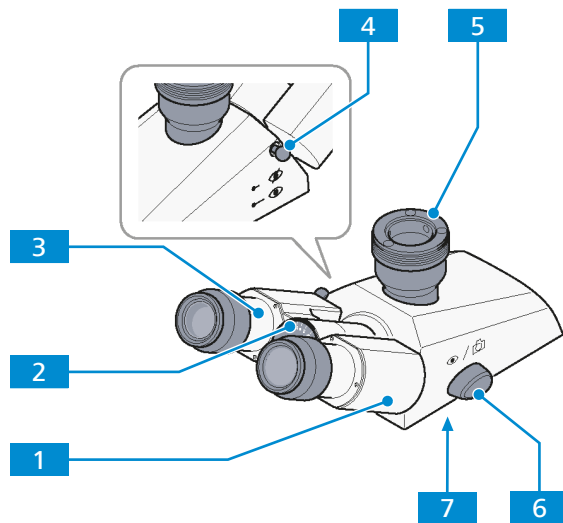


Fig. 15: Binocular Photo Tube 30°/23 (100:0/0:100)

- |   |  |
|---|--|
| <p><b>1</b> Binocular section</p> <p><b>3</b> Eyepiece socket</p> <p><b>5</b> Camera port</p> <p><b>7</b> Dovetail ring mount</p> | <p><b>2</b> Angle scale</p> <p><b>4</b> Eyepiece shutter</p> <ul style="list-style-type: none"> <li>▪ Push-pull rod pushed in: eyepiece shutter closed</li> <li>▪ Push-pull rod pulled out: eyepiece shutter opened</li> </ul> <p><b>6</b> Shift knob for selecting the graduation</p> <ul style="list-style-type: none"> <li>▪ Shift knob to front (eye symbol): 100% light to eyepieces</li> <li>▪ Shift knob to the rear (camera symbol): 100% light to camera</li> </ul> |
|---|--|

### 3.2.8.3 Binocular Photo Tube 20°/23 (100:0/0:100)

**Purpose** Binocular photo tubes are used to visualize the microscopic image by means of the eyepieces. Using the camera port, the microscopic image can be sent to a connected camera.

**Position** The binocular photo tubes are mounted on the top of the stand.

**Function** The pupil distance and the viewing height can be adjusted with the aid of the binocular section. The following features and controls are available:

- upright image
- camera port with toggle graduation (100:0/0:100)
- viewing angle 20°
- field of view 23 mm

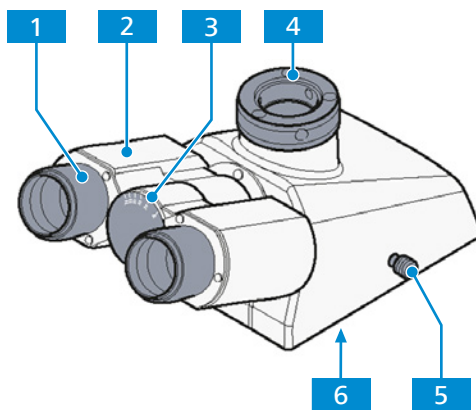


Fig. 16: Binocular Photo Tube 20°/23 (100:0/0:100)

- |  |                              |
|--|------------------------------|
| <b>1</b> Eyepiece socket                     | <b>2</b> Binocular section   |
| <b>3</b> Angle scale                         | <b>4</b> Camera port         |
| <b>5</b> Slider for selecting the graduation | <b>6</b> Dovetail ring mount |
- Slider pushed in: 100% light to eyepieces
  - Slider pulled out: 100% light to camera. 100% light to camera

### 3.2.8.4 Binocular Ergo Photo Tube -2° to 28°/23 (50:50), Reversed Image

**Purpose** Binocular photo tubes provide users with the convenience of accessing the sample images using eyepieces as well as monitors through camera at the same time. Users can change the angle from -2° to 28° to meet the ergonomic requirement of specific operators.

**Position** The binocular photo tubes are mounted on the top of the stand.

**Function** The pupil distance and viewing height can be adjusted by bending the binocular device up or down.

The following features and controls are available:

- Reverse image
- Camera port with splitting ratio 50:50
- Viewing angle -2° to 28°
- Field of view 23 mm

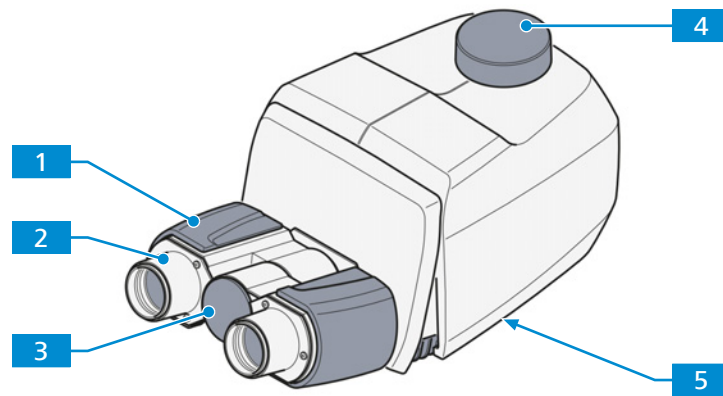


Fig. 17: Binocular ergo photo tube -2° to 28°/23 (50:50), reversed image

- |                              |                                |
|------------------------------|--------------------------------|
| <b>1</b> Binocular section   | <b>2</b> Eyepiece socket       |
| <b>3</b> Angle scale         | <b>4</b> Camera port (covered) |
| <b>5</b> Dovetail ring mount |                                |

### 3.2.8.5 Binocular Ergo Photo Tube -2° to 28°/25 (50:50), Reversed Image

**Purpose** Binocular photo tubes provide users with the convenience of accessing the sample images using eyepieces as well as monitors through camera at the same time. Users can change the angle from -2° to 28° to meet the ergonomic requirement of specific operators.

**Position** The binocular photo tubes are mounted on the top of the stand.

**Function** The pupil distance and viewing height can be adjusted by bending the binocular device up or down.

The following features and controls are available:

- Reverse image
- Camera port with splitting ratio 50:50
- Viewing angle -2° to 28°
- Field of view 25 mm

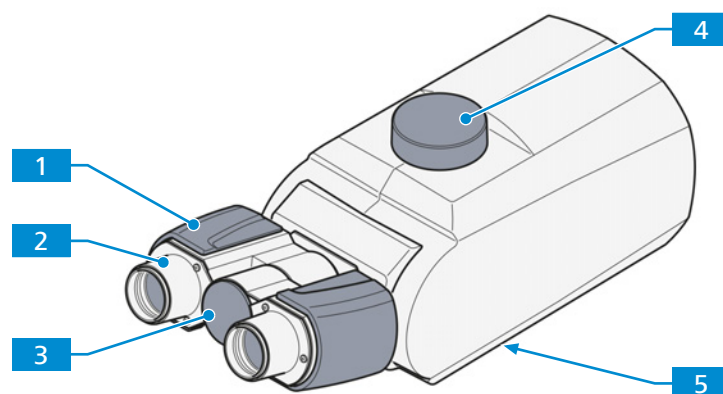


Fig. 18: Binocular ergo photo tube -2° to 28°/25 (50:50), reversed image

- |                              |                                |
|------------------------------|--------------------------------|
| <b>1</b> Binocular section   | <b>2</b> Eyepiece socket       |
| <b>3</b> Angle scale         | <b>4</b> Camera port (covered) |
| <b>5</b> Dovetail ring mount |                                |

### 3.2.9 Eyepieces

- Purpose** The eyepieces serve to observe the microscopic image.
- Position** The eyepieces are inserted into the eyepiece sockets of the binocular tube.
- Function** Both eyepieces are suitable for spectacle wearers. Additionally, they contain a focusing ring for compensation of defective vision. The provided diopter scale helps to find the correct setting. When using the microscope for fluorescence applications, the special eyecups with light protection can be used. However, they cannot be folded over and are not suitable for spectacle wearers.

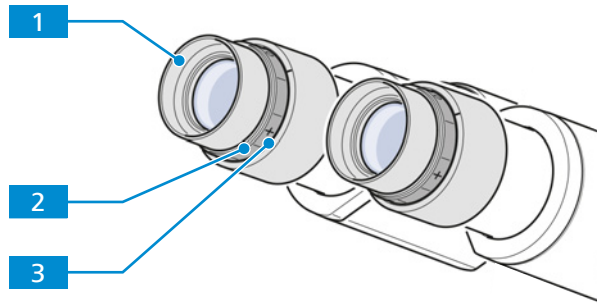


Fig. 19: Eyepiece

- |   |                        |
|---|------------------------|
| <b>1</b> Eyecup (e.g. Foldover rubber eyecup) | <b>2</b> Focusing ring |
| <b>3</b> Diopter scale                        |                        |

### 3.2.10 Eyepieces with Eyepiece Reticles

- Purpose** The eyepieces with eyepiece reticles serve to observe the microscopic image in special microscopy procedures.
- Position** The eyepieces with reticles are inserted into the tube.
- The eyepiece reticles must be inserted under dust-free conditions. This should be carried out only by ZEISS Service.

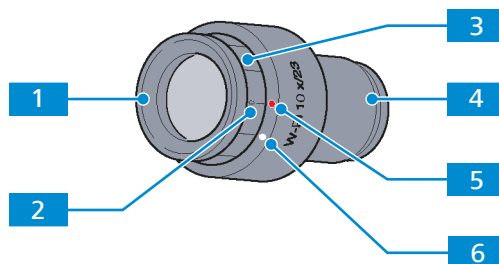


Fig. 20: Eyepiece with installed eyepiece reticle

- |  |   |
|--|---|
| <b>1</b> Eyecup, changeable  | <b>2</b> Diopter scale with zero point to facilitate finding the correct setting        |
| <b>3</b> Focusing ring for compensation of defective vision                          | <b>4</b> Mounting stop with inserted eyepiece reticle                                   |
| <b>5</b> Red dot, corresponds to the zero diopter setting when a reticle is inserted | <b>6</b> White dot, corresponds to the zero diopter setting when no reticle is inserted |

### 3.2.11 Nosepiece with Objectives

**Purpose** The nosepiece is used to hold the objectives and to swivel the desired objective into the beam path.

**Position** The nosepiece is mounted on the upper part of the stand.

The following features and controls are available:

- nosepiece with M27 mounting thread for five objectives
- one objective position is fixed and four positions can be centered with the aid of two screws each
- equipped with slot for 6x20mm sliders (compensators, analyzers, quarter plates or fluorescence protection shield)

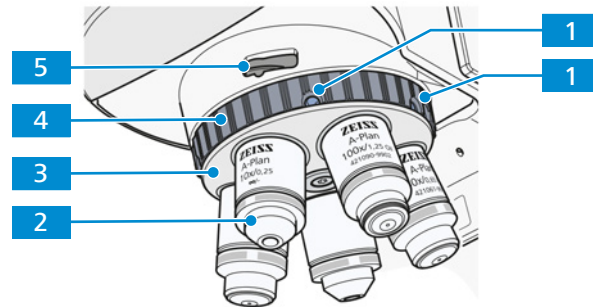


Fig. 21: Nosepiece with objectives

- |          |   |          |  |
|----------|---|----------|--|
| <b>1</b> | Centering screw, two for each centerable objective position | <b>2</b> | Objective                                  |
| <b>3</b> | Nosepiece   | <b>4</b> | Knurled ring for swivelling the nose-piece |
| <b>5</b> | Slot 6x20mm   |          |  |

### 3.2.12 Condenser Carrier

**Purpose** The condenser carrier is used to hold the condenser.

**Position** The condenser carrier is mounted onto the stage carrier.

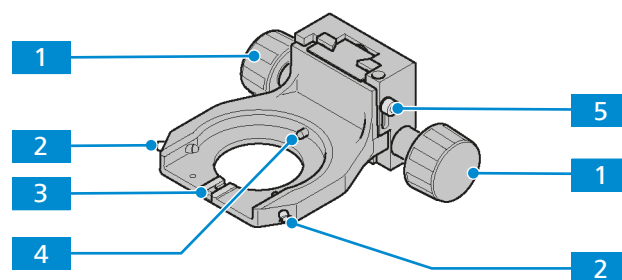


Fig. 22: Condenser carrier

- |          |   |          |   |
|----------|---|----------|---|
| <b>1</b> | Knurled knob for vertical adjustment (left/right) | <b>2</b> | Centering hex screw (left/right), optional: knurled screw |
| <b>3</b> | Orientation groove                                | <b>4</b> | Main spring   |
| <b>5</b> | Fastening screw for height stop                   |          |   |

### 3.2.13 Condensers

#### 3.2.13.1 Condenser 0.9/1.25 BF

**Purpose** Condensers are used to optimize the transmitted light illumination. The condenser 0.9/1.25 BF is usable for brightfield applications.

**Position** The condenser is mounted on the condenser carrier of the stand.

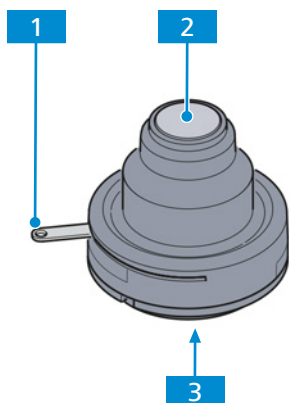


Fig. 23: Condenser 0.9/1.25 BF

- |          |  |          |            |
|----------|--|----------|------------|
| <b>1</b> | Lever for setting the aperture diaphragm | <b>2</b> | Front lens |
| <b>3</b> | Dovetail ring mount                      |          |            |

#### 3.2.13.2 Condenser 0.9/1.25 BF, DF, Ph1, Ph2, Ph3 with modulator disk

**Purpose** Condensers are used to optimize the transmitted light illumination. The condenser with modulator disk is usable for brightfield, darkfield and phase contrast applications.

**Position** The condenser is mounted on the condenser carrier of the stand.

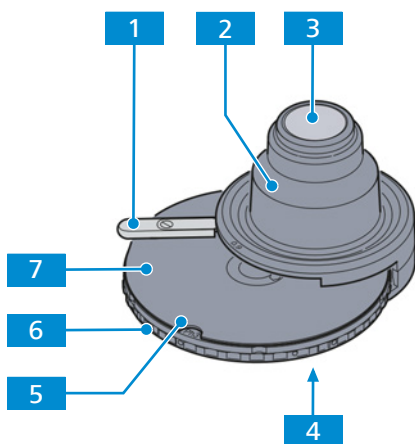


Fig. 24: Condenser 0.9/1.25 BF, DF, Ph1, Ph2, Ph3 with modulator disk

- |          |   |          |   |
|----------|---|----------|---|
| <b>1</b> | Lever for setting the aperture diaphragm                    | <b>2</b> | Front lens  |
| <b>3</b> | Condenser 0.9/1.25 BF, optionally condenser 0.9/1.25 BF Pol | <b>4</b> | Dovetail ring mount   |
| <b>5</b> | Display field of the adjusted modulator disk position       | <b>6</b> | Knurled ring for adjusting the position of the modulator disk |
| <b>7</b> | Modulator disk with 5 positions for condenser modules       |          |   |



### 3.2.13.3 Condenser 0.8 BF WD=5.8mm

**Purpose** Condensers are used to optimize the transmitted light illumination. The condenser 0.8 BF is usable for brightfield applications.

**Position** The condenser is mounted on the condenser carrier of the stand.

The following features and controls are available:

- NA of 0.8 in air and WD of 5.8 mm
- Suitable for brightfield observation with objectives 5x to 100x
- Compatible with low-power system for objectives 2.5x/4x

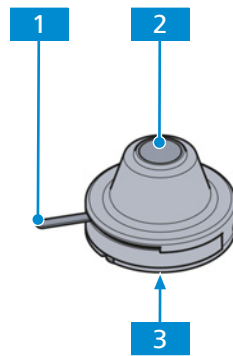


Fig. 25: Condenser 0.8 BF WD=5.8mm

**1** Lever for setting the aperture diaphragm

**2** Front lens

**3** Dovetail ring mount

### 3.2.14 Stages

A stage is a platform at right angles to the optical axis of the microscope, which carries the sample and which is often fitted with mechanical movements (as in a mechanical stage) to allow easy positioning of the object in the x- and y-axis, and movement along, and rotation about the z-axis.

#### 3.2.14.1 Rackless Mechanical Stage, 75x50 R

**Purpose** Mechanical stages are used for fixing and positioning the sample for examination.

**Position** The mechanical stages are mounted on the stage carrier of the stand.

**Function** The sample is fixed on the stage by means of the sample holder. For this purpose, the sample holder is equipped with a spring lever.

The sample is positioned in the beam path by means of the two coaxial drives in X and Y direction. The adjustment range can be read off the respective vernier scale.

The following features and controls are available:

- rackless stage
- coaxial drives in X and Y adjustment on the right (R), optionally on the left (L)
- travel range 75x50mm
- with hardcoat anodized surface

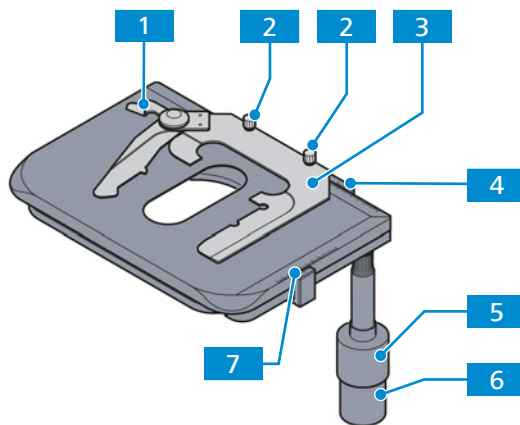


Fig. 26: Rackless mechanical stage, 75x50 R

**1** Spring lever

**3** sample holder for double slides 76x26

**5** Coaxial knurled knob for Y adjustment

**7** Vernier scale for display of the adjustment range in Y

**2** Knurled screw (2x) for fixing the sample holder to the stage

**4** Vernier scale for display of the adjustment range in X

**6** Coaxial knurled knob for X adjustment

### 3.2.14.2 Rotary Stage Pol 360° with Clamping Device

**Purpose** Rotary stages are used for fixing and positioning the sample for examination in polarized light.

**Position** The rotary stages are mounted on the stage carrier of the stand.

**Function** The sample is fixed on the stage by means of the clamping device.

The following features and controls are available:

- 360° rotation with lock
- click stop every 45°; enabled or disabled via the control knob

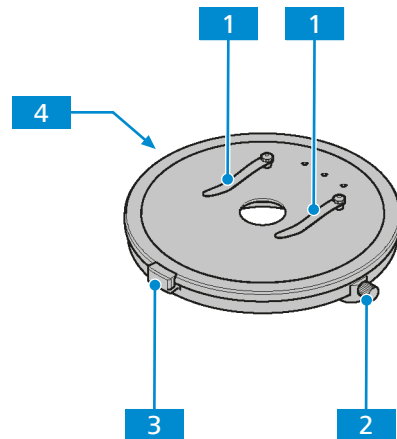


Fig. 27: Rotary Stage Pol 360° with clamping device

- |  |   |
|--|---|
| <p><b>1</b> Clamping device</p> <p><b>3</b> Click stop every 45°</p> | <p><b>2</b> Knurled screw for locking the rotation, 360° rotation possible</p> <p><b>4</b> Control knob to enable/disable the clip stop function (on the bottom side)</p> |
|--|---|

### 3.2.15 Reflector Turret

**Purpose** The reflector turret is used to hold the push-and-click (P&C) reflector modules and to swivel the desired reflector module into the beam path.

**Position** The reflector turret is mounted on the upper part of the stand above the nosepiece.

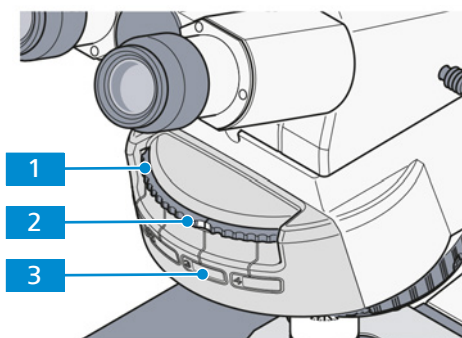


Fig. 28: 4-positions reflector turret

- |  |   |
|--|---|
| <p><b>1</b> Knurled ring to swivel the desired reflector module into the beam path</p> <p><b>3</b> Field for the supplied stickers, the stickers can be labeled with the filter combination data of the reflector module and pasted on the corresponding field</p> | <p><b>2</b> Marking on the knurled ring, indicates which reflector module is in the beam path</p> |
|--|---|

### 3.2.16 Storage Compartment

**Purpose** The storage compartment is used to store tools and the power cord (rolled up for transport).

**Position** The storage compartment is located at the back of the stand.

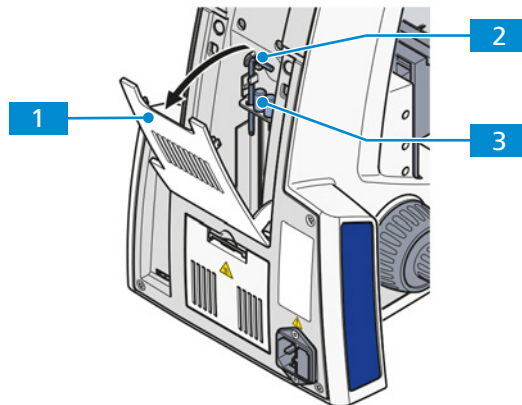


Fig. 29: Storage compartment

**1** Cover flap

**2** Angled hex key (3 mm)

**3** Two knurled hex keys (1.5 mm)

## 3.3 Light Manager Function

The Light Manager (LM) function allows users to save the favorable light intensity of specific objective and turret combinations. Afterwards, when switching between different objectives or reflector turrets, the saved light intensity will be automatically applied. Restarting the microscope will not vanish the saved value.

This value can be overwritten among different users with unlimited times.

Dazzle protection (ON by default) is used to switch off the light when switching objectives or reflector turrets such that the field of view is dark after first objective or turret leaves the position and before the second arrives in. Users can manually turn it off, so the field of view remains the brightness of the first objective or turret until the second one arrives in the position.

## 3.4 Microscopy and Contrast Techniques

The availability of microscopy and contrast techniques depends on the microscope type and the configuration.

### 3.4.1 Transmitted Light Brightfield Microscopy Using the KÖHLER Method

Transmitted light brightfield microscopy is the most common of all optical microscopy methods, since it can be used to quickly and easily examine high-contrast or stained samples (e.g. blood smears).

In order to obtain an image as close as possible to the object, not only the so-called direct beam bundles but also the indirect ones, i.e. the beam bundles diffracted and scattered at the preparation details, are of essential importance. According to ABBE, the larger the indirect beam components are, the more true to the object the microscopic image is.

The best performance of the microscope, and especially its objective, is achieved when the condenser, field diaphragm and aperture diaphragm are adjusted in accordance with the KÖHLER illumination principles.

### 3.4.2 Transmitted Light Darkfield Microscopy Using the KÖHLER Method

In the transmitted light darkfield microscopy you basically illuminate the sample with an illumination aperture which is higher than the one of the objective you are using.

In darkfield microscopy, only the diffracted and scattered light portions which are important for the imaging procedure get into the objective, whereas the indirect unaffected light beams are directed past the objective. Thus a resolution of fine structures can be achieved which is below the resolution capacity of a light microscope. The fine structures now appear bright and incandescent on a dark background.

Darkfield samples need to be kept impeccably clean, more so than samples for any other method. A fingerprint, dust or any dirt particle can have a negative effect, as they brighten the background and reduce the contrast of the object image.

### 3.4.3 Transmitted Light Phase Contrast Microscopy

The phase contrast method is ideal for examining thin uncolored samples, e.g. individual cells of cell cultures. Generally, the human eye cannot detect phase differences (variations in refractive index or thickness) within the different cell components.

The phase contrast method uses the optical modulators "annular phase diaphragm" and "phase ring" to convert the small phase differences in intensity differences which are visible to the human eye. The interference of different beams in the intermediate image is important for the generation of such images.

With the aid of the optically defined ring channel "annular phase diaphragm and phase ring", the bright direct light portions are attenuated and provided with a constant phase shift. The indirect light portions, however, which are diffracted by different cell particles, bypass this optical channel and their phase is affected by the difference in the sample's refractive index and thickness.

In the intermediate image plane, the partial beams are thus differently affected and achieve interference and strengthen or weaken each other (constructive and destructive interference) – depending on their phase. As a result, these interferences create image contents with intensity differences visible to the human eye.

### 3.4.4 Transmitted Light Polarization

The transmitted light polarization method is used for samples which change the polarization of the light. Such samples are called birefringent. Examples include crystals, minerals or polymers. If such birefringent substances are observed between crossed polarizers, the birefringent portion of the sample appears bright while its surroundings remain dark.

#### 3.4.4.1 Detecting Birefringence

A birefringent substance can be recognized by turning the sample by 360° between crossed polarizers. The sample should show four bright and four dark appearances during the turning procedure. During the turning procedure, interference colors appear that range from gray (mostly for biological samples) through white, yellow and red until blue, depending on birefringence, thickness as well as sample orientation. The interference colors may be of the first or of a higher order.

#### 3.4.4.2 Determination of the Polarization Direction

The determination of the polarization direction of  $n_v$  or  $n_v'$  respectively (polarization direction with the absolute or relative largest index of refraction) and  $n_a$  or  $n_a'$  respectively (polarization direction with the absolute or relative smallest index of refraction) relative to the morphological directions, e.g. of crystal surfaces, crystal needles or fibers, provide an important signature of the material. This method is also used in the diagnosis of bio-crystals (e.g. gout and pseudo-gout).

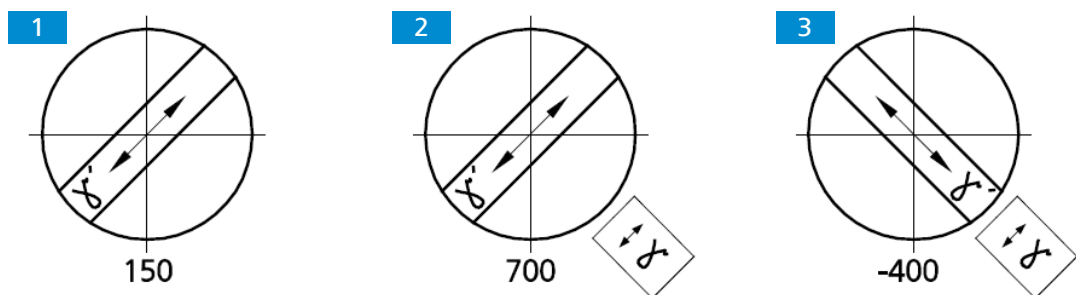


Fig. 30: Determining the polarization direction  $n_v$ , using a synthetic fiber as an example

When the lambda compensator is put in, the sample changes its color depending on its orientation (northeast-southwest or northwest-southeast). Like the sample, the lambda compensator is a birefringent object, but it has a defined path difference of 550 nm and a maximum oscillation direction  $n_v$  pointing strongly to northeast-southwest.

The changes in color are based on optical interference. It is necessary to compare the interference colors (path differences) in both diagonal positions (northeast-southwest and northwest-southeast).

The path difference results from the interference of the polarization of the sample and the polarization of the lambda compensator.

The largest path difference occurs when the polarization direction of the sample or the absolute or relative largest index of refraction ( $n_v$  or  $n_v'$ ) is parallel to the largest polarization direction of the lambda compensator. The sample appears then e.g. in blue-green **2**.

The smallest path difference occurs when the direction of polarization of the sample with the absolute or relative smallest index of refraction ( $n_a$  or  $n_a'$ ) is perpendicular to the polarization direction of the lambda compensator. The sample then appears e.g. yellow **3**.

The gray-white color appearing first in the bright position in the above example **1** corresponds to a path difference of 150 nm according to the Michel-Lévy color chart).

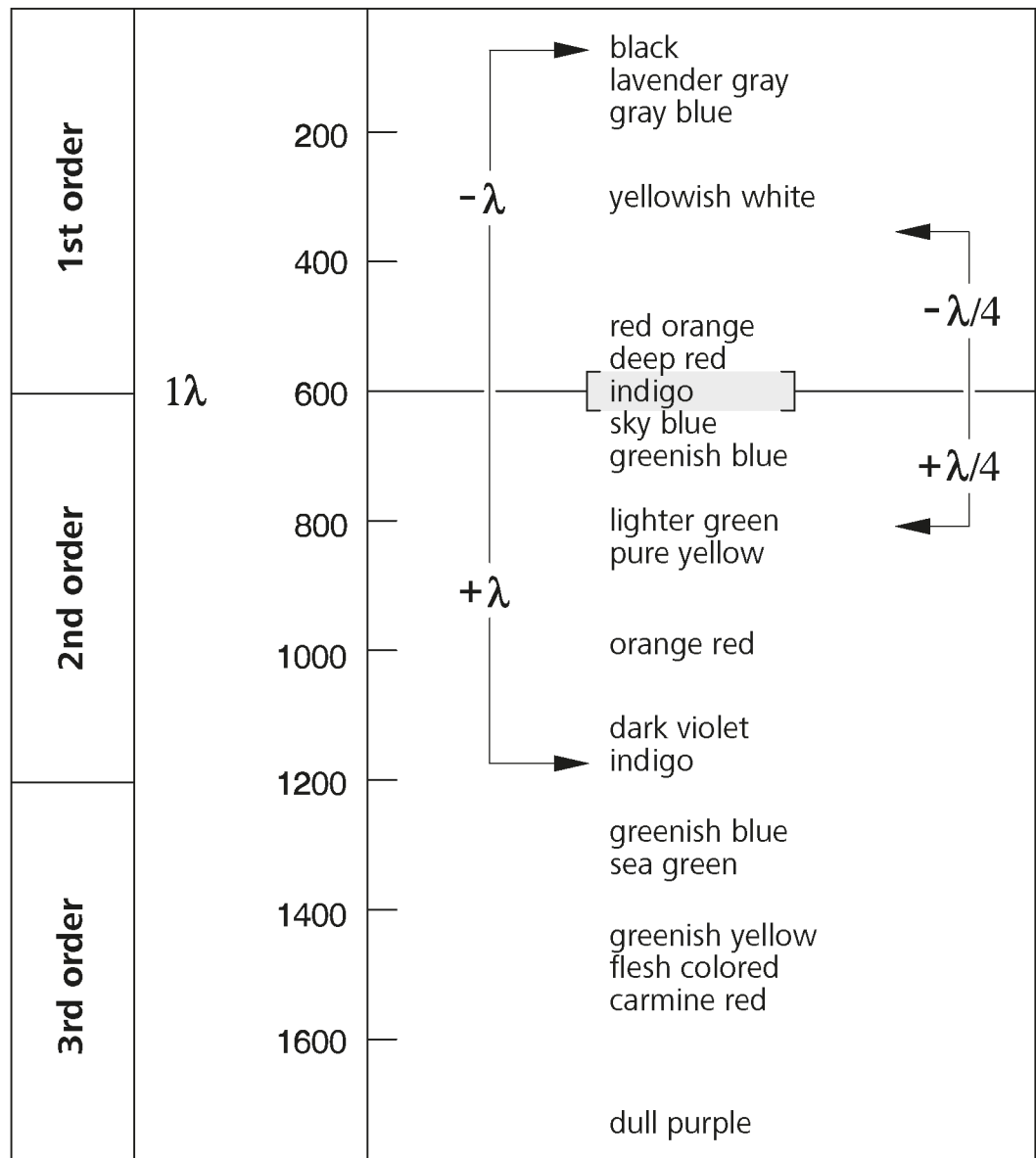


Fig. 31: Schematic diagram of the color charts developed by Michel-Lévy

When the lambda compensator is brought into the beam path, the non-birefringent "surroundings" of the synthetic fiber appear dark red, which corresponds to the path difference of the compensator of 550 nm (1st order interference color for the path difference of 550 nm corresponds to  $1\lambda$ ).

If the polarization direction ( $n_v$  or  $n_v'$ ) of the birefringent sample to be examined is parallel to the principal polarization direction ( $n_p$ ) of the lambda compensator, i.e. in the northeast-southwest direction, the path difference of the sample (e.g. gray-white: 150 nm) and the path difference of the lambda compensator (red: 550 nm) add up. This results in a color change of the sample from grayish white to greenish-blue (resulting path difference = 700 nm).

If the polarization direction of the birefringent sample to be examined is perpendicular to the principal polarization direction of the lambda compensator, i.e. in the northwest-southeast direction, the path difference of the sample (e.g. gray-white: 150 nm) is subtracted from the path difference of the compensator (red: 550 nm). In this case, the interference color of the sample visibly changes from gray-white to orange (resulting path difference = 400 nm).

### 3.4.4.3 Measuring Path Differences

Measuring compensators are required for the accurate measurement of path differences,. These compensators reset, i.e. compensate the path difference produced by the sample to zero (first order black). While the addition position as well as the subtraction position are of interest for the methods described above, only the subtraction position is relevant for measurements. Path differences in the sample can assume very small values ( $1/50 \lambda$  or 10 nm) and very large values (greater than  $10 \lambda$  or approx. 5500 nm and more) and thus determine the compensator appropriate for the measurement.

The suitable compensator is determined as follows:

- If more or less strong interference colors appear on the sample, the path difference ranges approximately between  $1/2 \lambda$  and  $5 \lambda$ .

The suitable compensator is:

B 0-5  $\lambda$  tilting compensator

- If the sample-side color changes from light gray/white to a strong interference color, when a lambda compensator (473704-0000-000) is inserted in the compensator slot, the path difference is  $(1/4 - 1/2) \lambda$ .

**NOTICE** A prerequisite for the occurrence of the color change effect may be the evaluation in two sample positions rotated at an angle of  $90^\circ$  from one another. For this purpose, rotate the centered stage (by 2 click stops).

The suitable compensator is:

B 0-5  $\lambda$  tilting compensator

or the DE SENARMONT compensation method up to  $1 \lambda$  using the 546/4 nm SENARMONT compensator.

**NOTICE** The DE SENARMONT compensation method requires the use of the rotatable analyzer.

- After insertion of the lambda compensator and rotation of the sample by  $90^\circ$ , the interference color remains white; in this case, however, it is a "higher-order white" and thus the path difference is  $> 5 \lambda$ .  
The suitable compensator is:  
K 0-30  $\lambda$  tilting compensator
- A dark gray appearing as the interference color indicates a very small path difference ( $\lambda/10$  or 54.6 nm).

### 3.4.4.4 Circular Polarization Contrast

Unlike standard polarization contrast, circular polarization contrast does not show any dark (extinction) positions that depend on the angle of rotation (azimuth) of the sample relative to polarizer or analyzer. This means that, while you are rotating the stage, the image will always look the same, as there are no bright and dark positions. With optical anisotropy all transparent samples show the interference colors that are characteristic to them.



### 3.4.4.5 Determining the Optical Character of Crystals

For the classification (and thus identification) of crystalline material – instead of the observation of the sample itself – the analysis of an interference image in the objective pupil provides the more valuable information. This image is visible in the eyepiece when an additional lens (so-called Bertrand lens) is switched on. Alternatively, the auxiliary microscope or a diopter may be used to view the interference image.

In contrast to orthoscopy, this is referred to as conoscopy, because the illumination is ideally provided by a wide open cone. In practice this means that the aperture diaphragm is fully open and the objective should likewise have a high aperture.

Crystal analysis is used to determine the optical character of transparent and weakly absorbent crystals. This method is also referred to as conoscopy.

Its main application is classic mineral microscopy. However, synthetic crystals, industrial minerals and plastics (e.g. films) can also be identified and characterized.

In the case of uniaxial crystals, the most favorable orientation for conoscopic viewing is obtained with those sample features (e.g. of a thin section) that in orthoscopic viewing change the brightness as little as possible. In this case, the direction of viewing and the optical axis are parallel. The same applies to biaxial crystals if viewed in or approximately in the direction of one of the two optical axes.

Crystalline anisotropic samples can be separated into optical uni- and biaxial samples, in each case with an "optically positive" or "negative" character.

**Uniaxial** crystals display a **black cross** when the optical axis is parallel to the direction of view. Depending on the level of birefringence and sample thickness, concentrically arranged colored interference rings (so-called isochromes) may appear (see also table below, second row).

The lines of this black cross remain closed when the stage is rotated. Depending on the section it may lie within or outside the displayed objective pupil.

With optically biaxial crystals, the cross resolves into two dark hyperbola branches (the so-called isogyres) depending on stage rotation, which are surrounded by colored interference patterns depending on the amount of birefringence and sample thickness (suggestive of the figure "8").

Inserting a lambda compensator or lambda/4 or a 0-4 lambda wedge compensator in the compensator slot with the initial state of the axial figure being as illustrated in the table below results in the following changes in color (shown schematically as blue and yellow areas) to the axial figure, thus allowing differentiation in "optically positive" and "optically negative".

	Optically uniaxial		Optically biaxial		
	Positive	Negative	Positive	Negative	
Lambda plate (white → blue → yellow)					+ = Blue – = Yellow
Quartz wedge (direction of motion at insertion)					Arrow = Direction of movement
Lambda/4 plate (position of black spots)					

Tab. 1: Determining optical character

In the case of less favourable sections in which the cross-hair center is optically uniaxial or the isogyres are optically biaxial samples outside the objective pupil, an assessment is possible as follows:

- If the black isogyres are **straight** and they run parallel to the pupil (in relation to the cross-hairs), the sample is **optically uniaxial**.
- If the black isogyres are **curved lines** which wander on a circular path through the pupil, the sample is **optically biaxial**.

#### Info

Axial figures can often be better displayed with circular polarization. In particular, the axis angle of optically biaxial samples (quasi the distance between the isogyres) can be determined more clearly. The optical character can also be determined. The lambda compensator (6x20mm), arranged in the compensator slot, is used for this purpose.

#### Info

Two storage compartments for compensators (6x20) are located on the back side of the conoscopy stand.

### 3.4.5 Transmitted Light Polarization for Conoscopic Observation

The determination of the optical character of transparent and weakly absorbent crystals is used to diagnose crystals. This method is also termed conoscopy. Its main application is classical mineral microscopy. However, it also facilitates the identification and characterization of synthetic crystals, industrial minerals and plastics (e.g. films).

For the classification (and thus identification) of crystalline matter, the examination of the interference image in the objective pupil delivers more valuable information than that obtained by viewing the sample itself. The interference image becomes visible in the eyepiece if an additional optical system (fixed or focusing Bertrand lens or, in the basic version, the auxiliary microscope or diopter) is used.

In contrast to orthoscopy, this technique is called conoscopy, because here the sample is ideally illuminated through a wide-open cone. In practical microscopic work, this means that the condenser front lens (0.9) must be in the light path, the aperture diaphragm fully open, and the objective, too, should be a high-aperture type.

### 3.4.6 Reflected Light Brightfield Microscopy Using the KÖHLER Method

Reflected light brightfield microscopy is the easiest and most commonly used RL-microscopy method. It is used to examine optically opaque samples or samples as e.g. cut, polished, etched metal or ores.

In order to obtain an image as close as possible to the object, not only the so-called direct beam bundles but also the indirect ones, i.e. the beam bundles diffracted and scattered at the preparation details, are of essential importance. According to ABBE, the larger the indirect beam components are, the more true to the object the microscopic image is.

The cone of light emerging from the reflected light light source is reflected on a color-neutral beam splitter before it passes through the objective which is focused on the sample surface (so-called condenser function). The objective collects the light reflected on the sample and creates, with the tube lens, the microscopic intermediate image. This image can then be examined visually or documented using a camera.

### 3.4.7 Reflected Light Darkfield Microscopy Using the KÖHLER Method

The reflected light darkfield method is applied when samples are examined, which do not have areas with different reflectivity (ideal brightfield samples), but which show deflections (as scratches, cracks, dust particles etc.) on the plane surface. All such light-scattering details appear bright in the darkfield, while the reflective plane areas remain dark.

### 3.4.8 Reflected Light Polarization Microscopy

Reflected light polarization is a contrasting method suited for cut, polished surfaces of mineral ore, coal, ceramics, special metals and alloys. Depending on the orientation of the crystals and the sample details, the cut surfaces often react differently when reflected in linearly polarized light.

The illumination light is polarized by the polarizer before passing through the objective onto the sample surface where it is reflected. Then the beam parts experience path differences depending on the structure and polarization of optical rotations which, when passing through the analyzer, are represented by different shades of gray. With the aid of a compensator with a  $\lambda$ -plate the gray contrast can be converted into a color contrast.

Even when examining "dark" sample surfaces, a rotatable  $\lambda/4$  plate in front of the objective (antireflective cap) helps eliminate the reflections which are inevitable when working with objectives with very low magnification.

A sample is birefractant when the sample details show differences in brightness and color which change when the direction of vibration of the polarizer or the stage is rotated. For samples with low birefractance using the analyzer with a rotatable lambda plate is recommended.

### 3.4.9 Reflected Light Fluorescence Microscopy

The reflected light fluorescence method is used to show fluorescent substances in typical fluorescent colors in high contrast. The light originating from a high-performance light source in a reflected light fluorescence microscope passes through a heat protection filter onto an excitation filter (bandpass). The filtered short-wave excitation radiation is reflected by a Dichroic Beam splitter and is focused on the sample through the objective. The sample absorbs the short-wave radiation before emitting longer-wave fluorescence radiation (Stokes' Law). This radiation is then captured from the image side by the objective and passes through the Dichroic Beam splitter. Last, the beams pass through a emission filter (longpass/bandpass) and only the long-wave radiation emitted by the sample passes.

The spectra of the excitation and the emission filter must match very closely. They must be inserted in a Reflector Module FL EC P&C reflector module together with the according Dichroic Beam splitter.

## 4 Installation


Perform only the installation work described in this document. All other installation work not described may only be carried out by an authorized ZEISS service representative.

### 4.1 Unpacking and Setting up the Microscope

- Procedure**
1. Open the packaging.
  2. Take the microscope, all assemblies, and accessories out of the packaging.
  3. Check them for completeness as per delivery note.
  4. Check all parts for damaging.
  5. Place the microscope on a vibration-free, level, and non-inflammable surface.

It is recommended to keep the original packing and store it away for later use, e.g. for stowing the microscope during periods of non-use or for returning the microscope to the manufacturer for repair.

## 4.2 Assembling the Binocular Tube

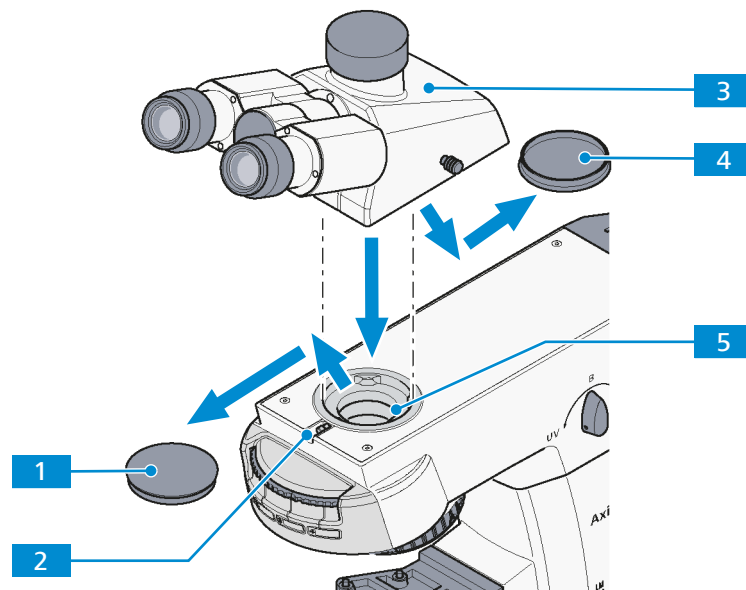


**Parts and Tools**  Hex key, 3.0 mm

### Procedure

1. Loosen the clamping screw **2**.
2. Remove the dust cap **1** from the dovetail ring mount on the stand side.
3. Remove the dust cap **4** from the underside of the binocular tube **3**.
4. Hold the binocular tube at an angle, insert it with the dovetail ring into the stand mount **5** and turn into a horizontal position.
5. Rotate the binocular tube into the desired observation position.
6. Re-tighten the clamping screw with the hex key.

### 4.3 Assembling the Binocular Tube with an Intermediate Plate

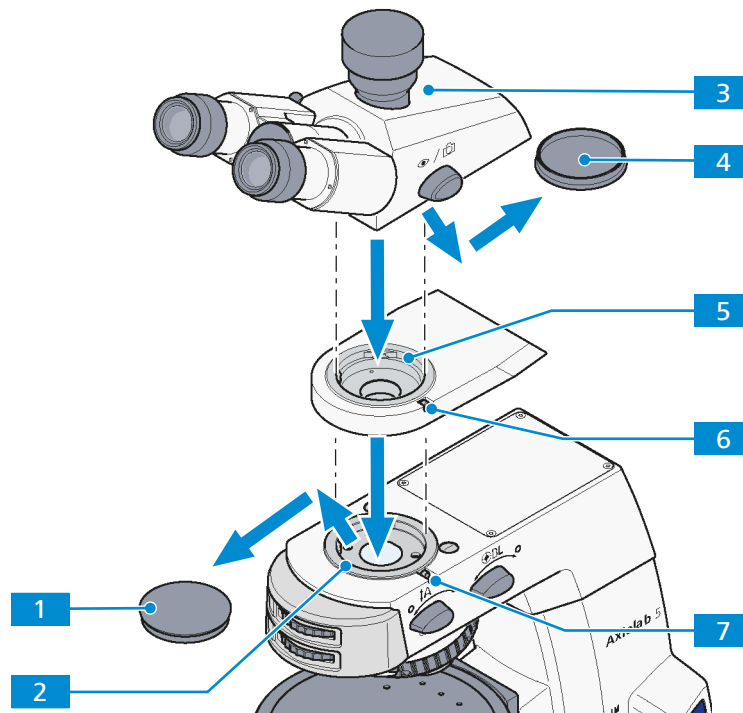

**Parts and Tools**  Hex key, 3.0 mm


#### Procedure

1. Loosen the clamping screw of the stand **7**.
2. Remove the dust cap **1** from the dovetail ring mount on the stand side.
3. Insert the intermediate plate **5** with its dovetail ring into the stand mount **2**.
4. Align it to the stand and tighten the clamping screw of the stand **7**.
5. Remove the dust cap **4** from the underside of the binocular tube **3**.
6. Hold the binocular tube **3** at an angle, insert it with the dovetail ring into the intermediate plate **5** and turn into a horizontal position.
7. Rotate the binocular tube into the desired observation position.
8. Re-tighten the clamping screw of the intermediate plate **6** with the hex key.

## 4.4 Assembling Components into the Binocular Tube

The following components can be inserted into the tube:

- eyepieces
- auxiliary microscope
- pinhole diaphragm

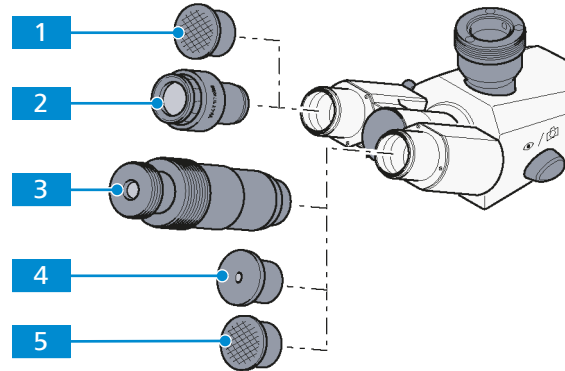


### Procedure

1. Remove both dust caps **1** / **5** from the tube.
2. Remove both eyepieces **2** from the box and insert them into the eyepiece socket of the tube to the stop.  
**NOTICE** Before inserting Pol eyepieces with tubes without upright reticles, the orientation screw on the reverse side of the eyepieces must be unscrewed. The eyepieces cannot otherwise be fully inserted.
3. Instead of an eyepiece insert an auxiliary microscope **3** or pinhole diaphragm **4** in one eyepiece socket.


Proceed in the reverse order for removal.

## 4.5 Assembling the Objectives

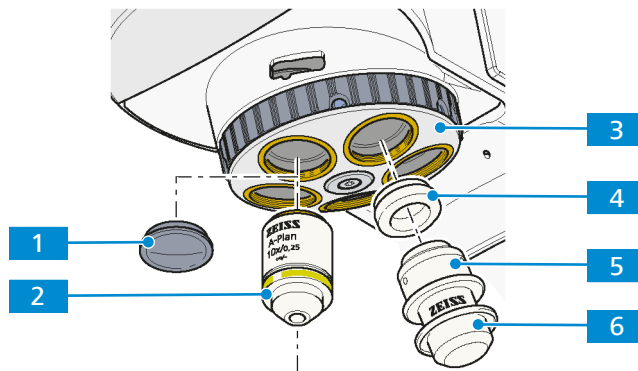

**Prerequisite** ✓ Adapter W0.8/M27 is required when using W0.8 objectives or the sample marker.

**Procedure** 1. Move mechanical stage to lower stop using the focusing drive.

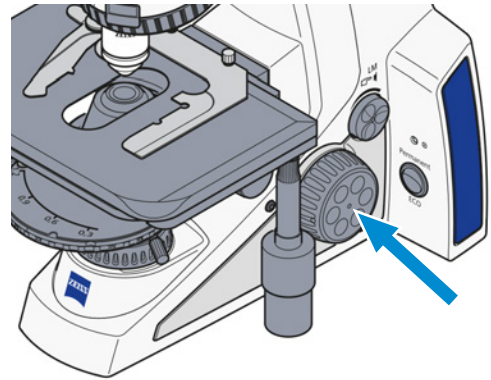


2. Remove the dust protection caps **1** from the appropriate openings in the nosepiece.
3. Remove objectives **2** from the case and screw them into the nosepiece **3**.
4. Carefully screw the objective into the opening. Start with the smallest magnification factor (set up clockwise) in nosepiece position 1.
5. Make sure it engages properly in the nosepiece's thread.
6. Instead of an objective, the sample marker **5** with an adapter W0.8/M27 **4** can be screwed on in any desired nosepiece position.
7. Apply the protective cap **6** to prevent sample marker from drying out.
8. Always replace the dust protection caps on any empty positions on the nosepiece.



## 4.6 Assembling Condensers


### 4.6.1 Assembling the Condenser 0.9/1.25 BF, DF, Ph1, Ph2, Ph3 with Modulator Disk

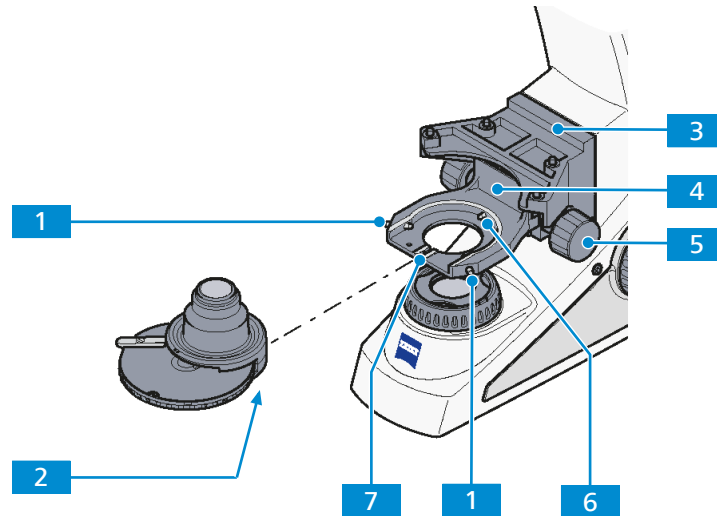

#### Procedure


1. Carefully move the stage carrier **3** with the focusing drive to the upper stop position.

**NOTICE** Make sure that the stage does not collide with the objective.

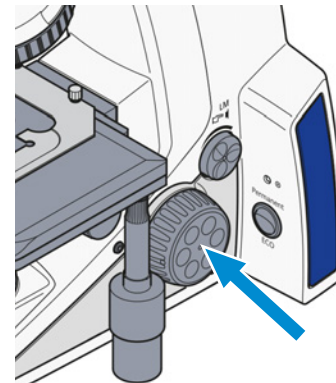


2. Unscrew both centering screws **1** on the condenser carrier until their ends are no longer visible.
3. Using the knurled knob for vertical adjustment **5**, push the condenser carrier down as far as it will go.  
**NOTICE** If using a low-power system, make sure that this does not come to rest on the luminous-field diaphragm.
4. Insert the condenser between the condenser carrier **4** and the stage carrier **3**. In doing so, align the screwed stud bolt **2** on the underside of the condenser with the groove **7** of the condenser carrier.
5. Press the condenser with the dovetail ring against the mainspring **6** of the condenser carrier until the condenser sits horizontally on the condenser carrier.
6. Screw in the centering screws until they engage with the dovetail ring of the condenser.

### 4.6.2 Assembling the Dry Darkfield Condenser

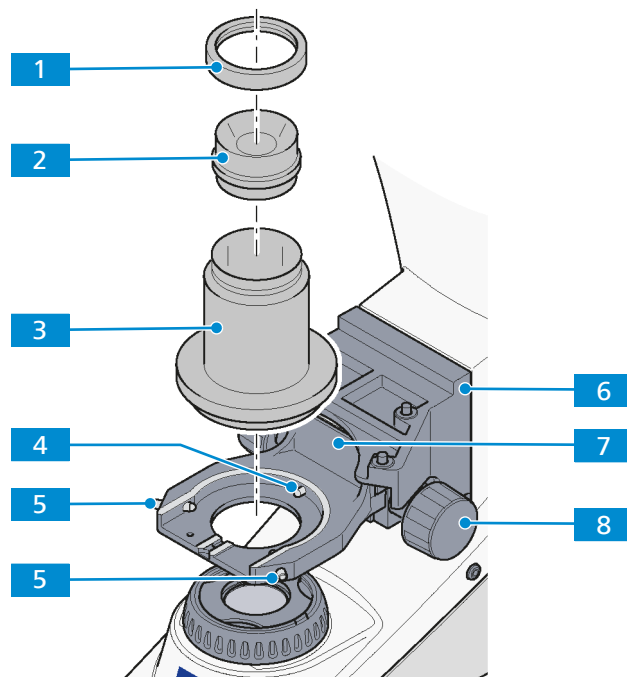

- Procedure**
1. Carefully move the stage carrier **6** with the focusing drive to the upper stop position.

**NOTICE** Make sure that the stage does not collide with the objective.

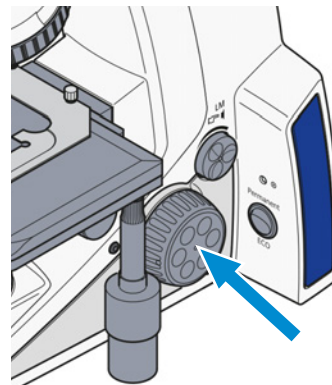

2. Using the knurled knob **8** for vertical adjustment, push the condenser carrier down as far as it will go.  
**NOTICE** If using a low-power system, make sure that this does not come to rest on the luminous-field diaphragm.
3. Insert the darkfield condenser **2** in the condenser holder Z **3**.
4. Fix the darkfield condenser with the fastening ring **1**.
5. Unscrew both centering screws **5** on the condenser carrier **7** until their ends are no longer visible.
6. Press the condenser holder Z with the dovetail ring against the mainspring **4** of the condenser carrier until the condenser holder Z sits horizontally on the condenser carrier.
7. Screw in the centering screws until they engage with the dovetail ring of the condenser holder Z.

### 4.6.3 Assembling the Condenser, Chromatic-aplanatic 0.9 BF DF PhC DIC

#### Info

If an additional component, e.g. a polarizer, has been mounted beneath the condenser carrier, the stage carrier should be removed before installing the condenser.


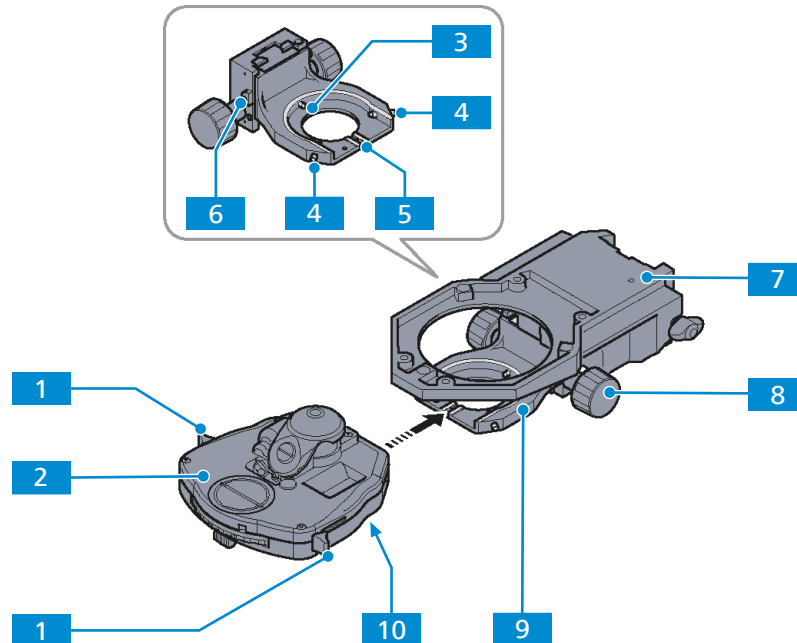


Fig. 32: Installing the condenser, achromatic-aplanatic 0.9 BF DF PhC DIC

#### Procedure

1. Carefully move the stage carrier **7** to the upper stop position. Use the focusing drive.

**NOTICE** Make sure that the stage does not collide with the objective.

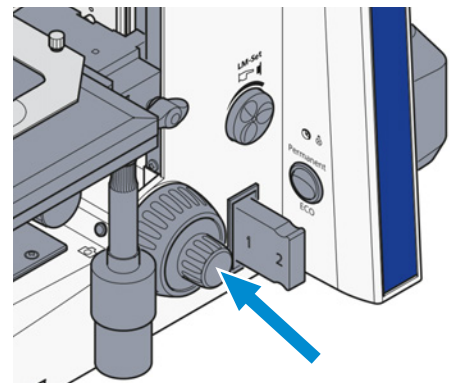


2. Swivel out the front lens on the condenser **2** using the lever **1**.
3. Unscrew both centering screws **4** on the condenser carrier **9** until their ends are no longer visible.
4. Loosen the clamping screw **6** of the condenser carrier until the maximum vertical adjusting range is usable.
5. Using the knurled knob **8** for vertical adjustment, push the condenser carrier down as far as it will go.  
**NOTICE** If using a low-power system, make sure that this does not come to rest on the luminous-field diaphragm.
6. Insert the condenser between the condenser carrier and the stage carrier **7**. In doing so, align the screwed stud bolt **10** on the underside of the condenser with the groove **5** of the condenser carrier.
7. Press the condenser with the dovetail ring against the mainspring **3** of the condenser carrier until the condenser sits horizontally on the condenser carrier.


8. Screw in the centering screws **4** until they engage with the dovetail ring of the condenser.
9. Screw in the clamping screw **6** without clamping the vertical drive.

Proceed in the reverse order for removal.

#### 4.6.4 Assembling the Modulator Disk in the Condenser 0.9 BF Pol

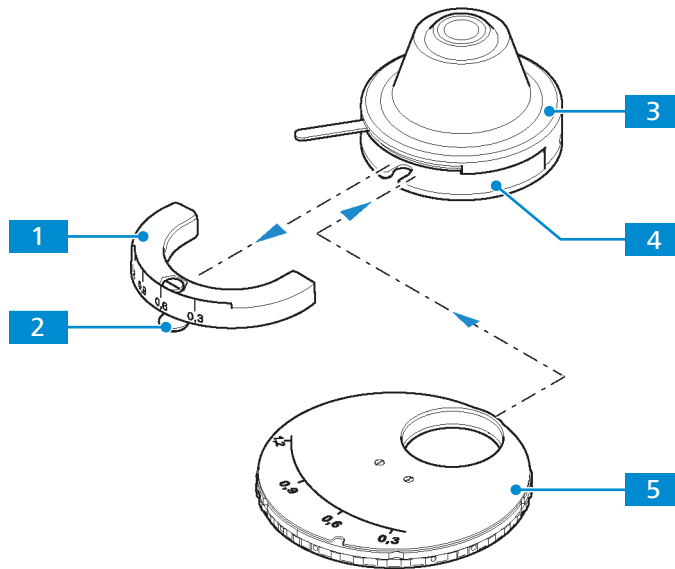

**Parts and Tools**  Hex key, 3.0 mm


**Prerequisite**  The condenser **3** is removed from the condenser carrier [► 59].

- Procedure**
1. Loosen the clamping screw **2** of the condenser's dial segment **1** and pull out the dial segment.
  2. Slide the modulator disk **5** with its two-pronged forked opening pointing forward into the condenser opening **4**.
  3. Make sure that the disk engages in the guide on both inner sides of the condenser.
    - The guide serves as a stop for the modulator disk. The pin of the disk's clamping screw must slide into the orientation groove of the condenser.
  4. Tighten the disk's clamping screw.
  5. Replace the condenser in its carrier.

## 4.7 Loading the Reflector Turret


The reflector turret with four positions is firmly installed in the reflected light illumination module for fluorescence or materials stands.

### 4.7.1 Assembling Reflector Modules

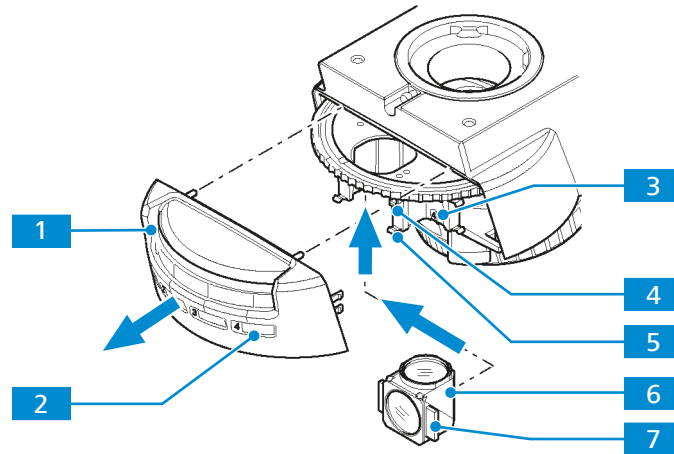

- Procedure**
1. Remove the cover cap **1** from the stand towards the front.
  2. Turn the reflector turret to the desired position **3**.
  3. Insert the reflector module **6** together with the retaining brackets **7** on the right and left diagonally from below into the upper spring clips **4** (right/left) of the reflector turret position.
  4. Apply pressure to the reflector module from below until it also securely engages with the lower spring clips **5** of the reflector turret position.
  5. Fit the cover cap onto the stand as straight as possible to avoid the knurled ring of the reflector turret becoming jammed and damaged.
  6. Apply pressure to the cover cap until the retaining brackets have engaged.
  7. Apply the supplied stickers with the filter combination data of the respective module to the corresponding field **2** of the cover cap (positions 1 to 4).

### 4.7.2 Removing Reflector Modules

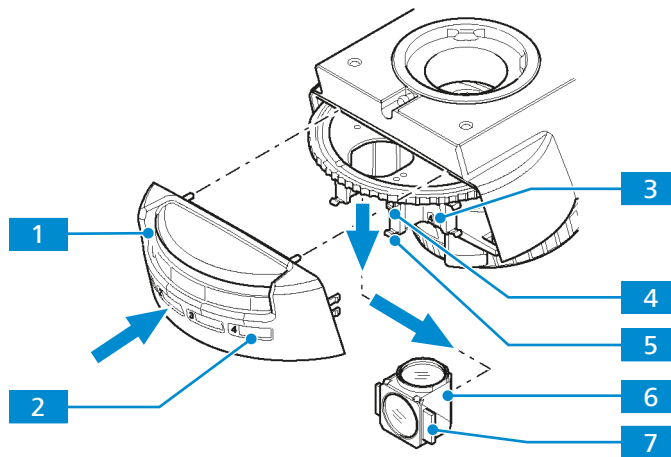


- Procedure**
1. Remove the cover cap **1** from the stand towards the front.
  2. Turn the reflector turret to the desired position **3**.
  3. Slightly tilt the reflector module **6** in order to detach it from the lower spring clips **5**, then from the upper spring clips **4** of the reflector turret using the retaining bracket **7**.
  4. Fit the cover cap onto the stand as straight as possible to avoid the knurled ring of the reflector turret becoming jammed and damaged.
  5. Apply pressure to the cover cap until the retaining brackets have engaged.
  6. Update the filter data on the sticker for the corresponding field **2** on the cover cap.

## 4.8 Connecting the Microscope to the Mains

**Prerequisite** ✓ The microscope is switched off.

- Procedure**
1. Connect the power cord to the mains socket of the stand.

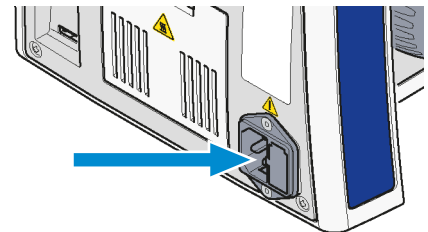

2. Connect the power cord to the mains.

## 5 Operation

This chapter describes switching on/off the microscope as well as the operating steps with the microscope.

### Info

For additional information and detailed descriptions, refer to further applicable documents or ask your ZEISS Sales & Service Partner.

### Info

Further information on the software and its operation is available in the software's online help.

### 5.1 Prerequisites for Commissioning and Operation

The following basic prerequisites are necessary for commissioning and operation:

- This document was read prior to commissioning or operation and kept for further use.
- The chapter **Safety** was read and understood.
- The operator is acquainted with the general Windows-based programs.
- If required: Basic training and safety briefing were successfully completed.

### 5.2 Switching On the Microscope

#### **WARNING**


##### **Skin or eye injury due to hazardous light emission**

The light source belongs to Risk Group 3 as specified in IEC 62471 and emits LED radiation and UV radiation. Skin or eye injury can result from the exposure.

- ▶ Avoid any eye and skin exposure to the light-emitting aperture of the light source.
- ▶ Avoid exposure of skin to the radiation. Use suitable protective equipment/protective clothing if required.
- ▶ Before installing or removing the light source always make sure it is switched off.

**Prerequisite** ✓ The microscope is connected to the mains.

**Procedure** 1. Switch the microscope on using the power switch **On/Off**.

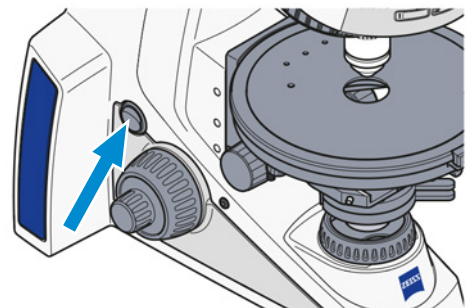


## 5.3 Adjusting

### 5.3.1 Adjusting the Position of the Eyepieces


#### Info

The adjustment of the interpupillary distance is correct when you see only one round image while looking through the two eyepieces.

- Procedure**
1. Set the interpupillary distance by rotating the eyepiece tubes symmetrically toward or away from one another.

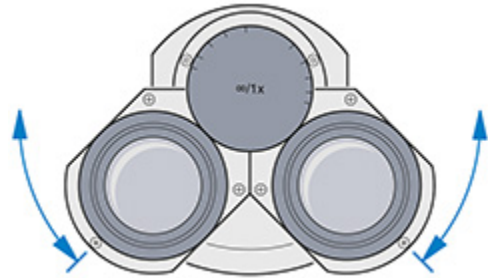

2. Set the viewing height by swivelling the whole eyepiece unit a full 180 ° upwards or downwards.

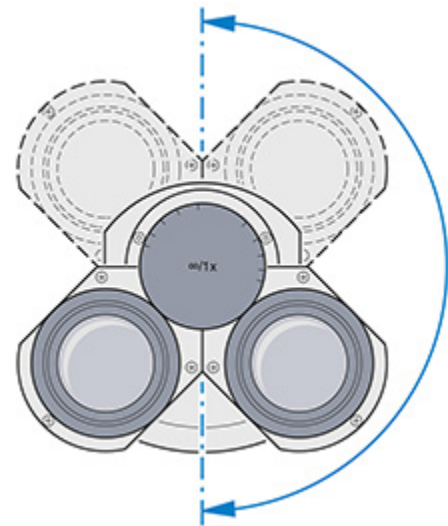

### 5.3.2 Adjusting for Ametropia when Using Eyepiece Reticles



- Prerequisite**
- ✓ Two adjustable eyepieces are installed
  - ✓ One eyepiece with reticle is installed.

- Procedure**
1. Focus on the line figure of the eyepiece reticle with the focusable eye lens of the adjustable eyepiece containing the eyepiece reticle.
  2. Focus on the microscopic image of a loaded sample with the focusing drive while observing with the eyepiece containing the eyepiece reticle.  
→ Both the microscopic image and the eyepiece reticle are in focus now.
  3. Focus the microscopic image for the second eye with the focusable eye lens of the second eyepiece.  
↳ Both microscopic images including the eyepiece reticle are thus in focus.  
From this point, use only the focusing drive for any subsequent focusing activity.




### 5.3.3 Adjusting the Height Stop on the Condenser Carrier


**Parts and Tools**  Hex key, 3.0 mm

**Prerequisite**  The microscope is operational.  
 A sample is positioned on the stage.


**Procedure** 1. Loosen the fastening screw of the height stop.

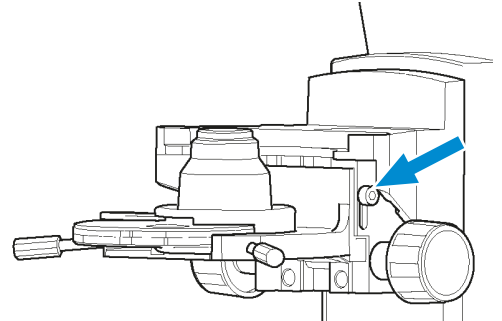


2. Focus the sample with the focusing drive.

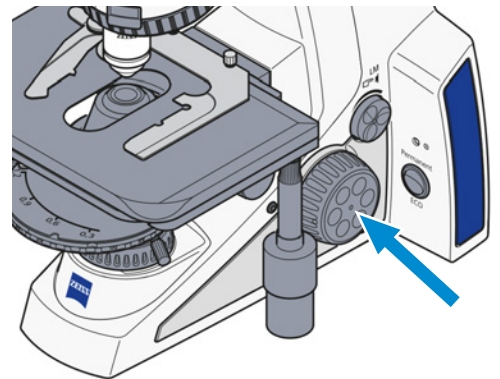

3. Close the luminous-field diaphragm.

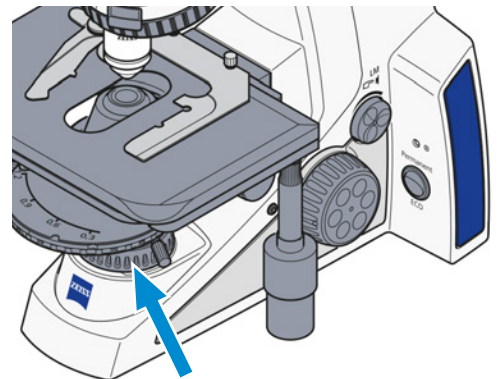


4. Adjust the condenser vertically until a sharp image is visible.

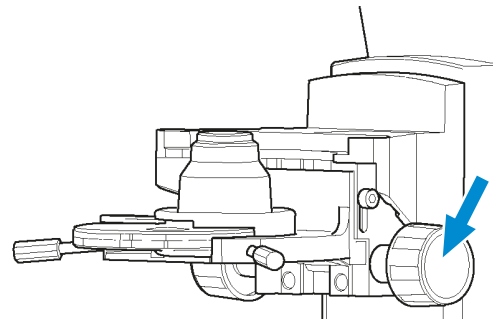


5. **NOTICE** The sample and the objective can be damaged when the sample is lifted out.

Carefully raise the condenser by a small amount without lifting out the sample.

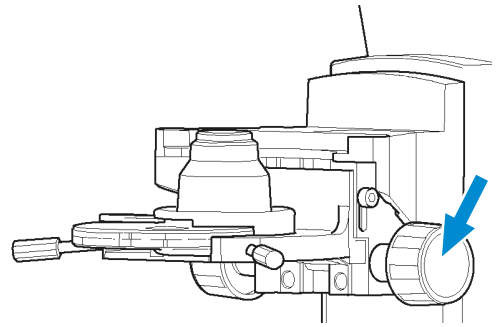

6. Tighten the fastening screw of the height stop.

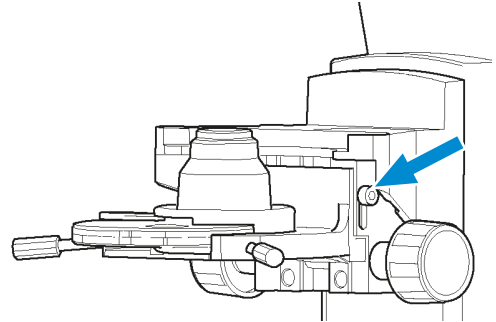


↳ The height stop is adjusted.

### 5.3.4 Using the Light Manager Function

#### 5.3.4.1 Switching On and Off the Light Manager Function

**Prerequisite** ✓ The microscope is operational.

- Procedure** 1. Press one of the **Snap** buttons and the **Intensity/LM** knob simultaneously for at least 1.5 seconds.

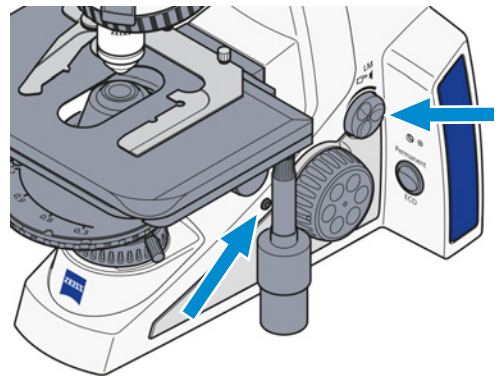


- The indicator light blinks in the following sequence, when the Light Manager is switched on: GREEN / GREEN / GREEN
- The indicator light blinks in the following sequence, when the Light Manager is switched off: GREEN / ORANGE / GREEN


### 5.3.4.2 Saving Light Intensity Values

- Prerequisite**
- ✓ The microscope is operational.
  - ✓ The Light Manager function is *enabled* [▶ 66].

- Procedure**
1. Switch to the first objective and/or reflector (if available) positions of interest using the knurled rings.

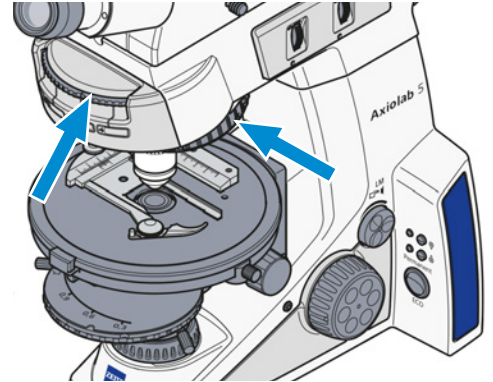

2. Set the desired light intensity by turning the **Intensity/LM** knob.

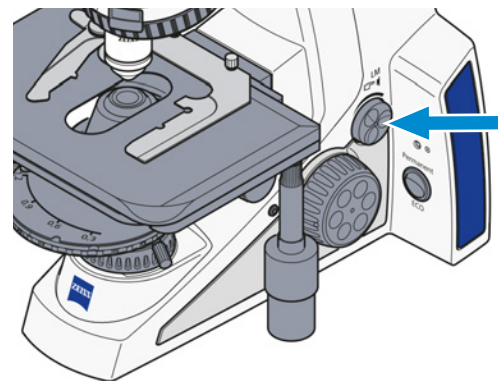

3. Press the **Intensity/LM** knob for at least 1.5 seconds.
  - The light intensity for this objective/reflector combination is saved.
  - When using LED as light source, the LED is switched off for 300 ms. This is visible through the eyepieces and serves as an indicator for the user.
4. Switch to the second objective/reflector position.
5. Press the **Intensity/LM** knob for at least 1.5 seconds.
  - Now a ratio between the first and the second objective/reflector combination is established.
6. Repeat the procedure to set the light intensity values for further objective/reflector combinations.
  - ↳ The light intensities for all set objective/reflector combinations are stored.
  - ↳ After switching on the microscope, the previous setting of the Light Manager will be restored.

### 5.3.4.3 Enabling/Disabling the Dazzle Protection Function


By factory default, the dazzle protection function is enabled.

- Procedure**
1. Press the left **Snap** button for at least 1.5 seconds to switch between enabled and disabled dazzle protection function.
    - Dazzle protection function disabled: The indicator light blinks ORANGE twice.
    - Dazzle protection function enabled: The indicator light blinks GREEN twice.

### 5.3.5 Setting the ECO/Permanent Mode

**Prerequisite** ✓ The microscope is operational.

**Procedure** 1. Select the ECO or Permanent mode for microscope illumination using the **ECO/Permanent** mode switch.

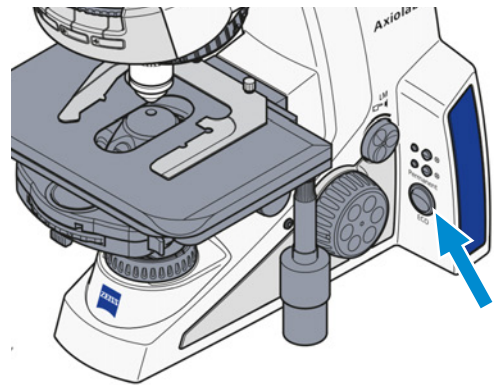

## 5.4 Setting Up for Transmitted Light Techniques


### 5.4.1 Setting Up for Transmitted Light Brightfield Microscopy

- High-contrast sample slide


The equipment of every microscope stand model enables the use of transmitted light brightfield microscopy. All available condensers (except special condensers such as darkfield condensers) can be used for transmitted light brightfield microscopy.

**Prerequisite** ✓ The microscope is operational.  
 ✓ The microscope is *adapted* [▶ 64] to the user.  
 ✓ The height stop of the condenser carrier is *adjusted* [▶ 65].

**Procedure** 1. If required, push the **TL** button for transmitted light illumination.

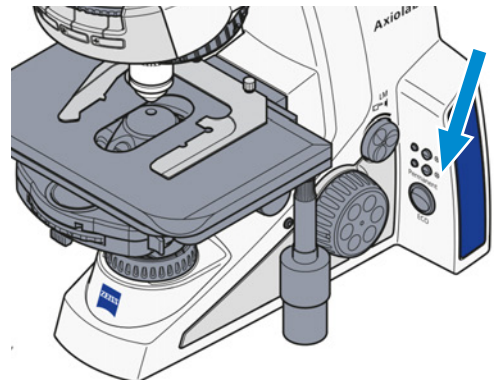


2. Adjust the image brightness using the **Intensity/LM** knob on the microscope stand.

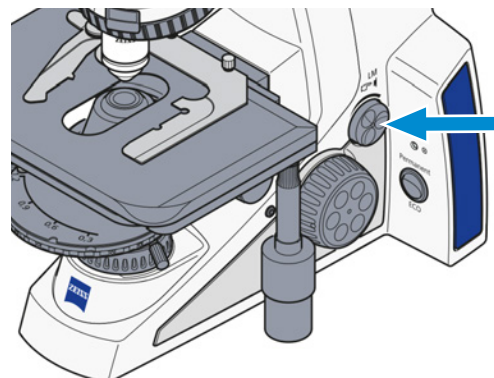


3. Insert the high-contrast sample into the sample holder of the stage.


4. Set position H (or B = brightfield), when using condensers with a turret/modulator disk and knurled ring.

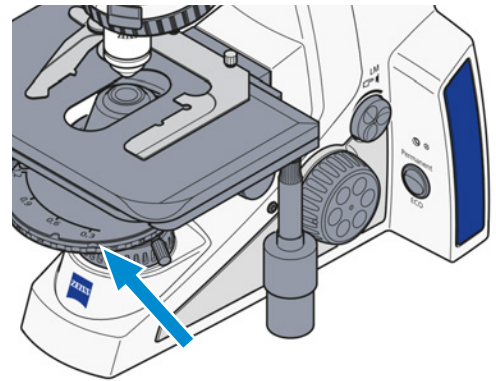


5. Swivel the front lens into the beam path with  $\geq 10\times$  objectives, if condensers with a swiveling front lens are used.

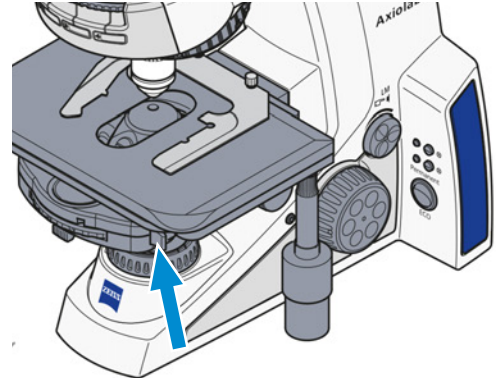

6. Set the condenser with the knurled knob for vertical adjustment to the upper stop.

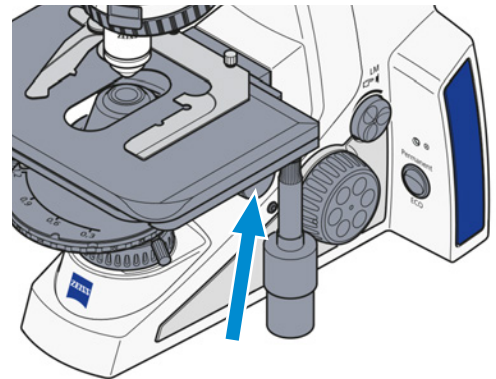


7. Swivel in the 10x objective on the nosepiece.

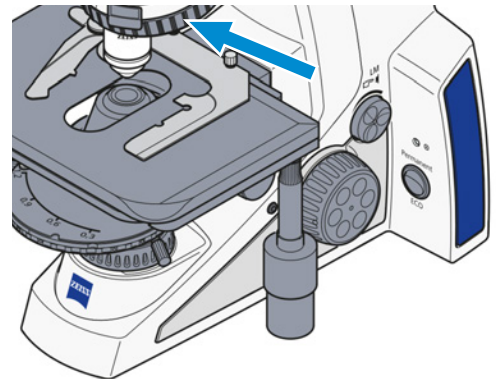




8. Focus the sample.

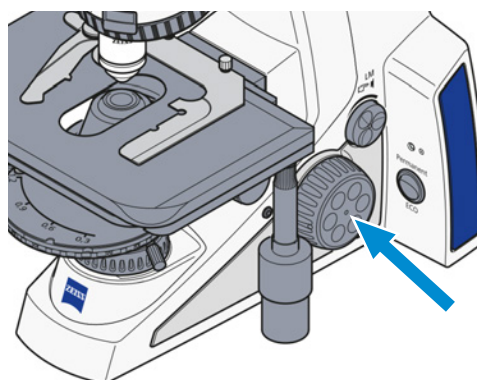


9. Close the luminous-field diaphragm until it is visible (even if not in focus) in the field of view

**A**

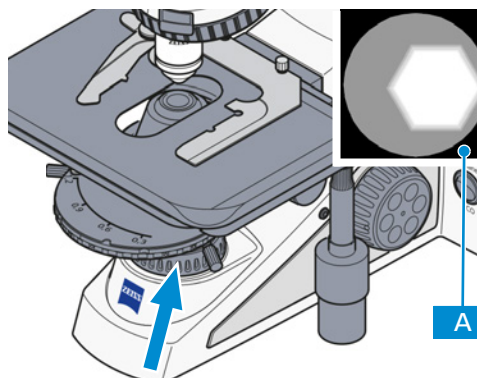


10. Turn the knurled knob for vertical adjustment to lower the condenser until the edge of the luminous-field diaphragm appears in focus

**B**

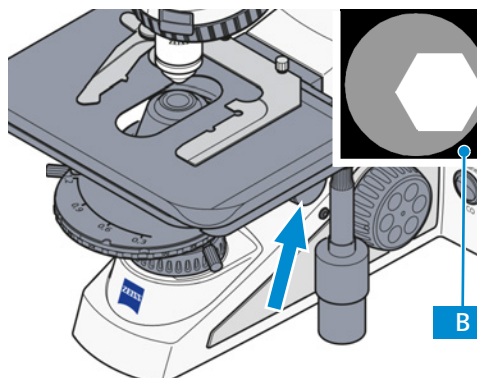


11. Center the luminous-field diaphragm using the two centering screws on the condenser carrier


**C**

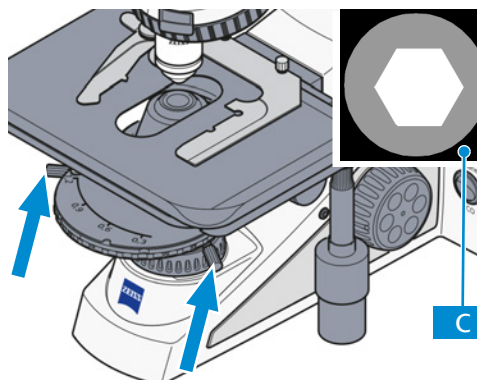

12. Open the luminous-field diaphragm until the edge of the diaphragm just disappears from the field of vision **D**.

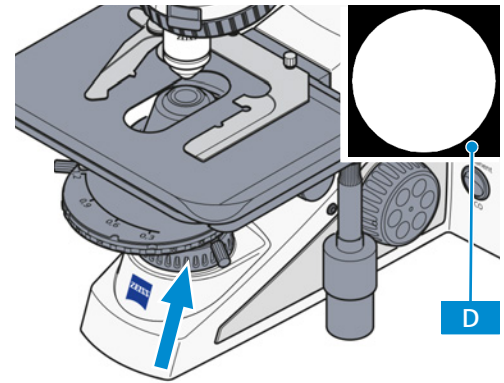

13. Remove an eyepiece from the binocular tube to adjust the aperture diaphragm (contrast).  
 14. Look into the tube with the naked eye.  
 15. Set the aperture diaphragm with the adjusting lever to between  $\frac{2}{3}$  -  $\frac{4}{5}$  of the diameter of the exit pupil of the objective **E**.

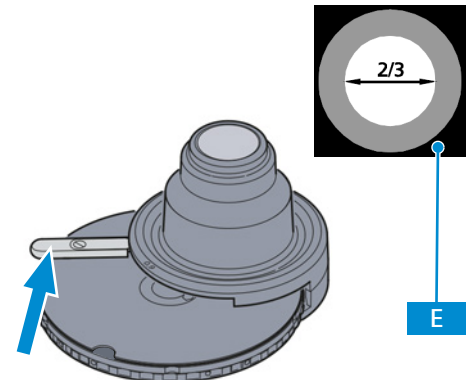

→ In most applications, this aperture diaphragm setting provides optimal contrast at almost ideal resolution, and is therefore the best compromise for the human eye.

16. Reinsert the eyepiece into the binocular tube.  
 17. Remove the high-contrast sample.  
 ↳ The illumination is now adjusted according to the KÖHLER method.

### Info


Every change of objective will result in a change in sample field size and objective aperture, together with a possible slight change in centering, so that for optimal results the luminous-field and aperture diaphragm adjustments must be repeated.

With objectives  $< 10\times$ , the front lens of the condenser (if swivelable) must be swivelled out of the beam path and the aperture diaphragm completely opened. For better contrast with such large object fields, the luminous-field aperture should be closed to a certain extent. Overclosing should be avoided so as not to impair the uniformity of the illumination of the field of view.


## 5.4.2 Setting Up for Transmitted Light Darkfield Microscopy

- Prerequisite**
- ✓ The microscope is operational.
  - ✓ The height stop of the condenser carrier is *adjusted* [▶ 65].
  - ✓ A suitable condenser for transmitted light darkfield microscopy is *installed* [▶ 57].
  - ✓ The illumination is adjusted for *transmitted light brightfield microscopy* [▶ 68].


**Procedure** 1. Set the modulator disk to position D (or DF = darkfield).

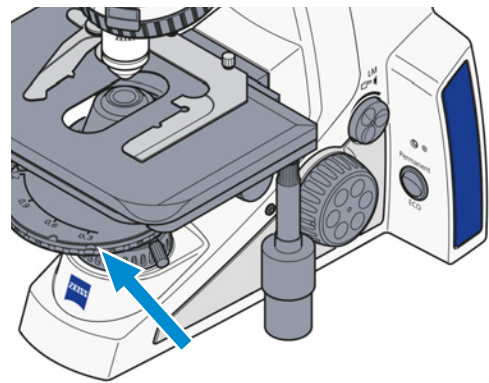

2. Turn the knurled knob for vertical adjustment of the condenser to the upper stop.

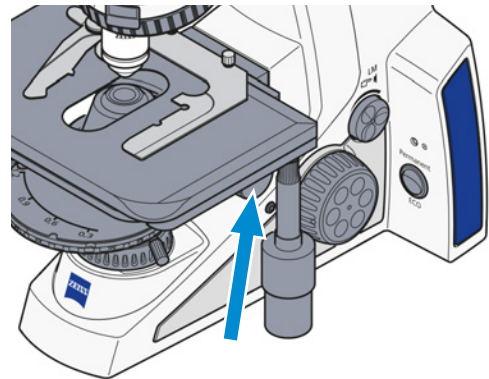


3. Swivel the objective with the highest possible aperture into position on the nosepiece.

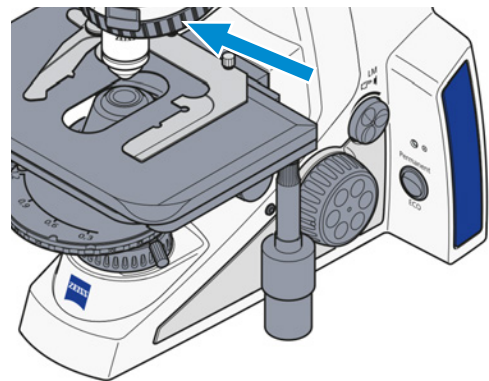


4. Place the sample on the stage.

5. Focus the sample.

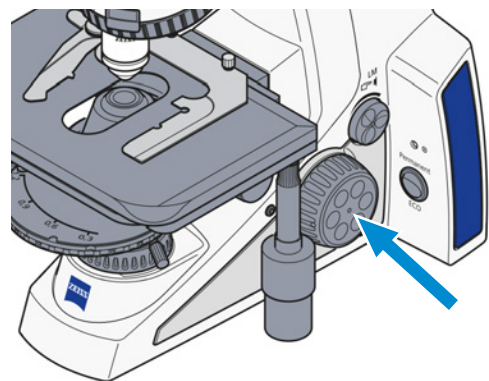





6. Close the luminous-field diaphragm enough to make it visible in the field of view (even if blurred).

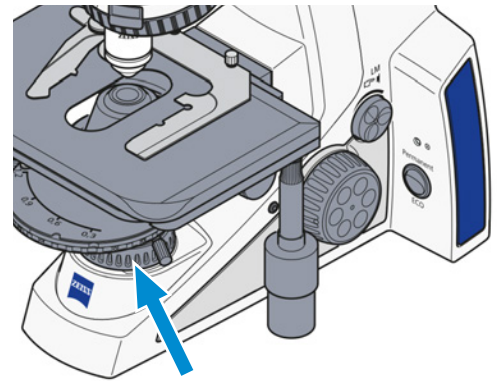

7. Lower the condenser until the edge of the field diaphragm appears sharp using the knurled knob for vertical adjustment.

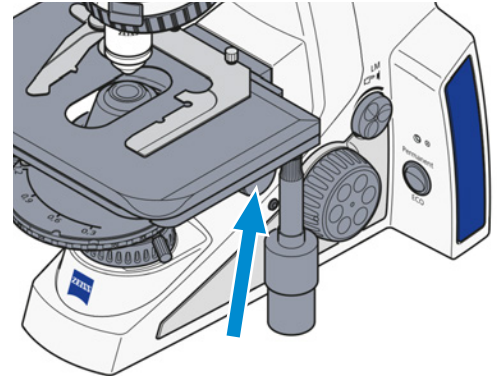

8. Center the luminous-field diaphragm on the condenser carrier using the adjustment screws.

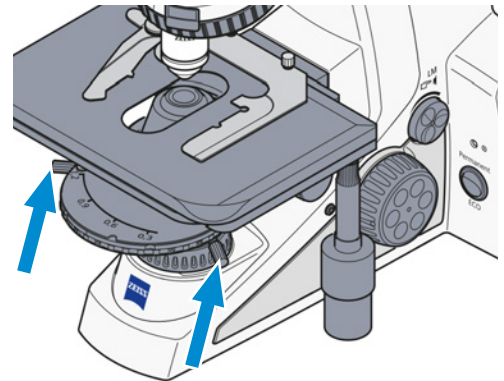

9. Open the luminous-field diaphragm enough to make the edge of the diaphragm disappear from the field of view.
  10. Remove one eyepiece or replace it with the auxiliary microscope.
  11. Check the centering of the darkfield diaphragm in the objective exit pupil.  
→ The objective exit pupil must appear homogeneously dark.
  12. If necessary, *center* [▶ 131] the darkfield diaphragm.
  13. if required, remove the auxiliary microscope.
  14. Insert the eyepiece.
  15. Adjust the condenser height using the knurled knob for vertical adjustment until no more lighter areas are visible in the field of view .
  16. Adjust the luminous-field diaphragm diameter to the size of the field of view.
- ↳ The illumination is now adjusted for transmitted light darkfield.


### Info

Darkfield microscopy requires samples to be considerably cleaner than in other techniques. In particular, fingerprints, dirt or dust particles have a negative effect, as they brighten the background of the field of view and decrease the contrast of the object image.

### 5.4.3 Setting Up for Transmitted Light Phase Contrast Microscopy

- Prerequisite**
- ✓ The microscope is operational.
  - ✓ Phase contrast objectives with the phase rings **PhC 1**, **PhC 2** or **PhC 3** are *installed* [► 56].
  - ✓ Condenser with modulator disk with centerable ring diaphragms **PhC 1**, **PhC 2** and **PhC 3** is *installed* [► 60].

- Procedure**
1. Swivel the phase contrast objective into the beam path (e.g. **Ph1**).
  2. Switch on the annular phase diaphragm on the condenser's revolver disk with the same labeling as the objective (e.g. **Ph1**).
  3. *Replace one eyepiece* [► 55] with an auxiliary microscope.
  4. With the adjusting fixture on the auxiliary microscope, focus the annular phase diaphragm and the phase ring in the objective exit pupil.
  5. Check the centering and the overlap of the lighter annular phase diaphragm (in the condenser) with the darker phase ring (in the objective). Both rings must be centered and overlapping

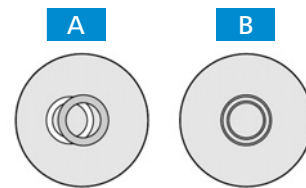

6. If the overlap is not exact **A**, *recenter the lighter annular phase diaphragm* [► 131].
7. Remove the auxiliary microscope and replace the eyepiece.

#### Info

To increase the image contrast, a green 32 x 4 interference broadband filter can be mounted on the field diaphragm or inserted into the color glass carrier (if available).

### 5.4.4 Setting Up for Transmitted Light Polarization

Polarization methods can be used in the transmitted light on the following microscope types:


- Axiolab 5 Pol-TL (430037-9130-000)
- Axiolab 5 Pol-TL/conoscopy (430037-9042-000)
- Axiolab 5 Pol-TL/RL (430037-9032-000)

The following requirements must be fulfilled:

- The microscope is operational.
- Strain-free objectives are installed in the *nosepiece* [► 56].
- The rotary stage is *installed* [► 120].
- A polarizer D (fixed or rotatable) is *installed* [► 128].
- Analyzer slider D, fixed, or lambda compensator or lambda/4 compensator is available.
- A depolarizer for avoiding unwanted polarizing effects is installed in the tube.
- The microscope is adjusted for *transmitted light brightfield microscopy* [► 68].
- The rotary stage is *centered* [► 121].
- The objectives are *centered* [► 122].

#### 5.4.4.1 Detecting Birefringence

For more information about the method, see chapter *Detecting Birefringence* [► 46].

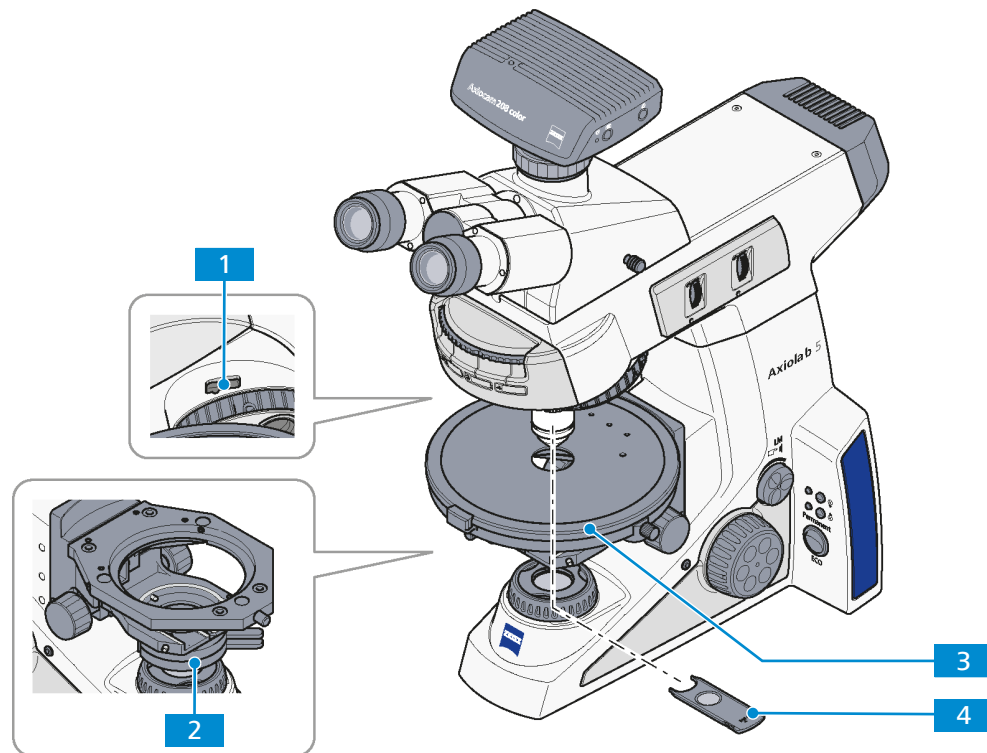

- Procedure**
1. Swing the polarizer **2** into the beam path.
  2. When using a rotatable polarizer, position it to 0°.
  3. Put the analyzer **4** into the slit for the compensator **1**.  
→ The field of view appears dark.
  4. Bring the sample into the field of view.
  5. With the rotary stage **3**, turn the sample.  
→ Normally, birefringent (anisotropic) objects will now show the interference color and intensity variations during rotation between crossed polarizers.  
Optically anisotropic substances may remain dark when an isotropic direction, e.g. from optically single-axle or double-axle crystals, is oriented parallel to the observation direction.

#### 5.4.4.2 Determination of the Polarization Direction

For more information about the method, see chapter *Determination of the Polarization Direction* [▶ 46].

- Prerequisite** ✓ An eyepiece with cross hair reticle is *installed* [▶ 55].  
 ✓ The Pol adjustment sample for polarization microscopy is available.

- Procedure**
1. Swing the polarizer into the beam path.
  2. When using a rotatable polarizer, position it to 0°.
  3. Put the analyzer into the slit for the compensator or swing analyzer module on the reflector turret/slider.  
 → The field of view appears dark.
  4. Place the Pol adjustment sample on the microscope.
  5. Turn the rotary stage until the adjustment sample appears dark.
  6. Remove the analyzer from the beam path.
  7. Align the reticle of the eyepiece along the split cracks of the adjusting sample.
  8. Return the analyzer into the beam path.
  9. Remove the adjustment sample.  
 → The forward direction of the polarizer and analyzer is parallel to the cross hair in the reticle (polarizer east-west, analyzer north-south).
  10. Turn the rotary stage with the sample, e.g. a synthetic fiber, until the sample reaches maximal darkness.  
 → The fiber is parallel to one of the two lines in the cross hair orientation.
  11. Turn the rotary stage by approx. 45° until the longitudinal axis of the fiber is pointing in northeast-southwest direction.  
 → The sample shows the strongest brightness (diagonal position). It can have any color in this position.
  12. Slide in the lambda compensator (possible only if used with screw-in analyzer in tube or intermediate plate).  
 → The sample changes its color depending on its orientation (northeast-southwest or northwest-southeast).

#### 5.4.4.3 Measuring the Path Differences


For more information about the method, see *Measuring Path Differences* [▶ 48].

- Prerequisite** ✓ The correct inter-pupillary distance in the binocular tube is set.

- Procedure**
1. Accurately position the sample to be examined on the center of the eyepiece reticle.
  2. Limit the aperture to a value of about 0.2.
  3. Turn the rotary stage until the sample is almost extinguished, i.e. **completely dark**, and set the 45° locking position.
  4. Rotate the stage **once** (by 45°) so that the sample is in a diagonal position (sample becomes bright).
  5. *Determine* [▶ 48] the suitable compensator.
  6. Insert the determined compensator into the slot as far as it will go.
  7. Use the enclosed operating instructions for measurement preparation and measurement procedure.

#### 5.4.4.4 Circular Polarization Contrast

For more information about the method, see *Circular Polarization Contrast* [▶ 48].

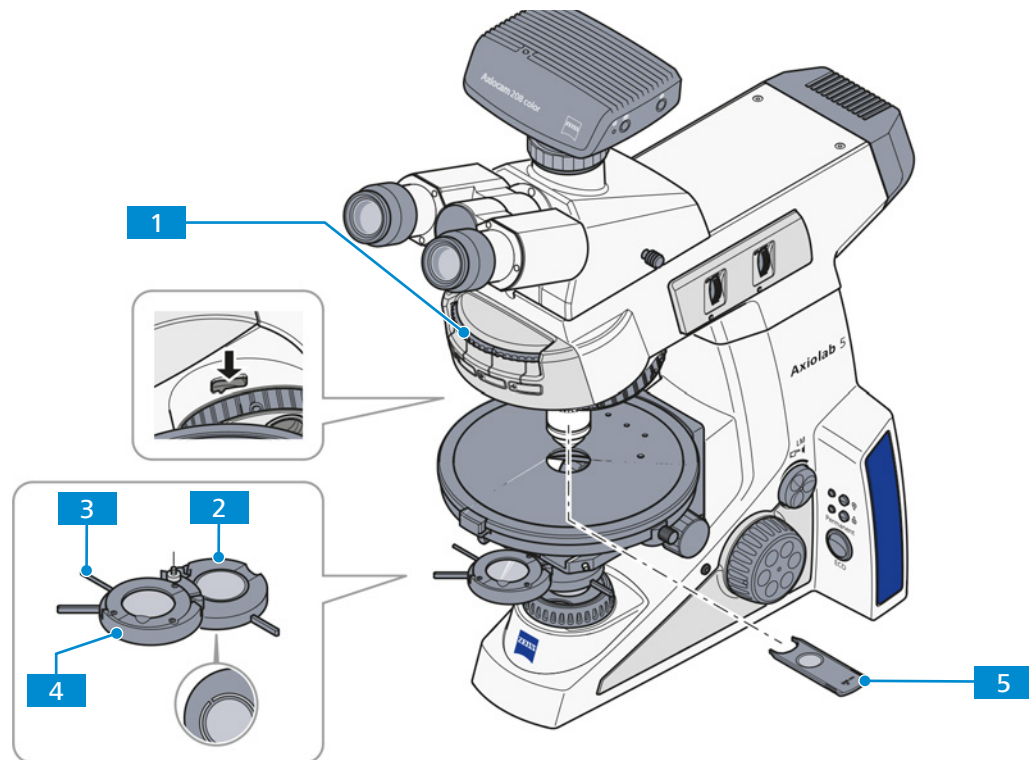

- Prerequisite**
- ✓ The circular polarizer D including the corresponding lambda/4 plate is installed.
  - ✓ The analyzer module is installed in the reflector turret or the stand is equipped with the intermediate plate for analyzer slider and the analyzer slider is available.
  - ✓ The lambda/4 compensator (6x20) is available.

- Procedure**
1. Remove the sample.
    - Initially do not use a sample for the further settings.
  2. Swivel the lower part **2** of the circular polarizer into the light path up to the click stop.
  3. Swing in the analyzer module on the reflector turret **1** or insert the analyzer slider into the intermediate plate.
  4. At full light intensity, assess the extinction (darkening) of the field of view without sample.
  5. Push the lambda/4 compensator **5** into the compensator slot above the nosepiece as far as it will go.
  6. Swivel the upper part **4** of the circular polarizer into the light path.
  7. Rotate the lever of the lambda/4 plate **3** of the D circular polarizer until the field of view appears dark gray.
    - The lever points 45° to the right.
    - The maximum extinction is obtained.
  8. Place the sample to be examined on the stage.
    - The samples appear constantly and independently of the stage rotation in their specific interference color, which depends on the material, sample thickness and orientation.
  9. For the detection of gout or pseudo-gout, select crystal needles that are oriented in the gamma direction (see marking on the lambda plate).
    - If the crystal needles oriented parallel to the gamma direction of the lambda plate are yellow, and the crystal needles lying at a right angle to the gamma direction are blue, the crystals are monosodium urate crystals (gout).

- If the crystal needles oriented parallel to the gamma direction of the lambda plate are blue, and the crystal needles lying at a right angle to the gamma direction are yellow, the crystals are calcium pyrophosphate crystals (pseudo-gout).

#### 5.4.4.5 Transmitted Light Polarization with the Conoscopy Stand

The present section applies to the following microscope type:

- Axiolab 5 Pol-TL/conoscopy (430037-9042-000)

#### NOTICE

##### Risk of mechanical damage

The movements of **rotary knobs A** and **BL** and the respective setting wheels are coupled with each other. Only one control element should therefore be operated at a time and the movement of the other should not be inhibited or blocked. Otherwise, mechanical damage may occur.


- ▶ Operate only one of the 4 controls (rotary knobs, setting wheels) at a time.

#### Info

If the **rotary knob BL** is set to the **On** position, **rotary knob A** is automatically moved into the **On** position if it is not there already. If, on the other hand, **rotary knob A** is set to the **Off** position, **rotary knob BL** is automatically moved into the **Off** position if it is not there already.

##### 5.4.4.5.1 Detecting Birefringence

For more information about the method, see chapter *Detecting Birefringence* [► 46].

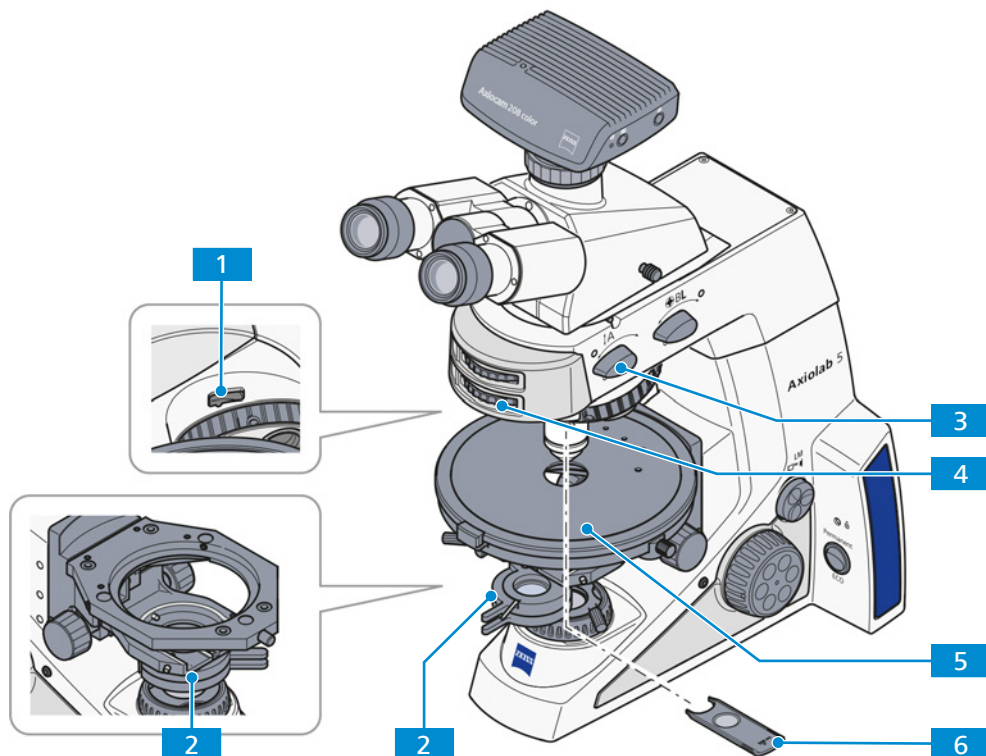


- Prerequisite** ✓ The illumination is adjusted for *transmitted light brightfield microscopy* [► 68] using the KÖHLER method.
- ✓ Lambda compensator or lambda/4 compensator is available.



- Procedure**
1. Swing the polarizer **2** into the beam path.
  2. When using a rotatable polarizer, position it to 0°.
  3. Swivel the analyzer into the beam path (on position) with rotary knob **A 3**.
  4. Adjust the analyzer using the setting wheel **4** until the field of view is dark.
  5. Bring the sample into the field of view.
  6. With the rotary stage **5**, turn the sample.
    - Normally, birefringent (anisotropic) objects will now show the interference color and intensity variations during rotation between crossed polarizers. Optically anisotropic substances may remain dark when an isotropic direction, e.g. from optically single-axle or double-axle crystals, is oriented parallel to the observation direction.
  7. If required, put the lambda compensator or lambda/4 compensator **6** into the slit for the compensator **1** for detecting birefringence.

#### 5.4.4.5.2 Determination of the Polarization Direction

For more information about the method, see chapter *Determination of the Polarization Direction* [▶ 46].

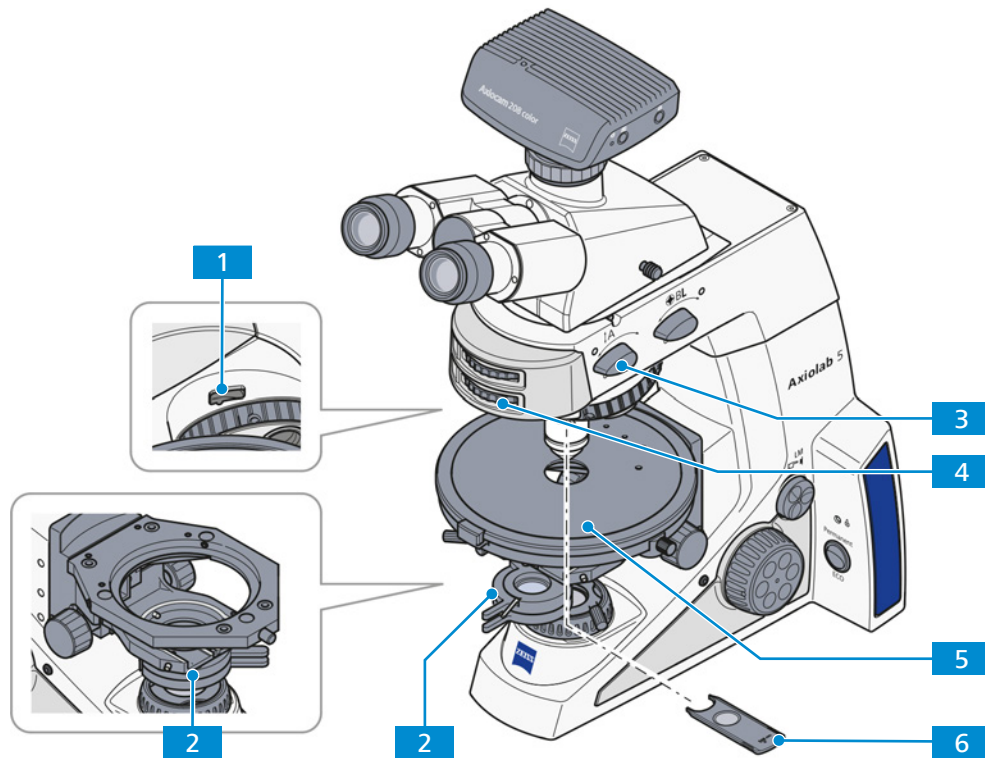

- Prerequisite**
- ✓ An eyepiece with cross hair reticle is *installed* [▶ 55].
  - ✓ The Lambda or lambda/4 compensator are available.
  - ✓ The Pol adjustment sample for polarization microscopy is available.

- Procedure**
1. Swing the polarizer **2** into the beam path.
  2. When using a rotatable polarizer, position it to 0°.
  3. Swivel the analyzer into the beam path (on position) with rotary knob **A 3**.
  4. Adjust it using the setting wheel **4** until the field of view is dark.
  5. Place the Pol adjustment sample on the microscope.
  6. Turn the rotary stage **5** until the adjustment sample appears dark.

7. Swivel out the analyzer **3**.
8. Align the reticle of the eyepiece along the split cracks of the adjusting sample.
9. Return the analyzer **3** into the beam path.
10. Remove the adjustment sample.
  - The forward direction of the polarizer and analyzer is parallel to the cross hair in the reticle (polarizer east-west, analyzer north-south).
11. Turn the rotary stage with the sample, e.g. a synthetic fiber, until the sample reaches maximal darkness.
  - The fiber is parallel to one of the two lines in the cross hair orientation.
12. Turn the rotary stage by approx. 45° until the longitudinal axis of the fiber is pointing in northeast-southwest direction.
  - The sample shows the strongest brightness (diagonal position). It can have any color in this position.
13. Slide in the lambda compensator or lambda/4 compensator **6** in the slit **1**.
  - The sample changes its color depending on its orientation (northeast-southwest or northwest-southeast).

#### 5.4.4.5.3 Measuring the Path Differences

For more information about the method, see *Measuring Path Differences* [► 48].


- Prerequisite**
- ✓ The illumination is adjusted for *transmitted light brightfield microscopy* [► 68].
  - ✓ The correct *inter-pupillary distance* [► 64] in the binocular tube is set.

- Procedure**
1. Accurately position the sample to be examined on the center of the eyepiece reticle.
  2. Limit the aperture to a value of about 0.2.
  3. Turn the rotary stage until the sample is almost extinguished, i.e. **completely dark**, and set the 45° locking position.
  4. Rotate the stage **once** (by 45°) so that the sample is in a diagonal position (sample becomes bright).
  5. *Determine* [► 48] the suitable compensator.
  6. Insert the determined compensator into the slot as far as it will go.
  7. Use the enclosed operating instructions for measurement preparation and measurement procedure.



#### 5.4.4.5.4 Determining the Optical Character of Crystals

For more information about the method, see chapter *Determining the Optical Character of Crystals* [▶ 49].

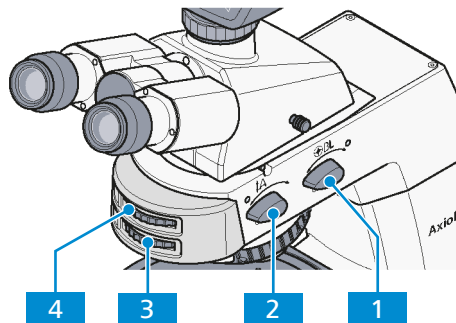


- Prerequisite**
- ✓ A strain-free objective is screwed in, recommended:
    - N-Achroplan 50x/0.8 Pol objective or
    - EC Plan-Neofluar 40x/0.9 Pol objective
  - ✓ A suitable condenser 0.9 Pol is installed.
  - ✓ The illumination is adjusted for *transmitted light brightfield microscopy* [▶ 68].

- Procedure**
1. Place the sample on the stage.
  2. Focus the sample.
  3. Swivel the analyzer into the beam path (on position) with rotary knob **A** 2.
  4. If necessary, change the polarization direction of the analyzer using the setting wheel 3.
  5. Place a selected crystal in the center of the crossline reticle.
  6. Swivel in the N-Achroplan 50x/0.8 Pol objective or EC Plan-Neofluar 40x/0.9 Pol objective and focus using the focusing drive.
  7. If necessary, close the luminous-field aperture to avoid superimposition of axial figures of neighboring crystals on the axial figure to be examined.
    - The smallest crystal value that can be faded out is 170 µm.
  8. If necessary, switch on Bertrand lens **BL** 1.
    - The axial figure will appear in the field of view.
  9. Bring the axial figure into focus using the setting wheel 4.
  10. Determine the optical character of the crystal.

#### 5.4.4.5.5 Circular Polarization Contrast

For more information about the method, see *Circular Polarization Contrast* [▶ 48].

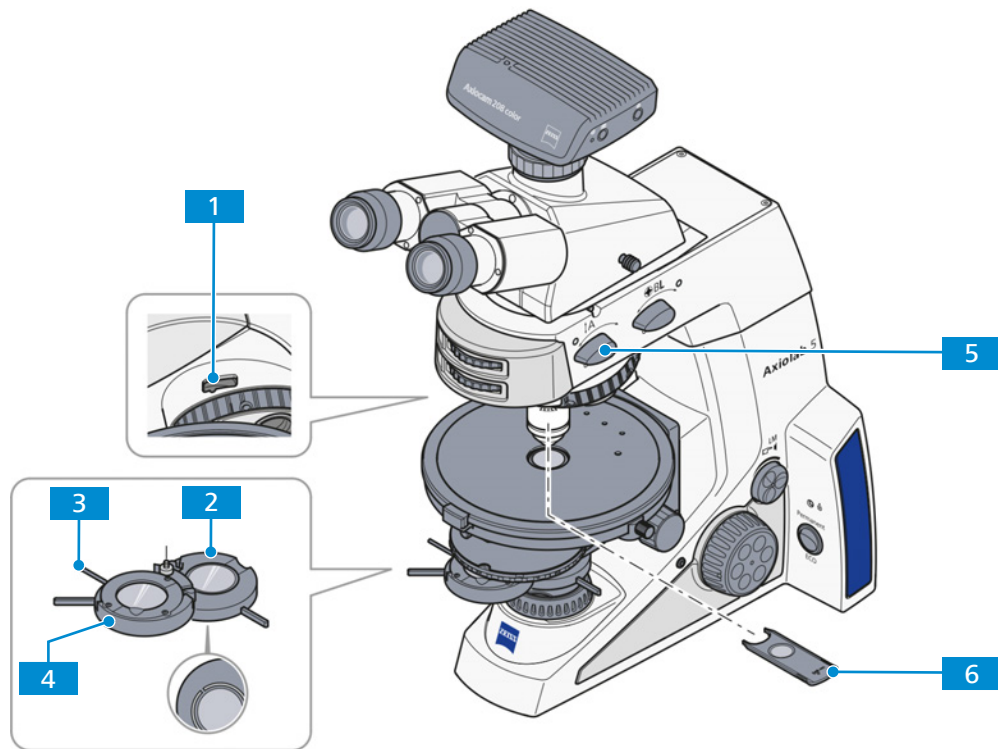

- Prerequisite**
- ✓ A strain-free objective is screwed in, recommended:
    - N-Achroplan 50x/0.8 Pol objective or
    - EC Plan-Neofluar 40x/0.9 Pol objective
  - ✓ The circular polarizer D including the corresponding lambda/4 plate is installed.
  - ✓ The lambda/4 compensator (6x20) is available.
  - ✓ The illumination is adjusted for *transmitted light brightfield microscopy* [▶ 68].

- Procedure**
1. Remove the sample.
    - Initially do not use a sample for the further settings.
  2. Swivel the analyzer with rotary knob **A** **5** into the beam path.
  3. Swivel the lower part **2** of the circular polarizer into the light path up to the click stop.
  4. At full light intensity, assess the extinction (darkening) of the field of view without sample.
  5. If this is not optimal, align the analyzer as necessary.
  6. Push the lambda/4 compensator **6** into the compensator slit **1** above the nosepiece as far as it will go.
  7. Swivel the upper part **4** of the circular polarizer into the light path.
  8. Rotate the lever of the lambda/4 plate **3** of the circular polarizer until the field of view appears dark gray.
    - The lever points 45° to the right.
    - The maximum extinction is obtained.
  9. Place the sample to be examined on the stage.
    - The samples appear constantly and independently of the stage rotation in their specific interference color, which depends on the material, sample thickness and orientation.
  10. For the detection of gout or pseudo-gout, select crystal needles that are oriented in the gamma direction (see marking on the lambda plate).

- If the crystal needles oriented parallel to the gamma direction of the lambda plate are yellow, and the crystal needles lying at a right angle to the gamma direction are blue, the crystals are monosodium urate crystals (gout).
- If the crystal needles oriented parallel to the gamma direction of the lambda plate are blue, and the crystal needles lying at a right angle to the gamma direction are yellow, the crystals are calcium pyrophosphate crystals (pseudo-gout).


## 5.5 Setting Up for Reflected Light Techniques

### 5.5.1 Setting Up for Reflected Light Brightfield Microscopy


For more information about the method, see chapter *Reflected Light Brightfield Microscopy Using the KÖHLER Method* [► 50].

- Prerequisite**
- ✓ A light source for reflected light is available on the stand.
  - ✓ In the reflector turret, an ACR P&C brightfield reflector module for reflected light is installed.
  - ✓ The microscope is *adapted* [► 64] to the user.
  - ✓ The microscope is operational.


- Procedure**
1. If required, push the **RL** button for reflected light illumination.

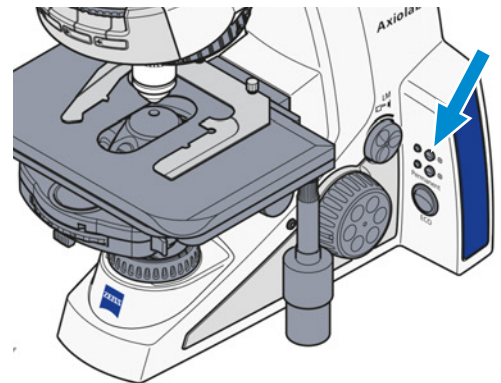


2. Adjust the image brightness using the **Intensity/LM** knob on the microscope stand.

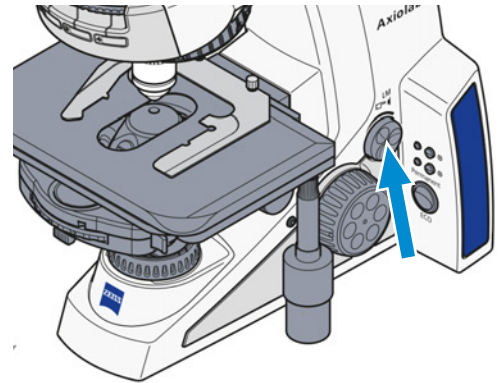


3. Place a high-contrast reflected light sample into the sample holder of the mechanical stage.
4. Swivel the 10x objective into the beam path.

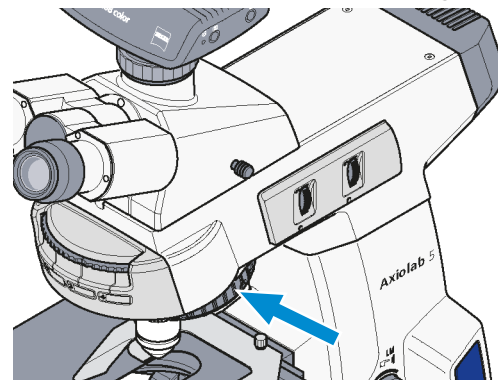


5. Swivel in the position with the brightfield reflector module on the reflector turret.

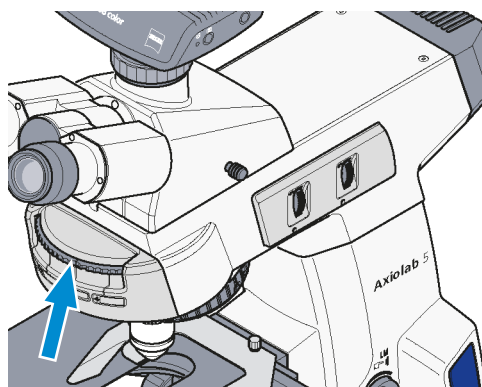


6. Focus the sample.  
Try to focus away from the sample to avoid any collision between the objective and sample.

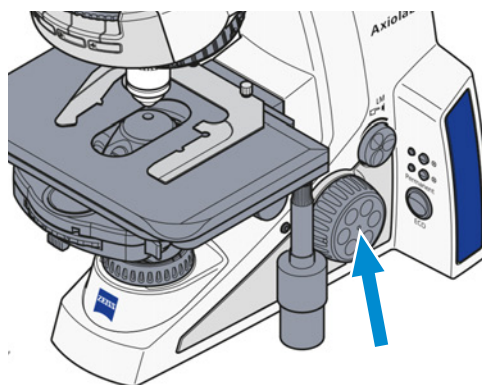


7. Set the knurled wheel of the aperture diaphragm **A** to a medium position.

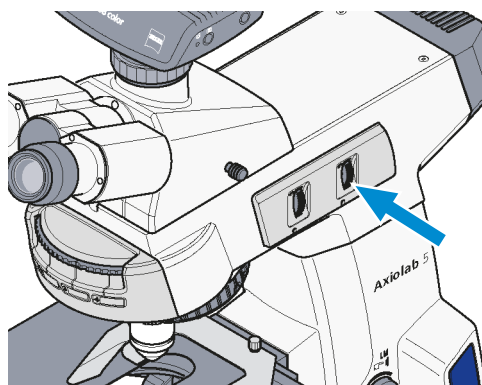


8. Adjust the knurled wheel of the luminous-field diaphragm **F** so that the field diaphragm becomes visible in the field of view.

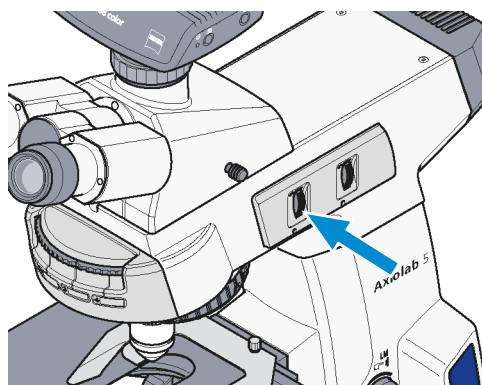


9. Focus on the edge of the luminous-field diaphragm using the focusing drive.

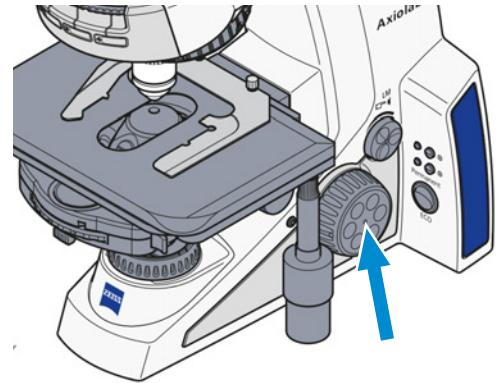

10. Open the luminous-field diaphragm enough to make the edge of the diaphragm disappear from the field of view.

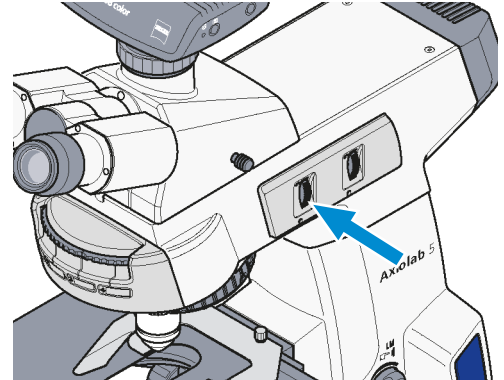

11. Remove an eyepiece from the binocular tube to adjust the aperture diaphragm (contrast).
12. Look into the tube with the naked eye or use the auxiliary microscope instead of the eyepiece.
13. Set the aperture diaphragm with the knurled wheel to between 2/3 and 4/5 of the exit pupil diameter of the objective.  
In most cases this aperture gives the best contrast at almost full resolution and is thus the best compromise for the human eye.

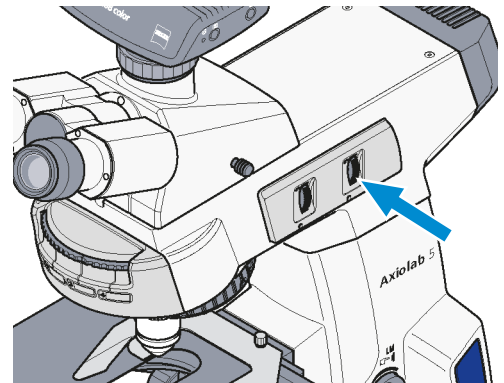

14. Replace the eyepiece.
  15. Remove the high-contrast reflected light sample.
  16. Re-adjust the focus using the coarse and fine focusing drives and set the image brightness according to the reflected light sample.
  17. Readjust the aperture diaphragm diameter after each objective change.
- ↳ The illumination is now adjusted according to the KÖHLER method.

### Info


Never use the aperture diaphragm for controlling image brightness. Use the **Intensity/LM** knob for illumination intensity!

### 5.5.2 Setting Up for Reflected Light Darkfield Microscopy


For more information about the method, see chapter *Reflected Light Darkfield Microscopy Using the KÖHLER Method* [► 51].

- Prerequisite**
- ✓ A light source for reflected light is available on the stand.
  - ✓ In the reflector turret, an ACR P&C brightfield reflector module for reflected light is installed.
  - ✓ A suitable objective for RL darkfield microscopy is installed.  
e.g. Epiplan-Neofluar, EC Epiplan-Neofluar, Epiplan objectives with the additional designation "HD"
  - ✓ The microscope is operational.
  - ✓ The microscope is adjusted for *reflected light brightfield microscopy* [► 83].
  - ✓ The displayed luminous-field aperture should be located slightly beyond the edge of the field of view.


- Procedure**
1. If used, remove the 6x20 mm compensator.
  2. Swivel in the objective position with the dark-field objective (HD) on the nosepiece.

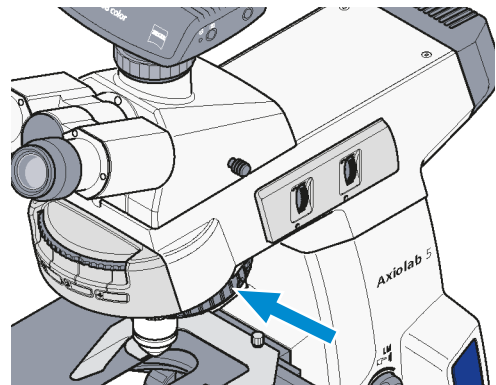

3. Swivel in the darkfield reflector module on the reflector turret.

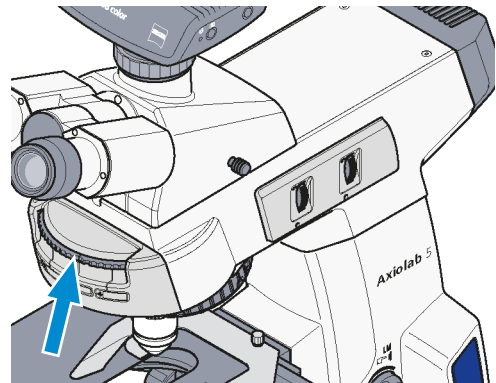


4. Completely open the aperture diaphragm **A**.

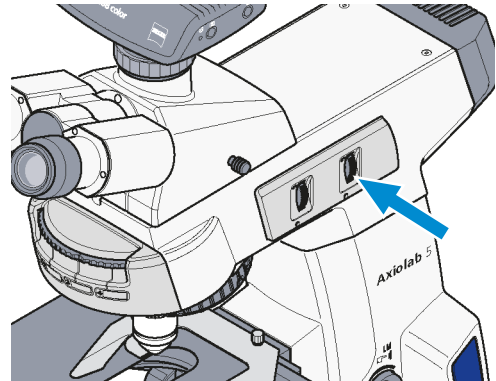

5. Switch off or remove the neutral filter as necessary.
6. Place the sample on the stage.



7. Focus the sample.

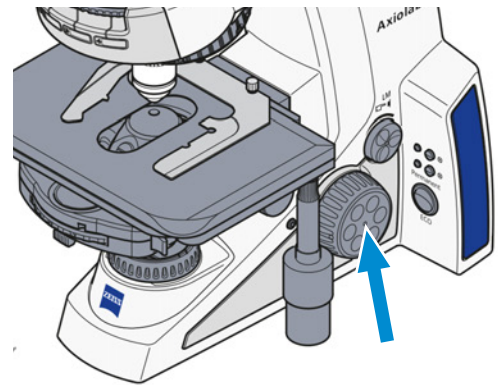


↳ The illumination is now adjusted for darkfield microscopy.

### 5.5.3 Setting Up for Reflected Light Polarization Microscopy


For more information about the method, see chapter *Reflected Light Polarization Microscopy* [► 51].

- Prerequisite**
- ✓ A light source for reflected light is available on the stand.
  - ✓ In the reflector turret, an ACR P&C DIC, C-DIC or Pol reflector module for reflected light is installed.
  - ✓ A suitable objective for RL polarization microscopy is installed.  
e.g. Epiplan-Neofluar Pol, EC Epiplan-Neofluar Pol, Epiplan Pol objectives with the additional designation "HD"
  - ✓ The microscope is operational.
  - ✓ The microscope is adjusted for *reflected light brightfield microscopy* [► 83].
  - ✓ A analyzer slider D with lambda plate or lambda compensator or lambda/4 compensator (6x20) is available.


- Procedure**
1. If required, push the **RL** button for reflected light illumination.

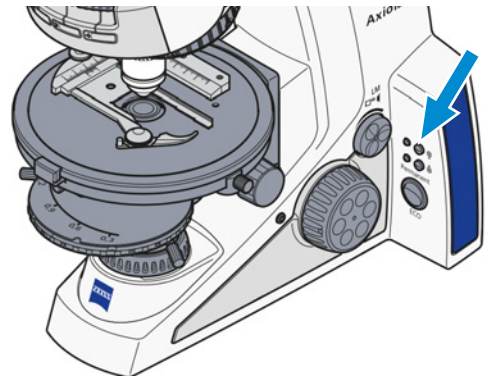


2. Swivel in the P&C (for DIC or Pol) reflector module on the reflector turret.

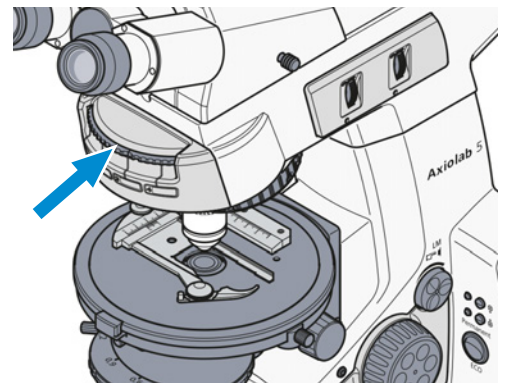


3. Insert the analyzer slider (or lambda compensator or compensator lambda/4) into the 6x20mm slot.

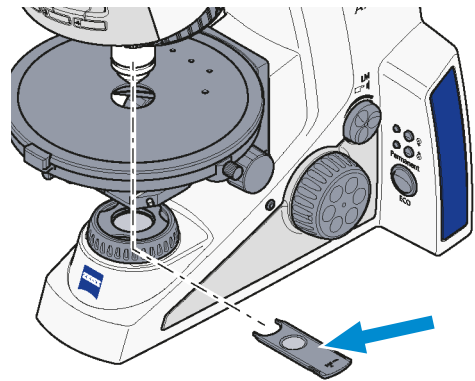

4. Place the sample on the stage.
5. Swivel in the Pol objective with the desired magnification level.

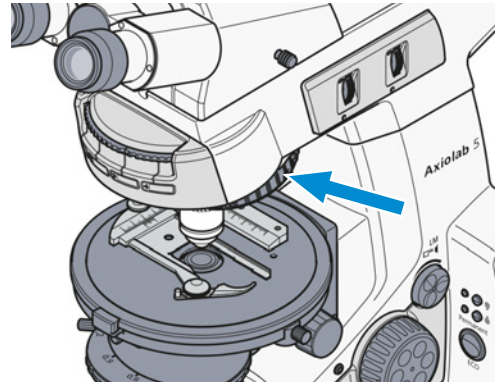

6. Focus the sample.

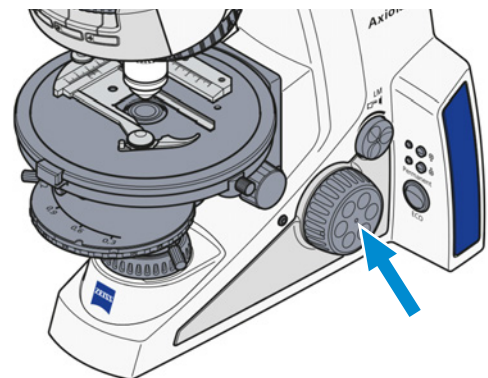

7. Observe the sample in the polarization contrast now present while turning the Pol rotary stage.

↳ The illumination is now adjusted for polarization microscopy.



### 5.5.4 Setting Up for Reflected Light Fluorescence Microscopy

The present section applies to the following microscope type:

- Axiolab 5 Bio-TL/FL (430037-9021-000, 430037-9120-000, 430037-9070-000)

For more information about the method, see chapter *Reflected Light Fluorescence Microscopy* [▶ 51].

#### WARNING

##### Skin or eye injury due to hazardous light emission

The light source belongs to Risk Group 3 as specified in IEC 62471 and emits LED radiation and UV radiation. Skin or eye injury can result from the exposure.

- ▶ Avoid any eye and skin exposure to the light-emitting aperture of the light source.
- ▶ Avoid exposure of skin to the radiation. Use suitable protective equipment/protective clothing if required.
- ▶ Before installing or removing the light source always make sure it is switched off.

#### NOTICE

##### Property damage due to heat emission

Microscope lamps emit a lot of heat which could damage the heat-sensitive fluorescence filters.


- ▶ Do not remove the heat protection filter when using a fluorescence filter.

#### Info


The adjustment of reflected light fluorescence is facilitated by starting with an objective of average magnification, e.g. EC Plan-Neofluar 20x/0.50, and a sample of high fluorescence. Demonstration samples can also be used for the start-up.

- Prerequisite**
- ✓ In the reflector turret, FL P&C reflector modules equipped with respective filter sets are installed.
  - ✓ Fluorescence protection shield is available.
  - ✓ A suitable objective for fluorescence microscopy is installed.  
e.g. EC Plan-Neofluar or Fluar (UV excitation)
  - ✓ The microscope is operational.
  - ✓ The microscope is adjusted for *reflected light brightfield microscopy* [▶ 83].


- Procedure**
1. Slide the fluorescence protection shield into the compensator slit above the nosepiece.

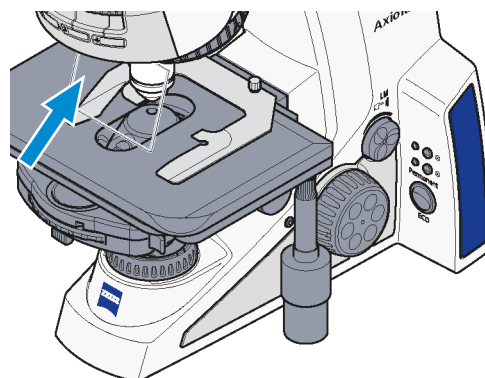


2. Swivel in the EC Plan-Neofluar objective on the nosepiece.

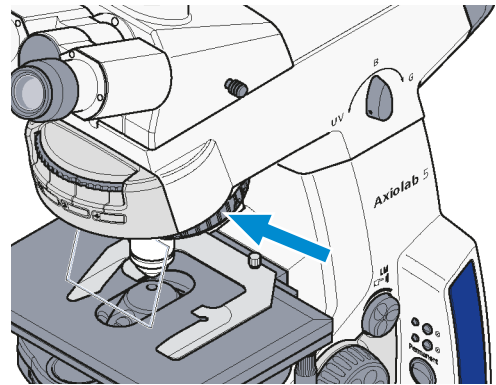


3. Initially set transmitted light illumination by pushing the **TL** button.

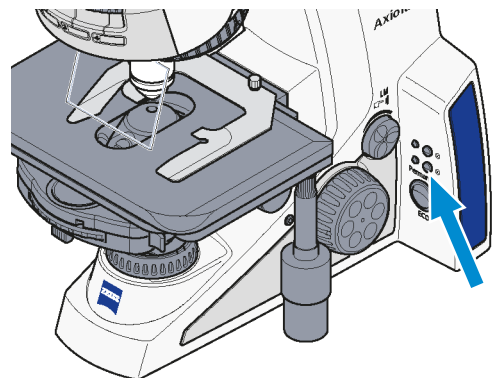


4. If necessary, turn the condenser turret to the **H** (B) position for transmitted light brightfield (or phase contrast if using a Ph objective).

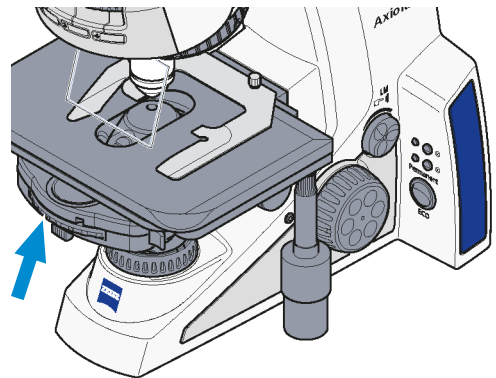


5. Seek the sample detail to be examined.
6. Adjust the light intensity by turning the **Intensity/LM** knob.

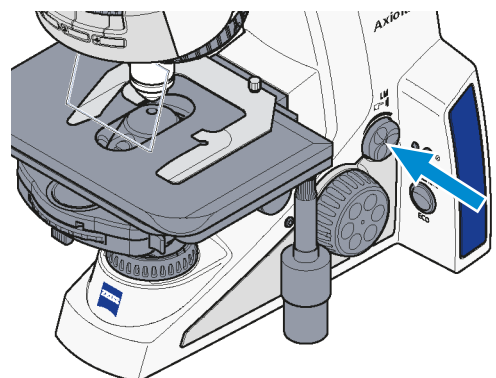

7. Focus the sample.

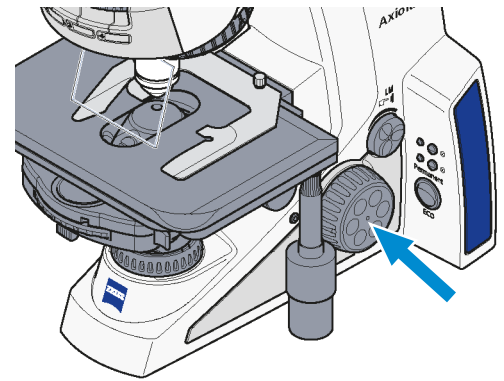


8. Swivel it in the FL P&C reflector module with the desired fluorescence filter combination (depending on the excitation mode).

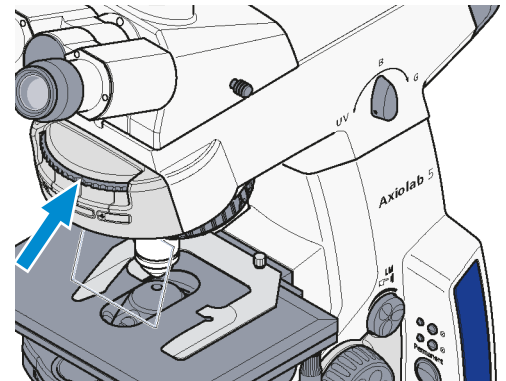


9. Swivel in the desired LED (UV, B or G) using the LED selection knob.

**CAUTION** To avoid dazzling when switching between the LEDs, the brightness should be reduced slightly beforehand.

**NOTICE** When switching between the three LEDs, the current brightness setting is adopted.

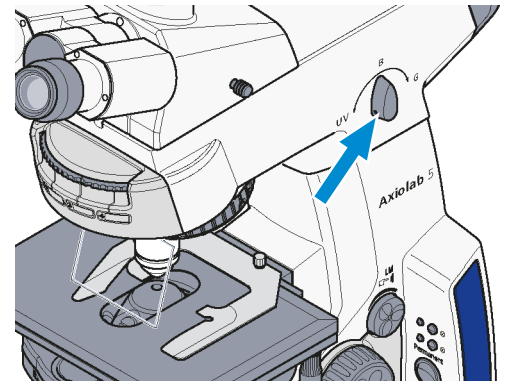

10. Set reflected light illumination by pushing the RL button.

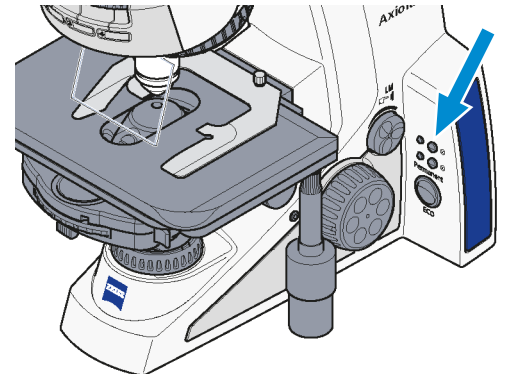


11. Adjust the light intensity for reflected light by turning the **Intensity/LM** knob.
12. Focus the sample.

↳ The illumination is now adjusted for fluorescence microscopy.

## 5.6 Switching Off the Microscope

**Procedure** 1. Switch the microscope off using the power switch **On/Off**.

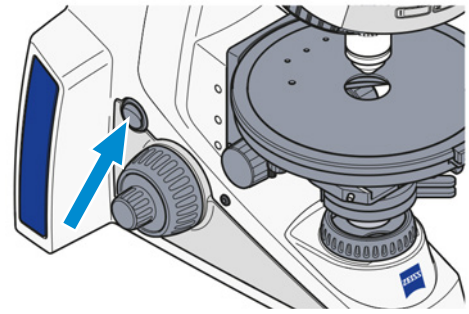

2. Cover the microscope with the dust cover.

## 6 Care and Maintenance

To ensure the best possible performance of the microscope and its components, maintenance must be performed on a regular basis. Keep the service logs of the microscope.

To maintain operational safety and reliability of the microscope, we recommend entering into a **ZEISS Protect Service Agreement**.

### Info

For additional information and detailed descriptions, refer to further applicable documents or ask your ZEISS Sales & Service Partner.

### 6.1 Safety during Cleaning and Maintenance

Only conduct preventive measures described here. All tasks of maintenance and cleaning not described may only be performed by an authorized ZEISS service representative.

Any unauthorized intervention or any operation outside the scope of the intended use can lead to injuries and property damage and voids all rights to warranty claims. Only original spare parts from ZEISS may be used.



### DANGER

#### Electric injury due to live parts

When the microscope and its components are still switched on, coming in contact with live parts can lead to electric shock or burn.

- ▶ Switch off the microscope and its components prior to opening or cleaning.
- ▶ Disconnect live parts from the power supply.

### NOTICE

#### Functional impairment due to dirt and moisture

Dirt, dust, and moisture can impair the microscope functionality and can cause short-circuits.

- ▶ Use the dust protection cover if the microscope is not used.
- ▶ The ventilation slots must be unobstructed at all times.
- ▶ Perform regular maintenance and cleaning according to the instructions in this document and according to the instructions in the applicable documents.
- ▶ Make sure that no cleaning liquid or moisture gets inside the microscope and its components.
- ▶ In case of damage, the affected parts of the microscope must be taken out of operation.

## 6.2 Maintenance Schedule

To maintain best possible performance of the microscope, it is essential to perform preventive maintenance on a regular basis. The recommended intervals depend on the total uptime of the microscope.

Interval	Part/Component	Activity
daily	Microscope	<p>Check the power cable and the plug for possible damage.</p> <p>If any damage is observed, turn the microscope off and secure it against inadvertent restarts immediately.</p> <p>Call in a qualified professional to remedy the problem.</p>

## 6.3 Maintenance Work

### 6.3.1 Cleaning an Optical Surface


#### NOTICE

##### Damage of optical surfaces due to improper cleaning

- ▶ Remove dust from the optical surface slowly and carefully.
- ▶ Remove dust on optical surfaces with a natural-hair brush or blow it off with a rubber bellows.
- ▶ Avoid touching optical surfaces with fingers.

- Parts and Tools**
- 🔧 Clean cloth
  - 🔧 Cotton swab
  - 🔧 Optical cleaning solution
  - 🔧 Lint-free cloth

- Procedure**
1. Moisten a cotton swab or a clean cloth with an optical cleaning solution, if necessary.
  2. Wipe optical surfaces in a circular motion towards the edge of the optics with slight pressure.

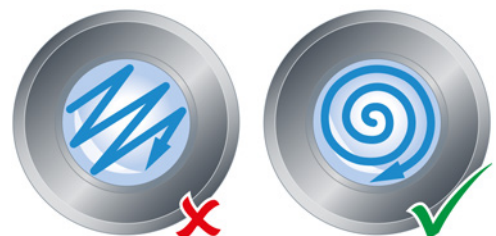

3. Dry with a lint-free cloth.

### 6.3.2 Removing Water-Soluble Contamination

#### NOTICE

##### Damage of the microscope due to wrong solutions or solvents

During cleaning, incorrectly used or wrong solutions/solvents can harm the microscope.

- ▶ Test clean solutions of unknown composition first on a less visible area of the unit.
- ▶ Use only tested solutions/solvents.

**Parts and Tools**

- 🔧 Clean cloth
- 🔧 Lint-free cloth

**Prerequisite** ✓ The microscope and its components are switched off and disconnected from the power supply.

- Procedure**
1. Remove dust and loose dirt particles with a soft brush or clean lint-free cloth.
  2. If necessary, moisten a clean cloth with water.
    - ➔ Stubborn dirt can be cleaned with all commercially available water solutions, benzine or alcohol (no solvent!). For cleaning coated parts, use a linen or leather cloth that is moistened with one of these substances.
  3. Wipe off the area with the cloth.
  4. Dry with a lint-free cloth.

**Info** Labels on the device may only be cleaned using a dry cloth.


### 6.3.3 Exchanging the Fuses T 15 A/H 250 V in the Stand

#### ⚠ DANGER

##### Electric injury due to live parts

When the microscope and its components are still switched on, coming in contact with live parts can lead to electric shock or burn.

- ▶ Switch off the microscope and its components prior to opening or cleaning.
- ▶ Disconnect live parts from the power supply.

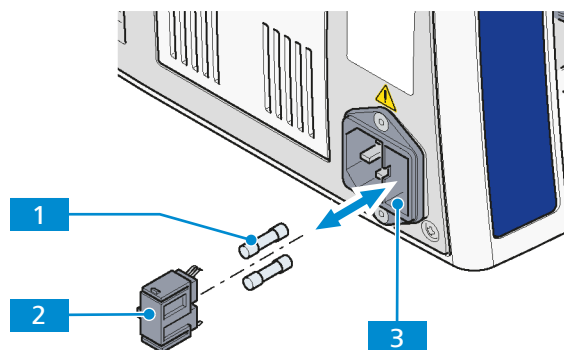

**Parts and Tools** 🔧 2x Fuse type T 15 A/H 250 V

- Prerequisite** ✓ The microscope is switched off.  
 ✓ The microscope is disconnected from the mains.

- Procedure**
1. If the fuses fail, first check the cause and remedy technical problems properly.
  2. Remove the fuse holder **2** on the rear side of the stand.
  3. Remove fuses **1** from the fuse holder.
  4. Insert new fuses.
  5. Push the fuse holder back into the fuse compartment **3** until it locks in place.
  6. Bring the microscope back into operation.


### 6.3.4 Mechanical Stages

Axiolab 5 stands are fitted with the respective mechanical stage at the factory according to customer requirements.

The friction adjustment of the coaxial knurled knobs is set at an average value at the factory.

Should the stage need to be replaced or the stage settings changed, proceed as follows.


#### 6.3.4.1 Adjusting the Drive Length on the Stage Drive

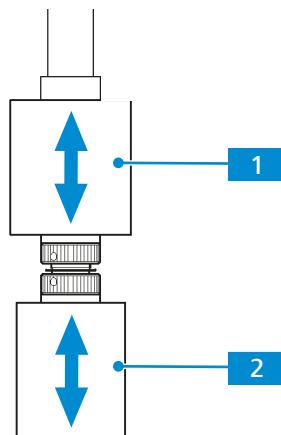

- Procedure**
1. Shift or lower the Coaxial knurled knob **1** for the Y adjustment within a range of approx. 15 mm.
  2. Shift or lower the Coaxial knurled knob **2** for the X adjustment within a range of approx. 15 mm.



### 6.3.4.2 Adjusting the Friction of the Coaxial Knurled Knobs on the Stage Drive

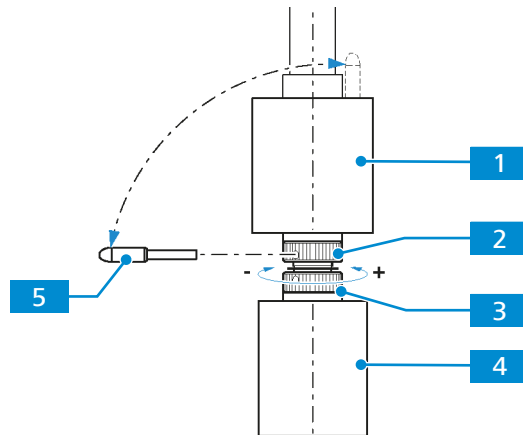

#### X drive

- Procedure**
1. Push the coaxial knurled knob for the X adjustment **4** all the way to the bottom.
  2. Remove the supplied adjusting pin **5** from the coaxial knurled knob for the Y adjustment **1**.
  3. Insert it into one of the holes of the lower hole nut **3**.
  4. Hold the coaxial knurled knob for the X adjustment **4** and turn the hole nut with the adjusting pin clockwise or counter-clockwise until the desired freedom of movement has been achieved.
    - Small friction adjustment: – (clockwise)
    - Large friction adjustment: + (counter-clockwise)
    - It should not be shifted more than one revolution.

#### Y drive

- Procedure**
1. Push the coaxial knurled knob for the Y adjustment **1** all the way to the top.
  2. Insert the supplied adjusting pin **5** into the hole of the upper hole nut **2**.
  3. Hold the coaxial knurled knob for the Y adjustment **1** and turn the hole nut with the adjusting pin clockwise or counter-clockwise until the desired freedom of movement has been achieved.
    - Small friction adjustment: – (clockwise)
    - Large friction adjustment: + (counter-clockwise)
    - It should not be shifted more than one revolution.
  4. Re-insert the adjusting pin into the coaxial knurled knob for the Y adjustment.

## 7 Troubleshooting

The following table provides hints for solving common problems. If you cannot solve the problem or if you are unsure about a certain technical difficulty, contact your local ZEISS service representative.

Symptom	Cause	Measure
No illumination light after switching on the microscope	Nosepiece and/or reflector turret are not engaged to defined positions.	Move the nosepiece and/or reflector turret to the left or right to engage the nosepiece and/or reflector turret to defined positions.
Shadows or inhomogeneous image brightness in the field of view; the field is not entirely visible	The vis/phot push-pull rod/shift knob on the photo tube is not in the correct functional position (intermediate position).	Move the vis/phot push-pull rod/shift knob to the correct functional position (end position).
	The nosepiece with objective is not engaged in the click-stop.	Push in nosepiece with objective until it engages.
	The condenser is not correctly adjusted.	<i>Set the condenser [▶ 68] correctly (adjustment, centering).</i>
	The aperture diaphragm is not correctly adjusted.	<i>Set the aperture diaphragm [▶ 68] correctly (opening).</i>
	The luminous-field aperture is not correctly adjusted.	<i>Set the luminous-field diaphragm [▶ 68] correctly (opening).</i>
Asymmetric image sharpness, e.g. one side is sharp, one side is blurred	The filter has not been inserted correctly in the filter mount.	Insert filter correctly in the filter mount.
	The condenser is not correctly adjusted.	<i>Set the condenser [▶ 68] correctly.</i>
	The nosepiece is not engaged in its locking position.	Engage the nosepiece in its locking position (click-stop).
Distinct focus differences when changing the objective	The sample has not been fixed on the mechanical stage.	Insert and fix the sample correctly in the sample holder.
	Focusable eyepieces are not adjusted correctly.	<i>Adjust the focusable eyepieces [▶ 64] according to the vision defect.</i>
	A Pol eyepiece was used in a binocular tube without upright image.	Use a matching eyepiece.
	The objective is not screwed in all the way.	Screw the objective in to the stop.
	The tube lens is not mounted, or mounted unnecessarily.	Mount the tube lens or remove it, as appropriate. If necessary, call in service personnel for check/repair.

Symptom	Cause	Measure
Low resolving power and poor image contrast	The aperture diaphragm is not correctly adjusted.	<i>Set the aperture diaphragm [▶ 68] as per 2/3 rule or the sample features.</i>
	The condenser is not correctly focused and the front lens 0.9 is not swivelled in and out correctly.	<i>Focus the condenser [▶ 68] and swivel front lens 0.9 in or out correctly.</i>
	A wrong cover glass thickness for transmitted light objectives corrected for 0.17 mm cover glass thickness is used.	Use standard cover glass with a thickness of 0.17 mm.
	The sample slide is placed up-side down.	Turn the sample slide over; the sample side should be on top.
	Immersion objectives are used with no or non-specified immersion oil.	Use immersion oil 518 N or 518 F from ZEISS.
	The immersion oil contains air bubbles.	Repeat the oiling procedure with fresh oil.
	The front lens of a dry objective is soiled with immersion oil.	Clean the front lens of the dry objective.
	Correction setting is not set to the proper thickness of the cover glass.	Adjust the correction setting ring to the correct cover glass thickness.
Left and the right field of view cannot be brought together in one image	Dirt or dust on the optical surfaces of objectives, eyepieces, condensers or filters.	Clean the respective optical components.
	The eyepiece distance (inter-pupillary distance) is not adjusted correctly.	<i>Adjust the inter-pupillary distance [▶ 64] correctly.</i>
Eye fatigue when using the microscope	The focusable eyepieces are not adjusted correctly.	<i>Adjust the focusable eyepieces [▶ 64] according to the vision defect.</i>
	The eyepiece distance (inter-pupillary distance) is not adjusted correctly.	<i>Adjust the inter-pupillary distance [▶ 64] correctly.</i>
	The focusable eyepieces are not adjusted correctly.	<i>Adjust the focusable eyepieces [▶ 64] according to the vision defect.</i>
	The image brightness is unacceptable.	Adjust the lamp brightness or insert a conversion filter.
	The binocular tube is optically/mechanically misaligned.	Call in service personnel for check/repair.

Symptom	Cause	Measure
Dirt or dust in the field of view	The condenser is not correctly focused and the front lens 0.9 is not swivelled in and out correctly.	<i>Focus the condenser [▶ 68] and swivel front lens correctly in or out.</i>
	The opening of the aperture diaphragm is too small.	<i>Set the aperture diaphragm [▶ 68] as per 2/3 rule or the sample features.</i>
	Dirt or dust on optical surfaces of objectives, eyepieces, condensers, filters or samples.	<i>Clean the optical surfaces [▶ 94] of the soiled components.</i>
LED/halogen lamp does not light up although the switch is in the On position	The power plug is not plugged into the mains outlet.	Insert the plug into the mains outlet. Ensure that the outlet and instrument have the same voltage.
	The lamp is not installed.	<i>Install lamp for transmitted light [▶ 132] or reflected light [▶ 136].</i>
	The lamp is defective.	<i>Replace lamp for transmitted light [▶ 132] or reflected light [▶ 136].</i>
	The fuses are defective.	<i>Replace fuses [▶ 95].</i>
	The installed electrical equipment may be defective.	Call in the service personnel to check components and replace if necessary.
	The power socket does not supply voltage.	Use a different mains power socket.
Halogen lamp flickers, illumination intensity is not stable	The halogen lamp is reaching the end of its service life.	<i>Replace lamp for transmitted light [▶ 132] or reflected light [▶ 136].</i>
	The power cable is not installed properly or is damaged.	Install the power cable properly or replace it.
	The plug of the LED/halogen lamp is not properly inserted in the socket.	<i>Insert the lamp plug for transmitted light [▶ 132] or reflected light [▶ 136] correctly.</i>
No light in eyepiece	The system is in ECO mode.	Turn the <b>Intensity/LM</b> knob clockwise to wake up the system.
	The light intensity is too low.	Turn the <b>Intensity/LM</b> knob clockwise to increase the light.
	The light was turned off by pressing the corresponding <b>RL/TL</b> button.	Press the <b>RL</b> or <b>TL</b> button according to the corresponding indicator (green).
	A wrong or missing reflector module for reflected light is in use.	Check the reflector turret and make sure the correct reflector is in use.

## 7.1 Resetting the Microscope to Factory Settings

### NOTICE


Use this function with caution as it will reset all existing configurations.

The default factory settings are:

- The Light Manager is enabled, but no light intensity values are saved.
- The light intensity is set to the initial minimum value.
- All stored configurations are cleared.

**Prerequisite** ✓ The microscope is operational.

**Procedure** 1. Press and hold down the **Intensity/LM** knob for 20 seconds.

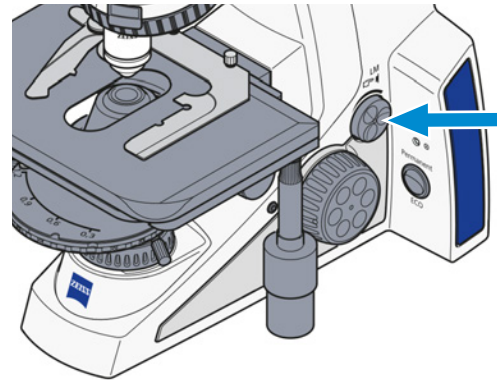

- While the **Intensity/LM** knob is held down from 3 s to 20 s, the indicator light blinks RED.
- It blinks GREEN after 20 s are reached.
- ↳ When the indicator light stops blinking and remains GREEN, the reset to the default factory setting is successful.

## 8 Decommissioning and Disposal

This chapter contains information on the decommissioning and disposal of the microscope and its expansions/components or accessories.

### 8.1 Decommissioning

If the microscope and its components are not used for an extended period of time such as several months, they should be shut down completely and secured against unauthorized access.

#### **DANGER**

##### **Electric injury due to live parts**

When the microscope and its components are still switched on, coming in contact with live parts can lead to electric shock or burn.

- ▶ Switch off the microscope and its components prior to opening or cleaning.
- ▶ Disconnect live parts from the power supply.

- Procedure**
1. Switch off the microscope.
  2. Pull the power supply plug.
  3. Protect microscope using a dust cover.

### 8.2 Transport and Storage

The following regulations must be observed before and during transport:

- The boxes must be secured during transport.
- Avoid rocking the boxes back and forth.
- Note the weight information on the package and on the shipping document.
- Where possible, the original packaging must be used for shipping or transport.
- Do not drop or bump the boxes during movement or storage. Acceleration must not exceed 10 g.
- Evaluate packaging shock and tilting sensors on delivery and after internal transport.

#### **Maximum shock resistance**

#### **Allowable temperature**

Allowable temperature during transport in packaging:

- Between -40 °C and +70 °C
- Relative humidity (without condensation) less than 75 % at 35 °C

Allowable temperature during storage:

- Between +10 °C and +40 °C
- Relative humidity (without condensation) less than 75 % at 35 °C

#### **Info**

**24 hours before installation** of the microscope it is required that the boxes are at recommended room temperature to avoid ingress of humidity, which is harmful to optical paths, and to ensure effective stability of the microscope during installation and testing.

### 8.3 Disposal

The microscope and its components must not be disposed of as domestic waste or through municipal disposal companies. They must be disposed of in accordance with applicable regulations (WEEE Directive 2012/19/EU). ZEISS has implemented a system for the return and recycling of devices in member states of the European Union that ensures suitable reuse according to the EU Directives mentioned.

ZEISS introduced a procedure for the return and recycling of the instruments within the member states of the European Union which ensures suitable recycling procedures conforming to the EU directives.

For more information on disposal and recycling please consult your ZEISS Sales & Service Partner. The microscope may not be disposed of in the household waste or through municipal waste disposal services. If the microscope is resold, the seller shall be obliged to inform the buyer that the microscope must be disposed of in accordance with the regulations.

The customer is responsible for decontamination.

### 8.4 Decontamination

A decontamination statement must be submitted before returning any used objects to the ZEISS location.

If reliable decontamination cannot be guaranteed, the hazard must be marked according to applicable regulations. In general, a well-visible warning sign must be affixed to the article itself and to the outside of the packaging, together with detailed information on the type of contamination.

## 9 Technical Data and Conformity

This chapter contains important technical data as well as information on the conformity.

### 9.1 Performance Data and Specifications

Due to continuous development, we reserve the right to change specifications without notice.

#### Info

Your ZEISS Sales & Service Partner will provide you with the detailed installation requirements.

#### Weight and Sizes

Main Components	Width (mm)	Depth (mm)	Height (mm)	Weight (kg)
Basic stand without tube	approx. 210	approx. 304	approx. 357.5	approx. 8 to 20

#### Location requirements

The microscope may only be operated in closed rooms. The microscope should not be installed near radiators or windows with direct sunlight. The microscope must be placed securely on the table surface to prevent slipping and falling.

Compliance with the installation requirements of the microscope and the availability of the requested supplies is the responsibility of the customer and has to be readily available at the time of installation.

Installation site	Exclusively inside buildings
Altitude	Max. 2000 m above sea level
Atmospheric pressure	Min. 800 hPa

#### Air Conditioning and Quality

Temperature for operation	+10 °C to + 40 °C
Relative humidity (without condensation)	< 75 %
Atmospheric pressure / altitude	800 to 1060 hPa / ≤ 2000 m above sea level
Pollution degree	2
Operational area	closed rooms

#### Mains connection Microscope stand

Protection class	I
Ingress protection rating	IP20 (IEC 60529)
Overvoltage category	II
Nominal AC voltage	100 to 240 V (AC), ±10 %
Nominal frequency	50 to 60 Hz
Power consumption	max. 100 VA
Main power plug	Local mains plug will be supplied.
Addition building PE	The system must be connected to a building earth point at all times.
Fuses in the stand	2x T 3.15 A/H, 5x20 mm (in compliance with IEC 127)



<b>Mains connection Dual observation</b>	Protection class	II		
	Overvoltage category	II		
	Nominal AC voltage (power adapter)	100 to 240 V (AC), $\pm 10\%$		
	Nominal frequency (power adapter)	50/60 Hz		
	Nominal input current	0.2 A		
	Power consumption (Dual observation)	5 VA (5 VDC, 1 A)		
<b>LED illumination TL/RL</b>	Power consumption	max. 10 VA		
	Adjustment of light source	continuous approx. 10 to 800 mA		
<b>Halogen illumination TL/RL</b>	Power consumption	max. 35 VA		
	Adjustment of light source	continuous approx. 0.5 to 12 V		
<b>LED illumination fluorescence TL</b>	Wavelengths optional	385, 470, 505, 565, 590, 625 nm		
<b>Stand specifications</b>	Focusing	manual stage focusing		
	Coarse focusing	approx. 4 mm/revolution		
	Fine focusing	approx. 0.4 mm/revolution; 2 $\mu\text{m}$ scale interval		
	Lifting range	depending on the stand, 15 mm/30 mm		
	Height stop	factory pre-set		
	Objective change	manual		
	Reflector module change	manual		
<b>Tube specifications</b>	<b>Type</b>	<b>Viewing angle</b>	<b>Adjustment</b>	<b>Viewing height* in mm</b>
	Binocular tube 30°/23	30°	- None -	449/485
	Binocular photo tube 30°/23 (50:50)	30°	- None -	449/485
	Binocular photo tube 30°/23 (100:100)	30°	- None -	449/485
	Binocular photo tube 20°/23 (100:100)	20°	- None -	442/481
	Binocular ergo tube 15°/23 (50:50), telescopic, height, upright image	15°	height, telescopic	410/509
	Binocular tube 20°/23	20°	- None -	442/481
	Binocular photo tube 20°/23 Pol (100:100)	20°	- None -	442/481

Type	Viewing angle	Adjustment	Viewing height* in mm
Binocular ergo tube 20°/23 (100/100), reverse image, 44 mm height	20°	height	457/574
Binocular Ergo Photo Tube -2° to 28°/23 (50:50)	-2° to 28°	height	356/507
Binocular Ergo Photo Tube -2° to 28°/23 (50:50)	-2° to 28°	- None -	392/537

\* Range between the lower and upper setting of the eyepieces, e.g. 442/481 → 442 mm to 481 mm

All specifications are for an inter-pupillary distance of 65 mm.

## 9.2 Applicable Standards and Regulations

Observe all general and country-specific safety regulations as well as applicable environmental protection laws and regulations.

The microscope is in compliance with the requirements of the following regulations and directives:

2011/65/EU and delegated directive (EU) 2015/863	Directive 2011/65/EU of the European Parliament and of the Council of 8 June 2011 on the restriction of the use of certain hazardous substances in electrical and electronic equipment (RoHS), amended by Commission Delegated Directive (EU) 2015/863 of 31 March 2015
EN 61010-1:2019	Safety requirements for electrical equipment for measurement, control, and laboratory use – Part 1: General requirements
EN IEC 61326-1:2021	Electrical equipment for measurement, control and laboratory use - EMC requirements - Part 1: General requirements

According to directive 2011/65/EU (RoHS) the microscope and its accessories have been classified as instrument category 9 (Monitoring and control instruments including industrial monitoring and control instruments). They also fall under 2012/19/EU (WEEE).

European and International Directives / Standards: For more information on ISO and CSA certificates or CE Declarations of Conformity, contact your ZEISS Sales & Service Partner.

Only applicable for Axiolab 5 materials

2014/30/EU	Directive 2014/30/EU of the European Parliament and of the Council of 26 February 2014 on the harmonization of the laws of the Member States relating to electromagnetic compatibility
2014/35/EU	Directive 2014/35/EU of the European Parliament and of the Council of 26 February 2014 on the harmonization of the laws of the Member States relating to the making available on the market of electrical equipment designed for use within certain voltage limits

Only applicable for Axiolab 5

(EU) 2017/746	Regulation (EU) 2017/746 of the European Parliament and of the Council of 5 April 2017 on in vitro diagnostic medical devices and repealing Directive 98/79/EC and Commission Decision 2010/227/EU
EN IEC 61010-2-101:2022	Safety requirements for electrical equipment for measurement, control, and laboratory use - Part 2-101: Safety requirements for in vitro diagnostic (IVD) medical equipment
EN IEC 61326-2-6:2021	Electrical equipment for measurement, control and laboratory use - EMC requirements - Part 2-6: Particular requirements - In vitro diagnostic (IVD) medical equipment

## 10 Accessories and System Expansions

Only the following accessories may be used with the microscope as their safe use has been confirmed by ZEISS. Only original parts from ZEISS may be used. Check in advance whether your microscope can be retrofitted with a system expansion or accessories.

After installation or conversion it must be carefully checked whether the microscope and its system expansions/accessories are in a safe operational state and whether unused ports are closed. For details and safety measures refer to the associated documents.

### Info


For additional information and detailed descriptions, refer to further applicable documents or ask your ZEISS Sales & Service Partner.

Name	Description/Info
Objectives	<p>The performance of the microscope objectives affects the image quality of your microscope like no other system component. Whether you work with histological samples, cell samples or entire organisms – the suitability of microscope objectives for your application depends on various factors.</p> <p>More detailed information on available and recommended objectives can be found at <a href="https://www.micro-shop.zeiss.com/de/de/shop/objectives">https://www.micro-shop.zeiss.com/de/de/shop/objectives</a> or ask your ZEISS Sales &amp; Service Partner.</p>
Sliders	<p>The following sliders are available:</p> <ul style="list-style-type: none"> <li>Analyzer slider D/A, with lambda plate, 360° rotatable</li> <li>Analyzer slider D/A, 360° rotatable</li> <li>Analyzer slider D/A, 360° fixed</li> <li>Analyzer slider D/A, with lambda plate, each rotatable +/- 10°</li> <li>Slider 12x46, with focusing Bertrand lens, for phase contrast and conoscopy</li> </ul>
Polarizers	<p>The following polarizers are available:</p> <ul style="list-style-type: none"> <li>Polarizer D, fixed, removable</li> <li>Polarizer D, 90°, rotatable, removable</li> <li>Polarizer, fixed, with lambda plate, rotatable</li> <li>Polarizer, rotatable, with color filter carrier</li> <li>Circular polarizer D</li> <li>Low-power system for objectives 2.5x/4x for condenser 0.9/1.25 H</li> <li>Color filter carrier 3x for filter d=32 mm</li> </ul>
Color filters	<p>The following color filters are available:</p> <ul style="list-style-type: none"> <li>Interference wide-band filter, green, d=32x4</li> <li>Heat absorbing filter KG 1, d=32x2</li> <li>Neutral density filter 0.06, d=32x2</li> <li>Neutral density filter 0.25, d=32x2</li> <li>Polarizing filter 32 mm</li> </ul>

Name	Description/Info
Eyepieces	<p>The following eyepieces and accessories are available:</p> <ul style="list-style-type: none"> <li>▪ Eyepiece E-PL 10x/23 GW, foc.</li> <li>▪ Eyepiece PL 16x/16 GW, foc.</li> <li>▪ Eyepiece PL 10x/23 GW, foc. POL with crossline graticule</li> <li>▪ Eyepiece PL 10x/23 GW, foc.</li> <li>▪ Auxiliary microscope</li> <li>▪ Pinhole diaphragm D= 30 mm</li> </ul>
Condensers	<p>The following condensers are available:</p> <ul style="list-style-type: none"> <li>▪ Ultra condenser 1.2/1.4 (0.75-1.0)</li> <li>▪ Dry darkfield condenser 0.8/0.95 (0.6-0.75)</li> <li>▪ Condenser 0.9/1.25 H</li> <li>▪ Condenser 0.9 H Pol</li> <li>▪ Condenser 0.8 BF WD=5.8mm</li> <li>▪ Condenser, achrom.-aplan. 0.9 BF</li> <li>▪ Condenser, achrom.-aplan. 0.9 BF DF PhC DIC</li> <li>▪ Condenser, achrom.-aplan. 0.9 BF Pol</li> </ul>
Stages	<p>The following stages are available:</p> <ul style="list-style-type: none"> <li>▪ Rackless mechanical stage, 75x30 R</li> <li>▪ Rackless mechanical stage, 75x50 R</li> <li>▪ Rackless mechanical stage, 75x50 L</li> <li>▪ Mechanical stage 75x30 R</li> <li>▪ Pol rotary stage, 360°, with clamping device</li> </ul>
Sample holders	<p>The following sample holders are available:</p> <ul style="list-style-type: none"> <li>▪ Specimen holder for single slide, one-hand operation with spring lever, left</li> <li>▪ Specimen holder for double slides 76x26</li> <li>▪ Attachable Pol object guide, 47x27 mm</li> </ul>
Light sources	<p>The following light sources are available:</p> <ul style="list-style-type: none"> <li>▪ LED module 625 nm for FL</li> <li>▪ LED module 505 nm for FL</li> <li>▪ LED module 470 nm for FL</li> <li>▪ LED module 385 nm for FL</li> <li>▪ LED module 565 nm for FL</li> <li>▪ Illuminator TL/RL LED 10 Axiolab</li> </ul>

Name	Description/Info
Tubes	<p>The following tubes are available:</p> <ul style="list-style-type: none"> <li>▪ Binocular ergophototube 20°/23 (100:0/0:100), reversed image</li> <li>▪ Binocular ergophototube -2° to 28°/23 (50:50), reversed image</li> <li>▪ Binocular ergophototube -2° to 28°/25 (50:50), reversed image</li> <li>▪ Binocular phototube, Pol, 20°/23 (100:0/0:100), upright image</li> <li>▪ Binocular tube 30°/23, reversed image</li> <li>▪ Binocular tube 30°/23, upright image</li> <li>▪ Binocular phototube, 30°/23 (50:50), reversed image</li> <li>▪ Binocular phototube, 30°/23 (100:0/0:100), reversed image</li> <li>▪ Binocular phototube, 20°/23 (100:0/0:100), upright image</li> <li>▪ Binocular phototube, 30°/23 (98:2), reversed image</li> <li>▪ Binocular phototube 30°/23 (100:0/30:70/0:100), reversed image</li> </ul>
Cameras	<p>The following cameras and accessories are available:</p> <ul style="list-style-type: none"> <li>▪ Axiocam 202 mono (see separate instruction manual)</li> <li>▪ Axiocam 208 color (see separate instruction manual)</li> <li>▪ Camera adapter 60N-C 2/3" 0.5x</li> <li>▪ Camera adapter 60N-C 2/3" 0.63x</li> <li>▪ Camera adapter 60N-C 1" 1.0x</li> <li>▪ Video adapter 60 C 1/3" 0.4x</li> </ul>

## 10.1 Assembling the Reversible Eyecups

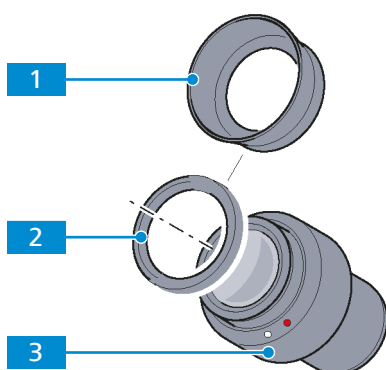

- Procedure**
1. Remove the eyeglass protection ring **2** from the eyepiece **3**.
  2. Mount the reversible eyecup **1**.

Sometimes the eyeglass protection rings are seated very tightly in the eyepiece groove, so you may need a blunt sample (wooden stick) to prod them off.

## 10.2 Binocular Tube

Various tubes with different inclination angles enable suitable eye levels to be selected for observation.

### 10.2.1 Binocular Tube 30°/23

**Purpose** Binocular tubes are used to visualize the microscopic image by means of the eyepieces.

**Position** The binocular tubes are mounted on the top of the stand.

**Function** The pupil distance and the viewing height can be adjusted with the aid of the binocular section. The following features and controls are available:

- Optionally with upright or reversed image
- viewing angle 30°
- field of view 23 mm

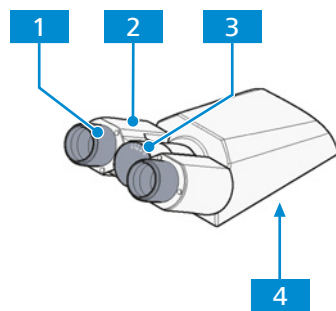


Fig. 33: Binocular Tube 30°/23

- |                          |                              |
|--------------------------|------------------------------|
| <b>1</b> Eyepiece socket | <b>2</b> Binocular section   |
| <b>3</b> Angle scale     | <b>4</b> Dovetail ring mount |

### 10.2.2 Binocular Photo Tube Pol 20°/23 (100:0/0:100)

**Purpose** Binocular photo tubes are used to visualize the microscopic image by means of the eyepieces. Using the camera port, the microscopic image can be sent to a connected camera.

**Position** The binocular photo tubes are mounted on the top of the stand.

**Function** The pupil distance and the viewing height can be adjusted with the aid of the binocular section. The following features and controls are available:

- upright image
- camera port with toggle graduation (100:0/0:100)
- viewing angle 20°
- field of view 23 mm



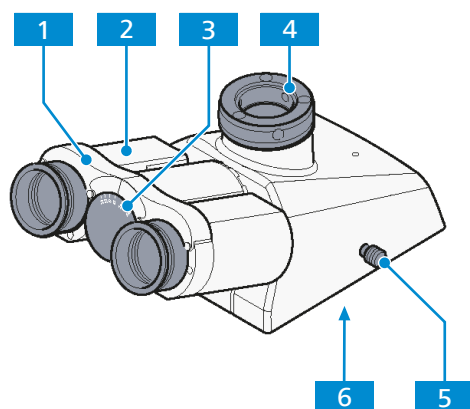


Fig. 34: Binocular Photo Tube Pol 20°/23 (100:0/0:100)

- |          |                                     |          |                     |
|----------|-------------------------------------|----------|---------------------|
| <b>1</b> | Eyepiece socket                     | <b>2</b> | Binocular section   |
| <b>3</b> | Angle scale                         | <b>4</b> | Camera port         |
| <b>5</b> | Slider for selecting the graduation | <b>6</b> | Dovetail ring mount |
- Slider pushed in: 100% light to eyepieces
  - Slider pulled out: 100% light to camera

### 10.2.3 Binocular Ergo Photo Tube 20°/23 (100:0/0:100)

**Purpose** Binocular photo tubes are used to visualize the microscopic image by means of the eyepieces. Using the camera port, the microscopic image can be sent to a connected camera.

**Position** The binocular photo tubes are mounted on the top of the stand.

**Function** The pupil distance and the viewing height can be adjusted with the aid of the binocular section.

The following features and controls are available:

- upright image
- camera port with toggle graduation (100:0/0:100)
- viewing angle 20°
- field of view 23 mm, usable 22 mm
- vertical adjustment of 44 mm with vertical scale

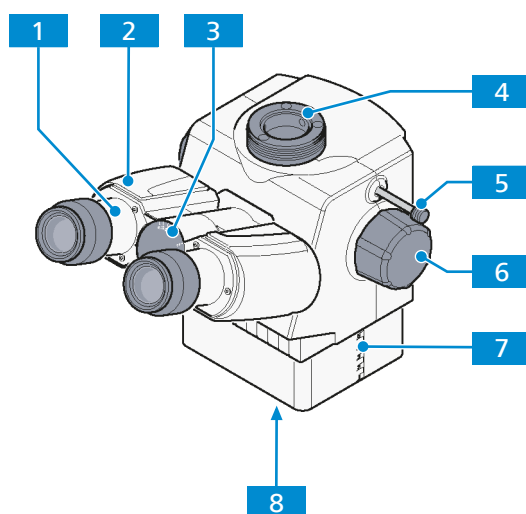


Fig. 35: Binocular Ergo Photo Tube 20°/23 (100:0/0:100)

- |  |   |
|--|---|
| <b>1</b> Eyepiece socket   | <b>2</b> Binocular section                                    |
| <b>3</b> Angle scale   | <b>4</b> Camera port  |
| <b>5</b> Slider for selecting the graduation <ul style="list-style-type: none"> <li>▪ Slider pushed in: 100% light to eye-pieces</li> <li>▪ Slider pulled out: 100% light to camera</li> </ul> | <b>6</b> Rotary knob for vertical adjustment (right and left) |
| <b>7</b> Vertical scale  | <b>8</b> Dovetail ring mount                                  |

#### 10.2.4 Binocular Photo Tube 30°/23 (98:2), Reversed Image

**Purpose** Binocular photo tubes provide users with the convenience of accessing the sample images using eyepieces as well as monitors through camera at the same time. This specific splitting ratio allow users to look at the sample with naked eyes even though light intensity is set to maximum level.

**Position** The binocular photo tubes are mounted on the top of the stand.

**Function** The pupil distance and viewing height can be adjusted by bending the binocular device up or down.

The following features and controls are available:

- Reverse image
- Camera port with splitting ratio 98:2
- Viewing angle 30°
- Field of view 23 mm

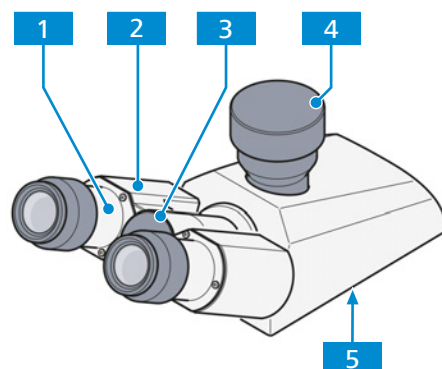


Fig. 36: Binocular photo tube 30°/23 (98:2), reversed image

- |                              |                            |
|------------------------------|----------------------------|
| <b>1</b> Eyepiece socket     | <b>2</b> Binocular section |
| <b>3</b> Angle scale         | <b>4</b> Camera port       |
| <b>5</b> Dovetail ring mount |                            |

### 10.2.5 Binocular Photo Tube 30°/23 (100:0/30:70/0:100), Reversed Image

**Purpose** Binocular photo tubes provide users with the convenience of accessing the sample images using eyepieces as well as monitors through camera at the same time. This specific splitting ratio allows users to freely switch between looking at the sample with naked eyes or using monitors or at simultaneously.

**Position** The binocular photo tubes are mounted on the top of the stand.

**Function** The pupil distance and viewing height can be adjusted by bending the binocular device up or down.

The following features and controls are available:

- Reverse image
- Camera port with splitting ratio 100:0/30:70/0:100
- Viewing angle 30°
- Field of view 23 mm

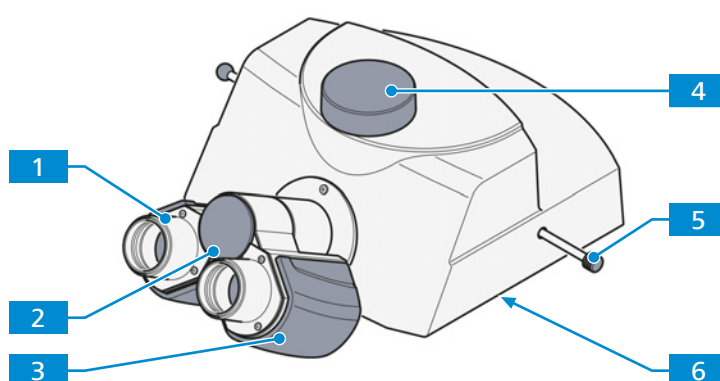



Fig. 37: Binocular photo tube 30°/23 (100:0/30:70/0:100), reversed image

- |   |                              |
|---|------------------------------|
| <b>1</b> Eyepiece socket                                      | <b>2</b> Angle scale         |
| <b>3</b> Binocular section                                    | <b>4</b> Camera port         |
| <b>5</b> Slider for selecting the graduation (right and left) | <b>6</b> Dovetail ring mount |
- Slider pushed in: 100% light to eyepieces
  - Slider pushed in to the middle: 30% light to eyepieces, 70% light to camera
  - Slider pulled out: 100% light to camera. 100% light to camera

## 10.3 Stages

A stage is a platform at right angles to the optical axis of the microscope, which carries the sample and which is often fitted with mechanical movements (as in a mechanical stage) to allow easy positioning of the object in the x- and y-axis, and movement along, and rotation about the z-axis.


### 10.3.1 Assembling a Mechanical Stage and Sample Holder

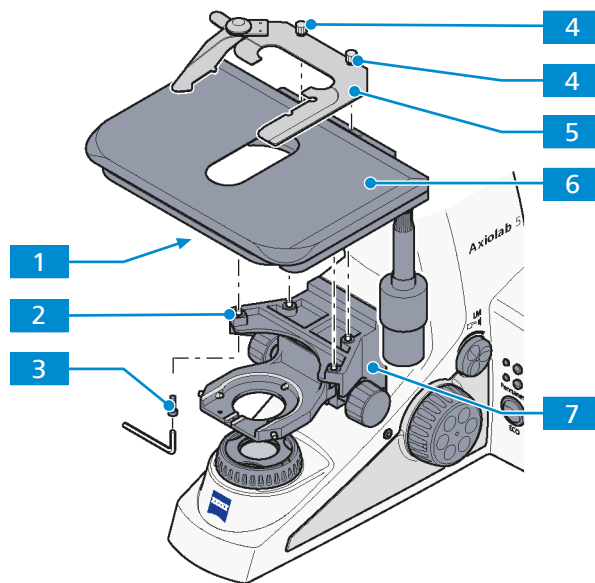

**Parts and Tools**    Hex key, 3.0 mm

- Procedure**
1. Place the stage **6** on the stage carrier **7** so that the threaded holes **1** on the bottom of the stage are positioned above the stage carrier openings **2** and hold the stage.
  2. Insert four fastening screws **3** through the stage carrier from below and screw them into the bottom of the stage.
  3. Turn the stage to orient it in an XY direction and tighten the fastening screws.
  4. Place the sample holder **5** on the stage and fasten the two clamping screws .

### 10.3.2 Removing a Mechanical Stage and Sample Holder

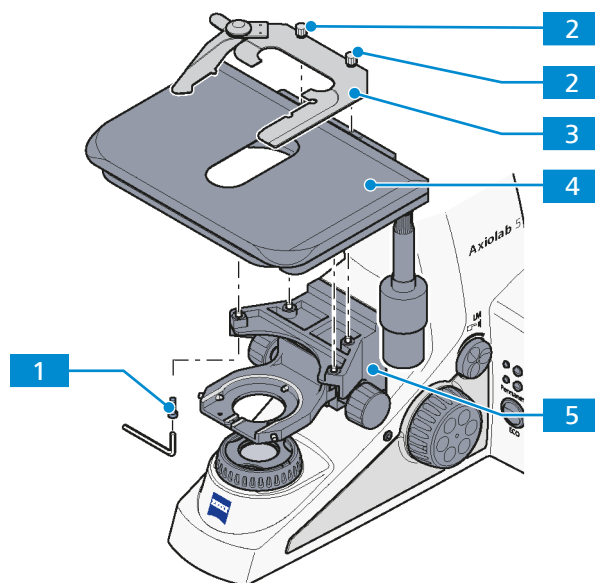

**Parts and Tools**  Hex key, 3.0 mm

- Procedure**
1. Loosen the two clamping screws **2** of the sample holder **3**.
  2. Remove the sample holder.
  3. Hold the stage **4** and remove the four fastening screws **1** on the stage carrier.
  4. Remove stage upwards from the stage carrier **5**.

### 10.3.3 Rackless Mechanical Stage, 75x30 R

**Purpose** Mechanical stages are used for fixing and positioning the sample for examination.

**Position** The mechanical stages are mounted on the stage carrier of the stand.

**Function** The sample is fixed on the stage by means of the sample holder. For this purpose, the sample holder is equipped with a spring lever.

The sample is positioned in the beam path by means of the two coaxial drives in X and Y direction. The adjustment range can be read off the respective vernier scale.

The following features and controls are available:

- rackless stage
- coaxial drives in X and Y adjustment on the right (R)
- travel range 75x30mm
- with hardcoat anodized surface

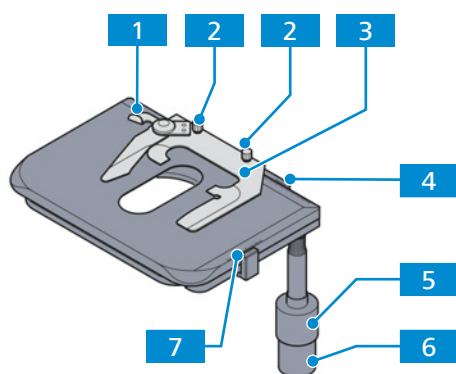


Fig. 38: Rackless mechanical stage, 75x30 R

- |   |   |
|---|---|
| <b>1</b> Spring lever   | <b>2</b> Knurled screw (2x) for fixing the sample holder to the stage |
| <b>3</b> Sample holder for one slide                            | <b>4</b> Vernier scale for display of the adjustment range in X       |
| <b>5</b> Coaxial knurled knob for Y adjustment                  | <b>6</b> Coaxial knurled knob for X adjustment                        |
| <b>7</b> Vernier scale for display of the adjustment range in Y |   |

### 10.3.4 Mechanical Stage, 75x30 R

**Purpose** Mechanical stages are used for fixing and positioning the sample for examination.

**Position** The mechanical stages are mounted on the stage carrier of the stand.

**Function** The sample is fixed on the stage by means of the sample holder. For this purpose, the sample holder is equipped with a spring lever.

The sample is positioned in the beam path by means of the two coaxial drives in X and Y direction. The adjustment range can be read off the respective vernier scale.

The following features and controls are available:

- stage with rack and pinion
- coaxial drives in X and Y adjustment on the right (R)
- travel range 75x30mm
- with hardcoat anodized surface

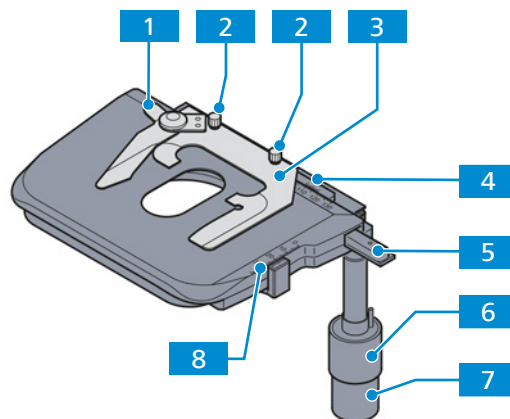


Fig. 39: Mechanical Stage, 75x30 R

- |   |  |
|---|--|
| <p><b>1</b> Spring lever</p> <p><b>3</b> Specimen holder for one slide</p> <p><b>5</b> Rack</p> <p><b>7</b> Coaxial knurled knob for X adjustment</p> | <p><b>2</b> Knurled screw (2x) for fixing the sample holder to the stage</p> <p><b>4</b> Vernier scale for display of the adjustment range in X</p> <p><b>6</b> Coaxial knurled knob for Y adjustment</p> <p><b>8</b> Vernier scale for display of the adjustment range in Y</p> |
|---|--|

### 10.3.5 Rotary Stage Pol 360° with Sample Guide

**Purpose** Rotary stages are used for fixing and positioning the sample for examination in polarized light.

**Position** The rotary stages are mounted on the stage carrier of the stand.

**Function** The sample is fixed on the stage by means of the sample guide. For this purpose, the sample guide is equipped with a spring lever.

The sample is positioned in the beam path by means of the two knurled knobs of the sample guide. The adjustment range can be read off the respective vernier scale.

The following features and controls are available:

- optionally equipped with: attachable sample guide for use of standard slides 45x25 mm and 75x25 mm (3"x1")
- 360° rotation with lock
- click stop every 45°; enabled or disabled via the control knob


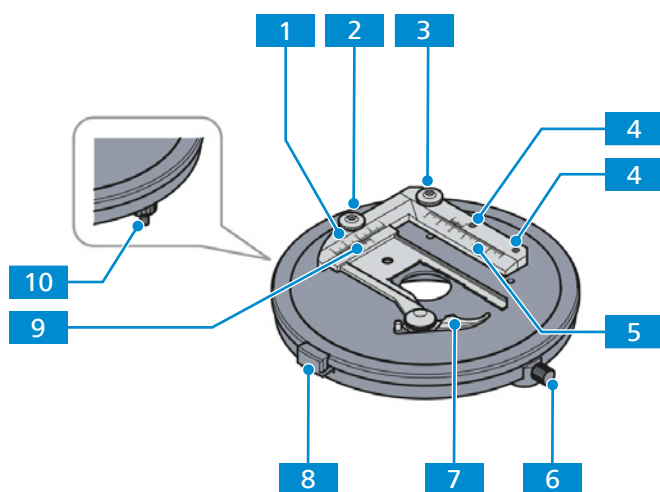


Fig. 40: Rotary Stage Pol 360° with sample Guide

- |   |   |
|---|---|
| <b>1</b> sample guide   | <b>2</b> Knurled knob for adjustment in Y direction                         |
| <b>3</b> Knurled knob for adjustment in X direction             | <b>4</b> Clamping screw (2x, AF 2) for fixing the sample guide to the stage |
| <b>5</b> Vernier scale for display of the adjustment range in X | <b>6</b> Knurled screw for locking the rotation, 360° rotation possible     |
| <b>7</b> Spring lever   | <b>8</b> Click stop every 45°   |
| <b>9</b> Vernier scale for display of the adjustment range in Y | <b>10</b> Control knob to enable/disable the clip stop function             |




### 10.3.5.1 Removing the Rotary Stage and Sample Guide or Clamping Device

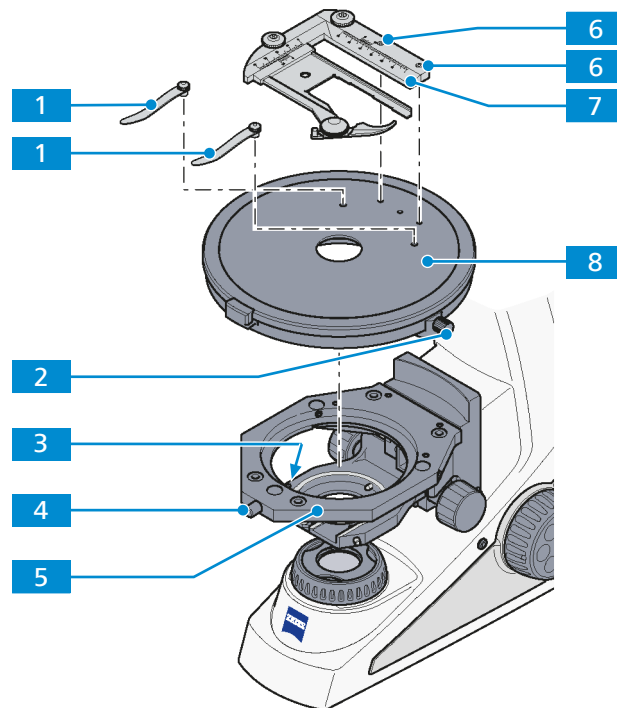

**Parts and Tools**  Hex key, 2.0 mm

- Procedure**
1. Loosen the screw cap **4** from the spring housing (about three rotations).
  2. Press the rotary stage **8** to the front against the spring-loaded pin **3**, lift it off the stage carrier **5** from the back and remove it upwards.
  3. Re-tighten the screw cap.
  4. If available, loosen the two clamping screws **6** on the sample guide **7**.
  5. Remove the sample guide by lifting it upwards.
  6. If available, remove the clamping device **1** from the rotary stage.

### 10.3.5.2 Assembling the Rotary Stage and Sample Guide or Clamping Device

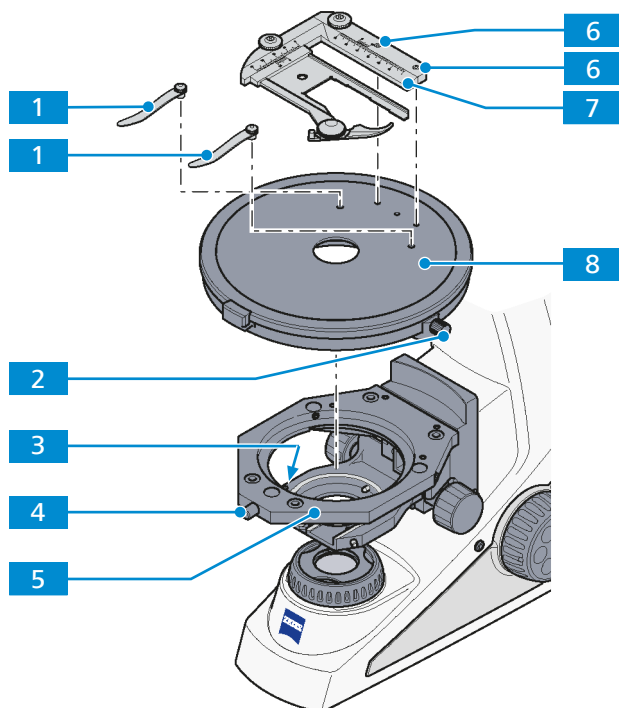


- Procedure**
1. Loosen the screw cap **4** of the spring housing with approx. three rotations.
  2. Place the rotary stage **8** with the groove of the dovetail (stage bottom) on the spring-loaded pin **3**.
  3. Attach the rotary stage with the knurled screw for locking the rotation **2** pointing to the front right.
  4. Press the rotary stage to the front against the spring-loaded pin and lower it towards the back into the stage carrier **5**, then release it.
  5. Re-tighten the screw cap.
  6. If required, insert the clamping device **1** into the holes provided.
  7. If required, insert the sample guide **7** with the two cylindrical pins on the underside into the holes provided and tighten the two clamping screws **6**. Use a 2 mm hex key.

### 10.3.5.3 Centering the Rotary Stage

With high-power objectives, centering can be exact only for one selected objective.

All stages are factory-precentered, i.e. while rotating the stage the sample detail set to the center will remain in the center. If the sample detail moves out of the center of the field of view while the stage is being rotated, the stage should be re-centered.

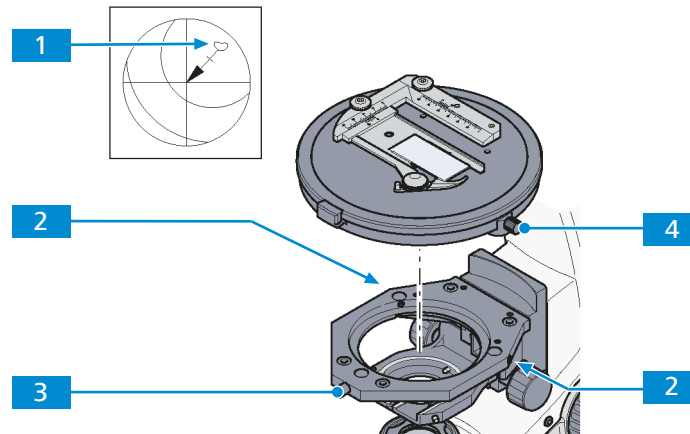

**Parts and Tools** Hex key, 1.5 mm


- Prerequisite**
- ✓ The rotary stage is installed in the stage carrier.
  - ✓ A high-contrast sample and an eyepiece with a crossline reticle are available.
  - ✓ The microscope is adjusted for *transmitted light brightfield microscopy* [▶ 68].

- Procedure**
1. Turn the nosepiece to swing the non-centering objective mount.
  2. Loosen the knurled screw for locking the rotation **4** and the screw cap **3** on the stage carrier.
  3. Rotate the stage to determine the position of maximum offset of the sample feature **1** (origin of arrow) from the center of the eyepiece reticle.
  4. Reset the two centering screws **2** on the stage carrier to move the sample detail by half an arrow length in the direction of the crossline center.
  5. Check whether the sample detail moves when the stage is rotated again; repeat the procedure, if required.
  6. Re-tighten the screw cap.

#### 10.3.5.4 Centering the Objectives of the Polarization Stand

**Purpose** Stage centering of the non-centering objective mount is necessary to ensure that a sample feature located in the center of the field of view does not drift out while rotating the stage. By centering of the remaining objectives, the sample feature remains in the center of the field of view even after changing the objective.

The nosepiece 5x Pol is equipped with one fixed and four centerable objective positions.

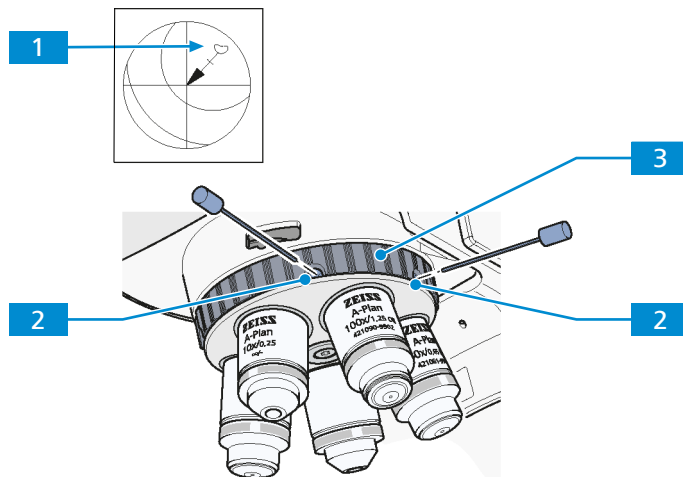

**Parts and Tools** 🔑 2x Hex key, 1.5 mm

- Prerequisite**
- ✓ The microscope is adjusted for *transmitted light brightfield microscopy* [▶ 68].
  - ✓ The *rotary stage* [▶ 121] is centered.
  - ✓ A high-contrast sample and an eyepiece with crossline reticle are available.

- Procedure**
1. Turn the nosepiece using the knurled ring **3** of the nosepiece to swing the non-centering objective position.
  2. Check if the rotary stage is centered to this objective.
  3. Turn the nosepiece to move a centering objective mount into the beam path.
  4. Rotate the stage to determine the position of maximum offset of the sample detail **1** (origin of the arrow) from the center of the eyepiece reticle.
  5. Reset the two centering screws **2** of the nosepiece position to move the sample detail by half the arrow length in the direction of the crossline center.
  6. Check whether the sample detail moves when the stage is rotated again; repeat the procedure if required.
  7. Center the other three objectives in the same manner.

## 10.4 Analyzer Sliders

### 10.4.1 Analyzer Slider TL/RL, Fixed

**Purpose** The analyzer slider is used to set the polarization contrast technique.

**Position** The analyzer slider is inserted into the 12x46 slot of the intermediate plate for analyzer slider 12x46 mounted between stand and tube.

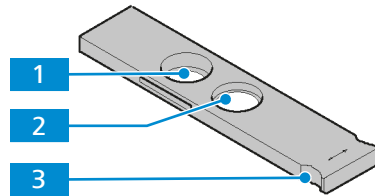


Fig. 41: Analyzer slider TL/RL, fixed

- |          |                |          |          |
|----------|----------------|----------|----------|
| <b>1</b> | Empty position | <b>2</b> | Analyzer |
| <b>3</b> | Handle         |          |          |

### 10.4.2 Analyzer Slider TL/RL, with Lambda Plate, 360° Rotatable

**Purpose** The analyzer slider is used to set the polarization contrast technique.

**Position** The analyzer slider is inserted into the 12x46 slot of the intermediate plate for analyzer slider 12x46 mounted between stand and tube.

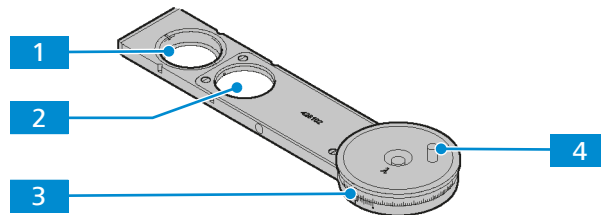


Fig. 42: Analyzer slider TL/RL, with lambda plate, 360° rotatable

- |          |                |          |                                   |
|----------|----------------|----------|-----------------------------------|
| <b>1</b> | Empty position | <b>2</b> | Analyzer and lambda plate         |
| <b>3</b> | Angle scale    | <b>4</b> | Handle to rotate the lambda plate |

### 10.4.3 Analyzer Slider TL/RL with Lambda Plate, each Rotatable +/- 10°

**Purpose** The analyzer slider is used to set the polarization contrast technique.

**Position** The analyzer slider is inserted into the 12x46 slot of the intermediate plate for analyzer slider 12x46 mounted between stand and tube.

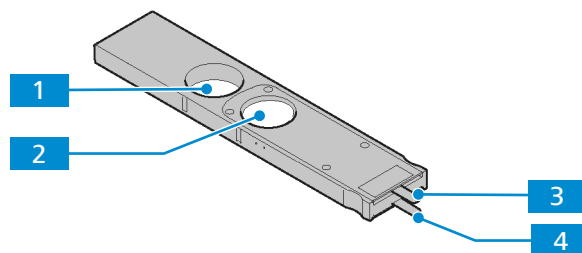


Fig. 43: Analyzer slider TL/RL with lambda plate, each rotatable +/- 10°

- |          |                                       |          |                                   |
|----------|---------------------------------------|----------|-----------------------------------|
| <b>1</b> | Empty position                        | <b>2</b> | Analyzer and lambda plate         |
| <b>3</b> | Handle for adjusting the lambda plate | <b>4</b> | Handle for adjusting the analyzer |

## 10.5 Polarizers

### 10.5.1 Polarizer D, Fixed, Removable

**Purpose** The polarizer for transmitted light is used to polarize the light of the transmitted light source. The polarizer can be swivelled into or out of the beam path using the handle.

**Position** The polarizer is mounted on the bottom of the condenser carrier.

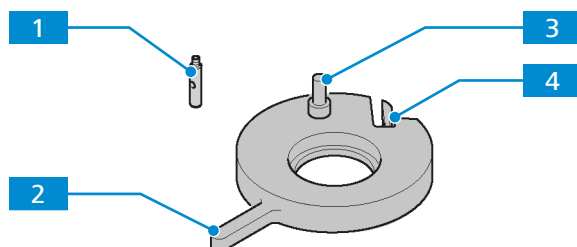


Fig. 44: Polarizer D, fixed, removable

- |                      |  |
|----------------------|--|
| <b>1</b> Locking pin | <b>2</b> Handle of the polarizer for swivelling in/out |
| <b>3</b> Holding pin | <b>4</b> Locking clamp                                 |

### 10.5.2 Polarizer D, 90°, Rotatable, Removable

**Purpose** The polarizer for transmitted light is used to polarize the light of the transmitted light source. The polarizer can be swivelled into or out of the beam path using the handle.

**Position** The polarizer is mounted on the bottom of the condenser carrier.

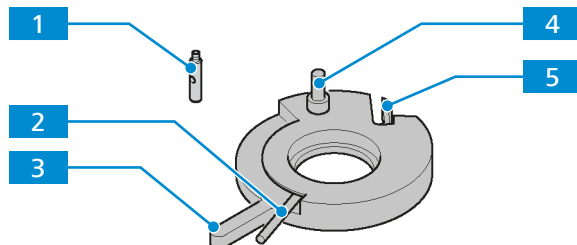


Fig. 45: Polarizer D, 90°, rotatable, removable

- |  |   |
|--|---|
| <b>1</b> Locking pin                                   | <b>2</b> Lever for rotating the polarizer |
| <b>3</b> Handle of the polarizer for swivelling in/out | <b>4</b> Holding pin                      |
| <b>5</b> Locking clamp                                 |   |

### 10.5.3 Polarizer, Fixed, with Lambda Plate, Rotatable

**Purpose** The polarizer for transmitted light is used to polarize the light of the transmitted light source. The polarizer can be swivelled into or out of the beam path using the handle.

**Position** The polarizer is mounted on the bottom of the condenser carrier.  
Polarizer and lambda plate can be swivelled in/out separately.

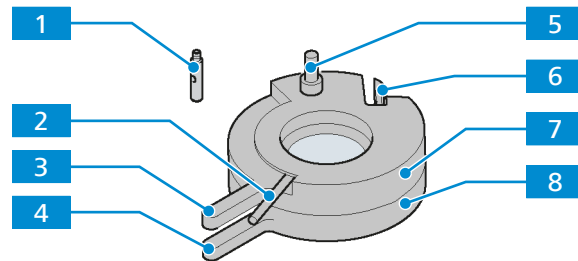


Fig. 46: Polarizer, fixed, with lambda plate, rotatable

- |   |  |
|---|--|
| <b>1</b> Locking pin                                      | <b>2</b> Lever for rotating the lambda plate           |
| <b>3</b> Handle of the lambda plate for swivelling in/out | <b>4</b> Handle of the polarizer for swivelling in/out |
| <b>5</b> Holding pin                                      | <b>6</b> Locking clamp                                 |
| <b>7</b> Lambda plate, rotatable by 90°                   | <b>8</b> Polarizer                                     |

### 10.5.4 Polarizer, Rotatable, with Color Filter Carrier

**Purpose** The polarizer for transmitted light is used to polarize the light of the transmitted light source. With the help of the color filter carrier optical filter elements can be placed in the beam path. The polarizer and the filter carrier can be swivelled into or out of the beam path using the handle.

**Position** The polarizer with filter carrier is mounted on the bottom of the condenser carrier.  
Polarizer and filter carrier can be swivelled in/out separately.

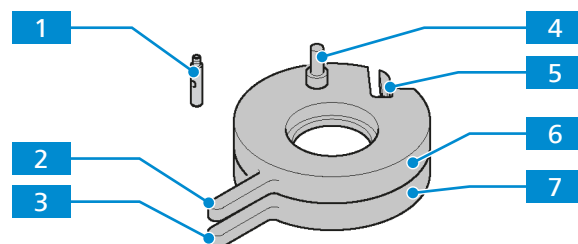


Fig. 47: Polarizer, rotatable, with color filter carrier

- |   |  |
|---|--|
| <b>1</b> Locking pin                        | <b>2</b> Handle of the polarizer for swivelling in/out |
| <b>3</b> Handle of the color filter carrier | <b>4</b> Holding pin                                   |
| <b>5</b> Locking clamp                      | <b>6</b> Polarizer                                     |
| <b>7</b> Color filter carrier               |  |



### 10.5.5 Circular Polarizer D

**Purpose** The polarizer for transmitted light is used to polarize the light of the transmitted light source. The polarizer can be swivelled into or out of the beam path using the handle.

**Position** The polarizer is mounted on the bottom of the condenser carrier.

Upper part and lower part of the polarizer can be swivelled in/out separately.

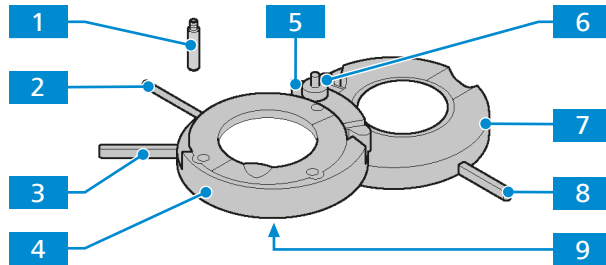


Fig. 48: Circular polarizer D

- |  |   |
|--|---|
| <b>1</b> Locking pin   | <b>2</b> Lever for rotating the lambda 4/plate, 90° possible                      |
| <b>3</b> Handle of the upper part of the polarizer for swivelling in/out | <b>4</b> Lambda/4 plate in the upper part of the circular polarizer               |
| <b>5</b> Locking clamp   | <b>6</b> Holding pin  |
| <b>7</b> Lower part of the circular polarizer                            | <b>8</b> Handle of the lower part of the circular polarizer for swivelling in/out |
| <b>9</b> Adjustment slit (2x)  |   |

### 10.5.6 Low-power System for Objectives 2.5x/4x

**Purpose** The low-power system is for full display field illumination when using an objective with a weak magnification factor (2.5x–4x) in combination with the Condenser 0.9/1.25 H.

**Position** The low-power system is mounted behind the condenser carrier.

**Function** It can be centered and remains swivelled into the beam path for as long as the respective objective is in use.

The illumination of weak objective magnifications can be centered with the centering screws. For this purpose, the condenser should be centered on the other objectives without the low-power system.

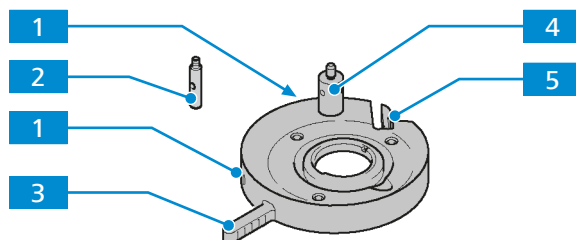



Fig. 49: Low-power system for objectives 2.5x/4x

- |   |                      |
|---|----------------------|
| <b>1</b> Centering screw (2x)                                 | <b>2</b> Locking pin |
| <b>3</b> Handle of the low-power system for swivelling in/out | <b>4</b> Holding pin |
| <b>5</b> Locking clamp  |                      |

### 10.5.6.1 Assembling and Centering the Low-power System

#### Info

The low-power system can be used only in combination with the condenser 0.9/1.25.

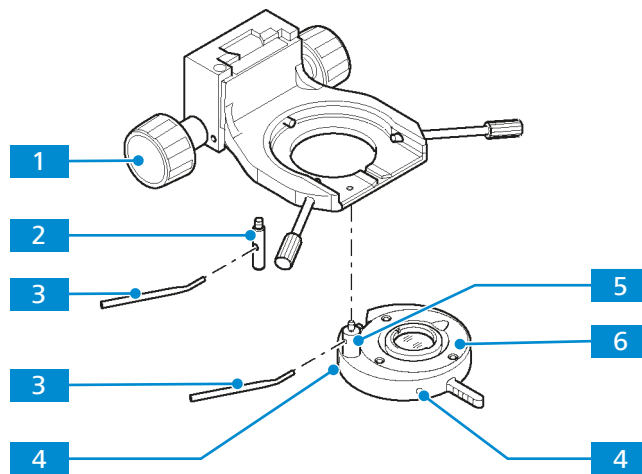

**Parts and Tools** 2x Hex key, 1.5 mm

#### Procedure

1. Lift the condenser carrier together with its knurled knob **1** upwards as far as it will go.
2. Hold the low-power system **6** parallel to the underside of the condenser carrier and screw the holding pin **5** of the low-power system with the angled adjusting lever **3** into the front left threaded hole below the condenser carrier as far as it will go.
3. Screw the locking pin **2** with the adjusting lever as far as it will go into the rear threaded hole of the condenser carrier.
4. Swivel the low-power system into the beam path using the handle.  
→ Make sure that it is securely engaged.
5. Switch on the microscope.
6. Set transmitted light illumination.
7. Open the aperture diaphragm and luminous-field diaphragm completely.
8. Adjust both adjustment screws **4** until the field of vision is optimally illuminated.

### 10.5.7 Color Filter Carrier 3x for Filter d=32mm

**Purpose** With the help of the color filter carrier optical filter elements can be placed in the beam path. The filter carriers can be swivelled into or out of the beam path using the handle.

**Position** The color filter carrier is mounted on the bottom of the condenser carrier. The three filter carriers can be swivelled in/out separately.

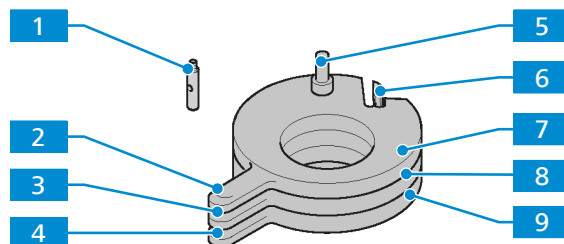



Fig. 50: Color filter carrier 3x for filter d=32mm

- |  |   |
|--|---|
| <b>1</b> Locking pin   | <b>2</b> Handle of the first filter carrier for swivelling in/out |
| <b>3</b> Handle of the second filter carrier for swivelling in/out | <b>4</b> Handle of the third filter carrier for swivelling in/out |
| <b>5</b> Holding pin   | <b>6</b> Locking clamp  |
| <b>7</b> First filter carrier                                      | <b>8</b> Second filter carrier                                    |
| <b>9</b> Third filter carrier                                      |   |

### 10.5.8 Assembling Polarizer or Color Filter Carrier onto the Condenser Carrier

The following polarizers or the filter carrier can be installed on the condenser carrier:

- Polarizer D, fixed, removable
- Polarizer D, 90° rotary, removable
- Polarizer, rotary, with color filter carrier
- Polarizer, fixed, with lambda plate, rotary
- Circular polarizer D, fixed, with rotary lambda/4 plate
- Color filter carrier 3x for filter d=32 mm

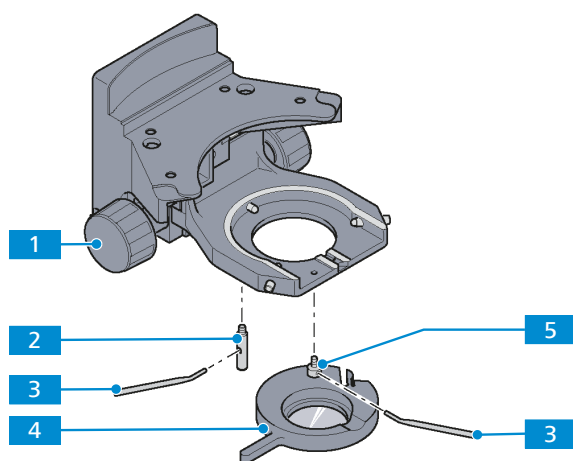

- Procedure**
1. Lift the condenser carrier upwards as far as it will go using the knurled knob **1**.
  2. Hold the polarizer **4** or filter carrier parallel to the underside of the condenser carrier and screw the holding pin **5** of the polarizer with the angled adjusting lever **3** into the front left threaded hole below the condenser carrier as far as it will go.
  3. Screw the locking pin **2** with the adjusting lever as far as it will go into the rear threaded hole of the condenser carrier.

Proceed in the reverse order for removal.

## 10.6 DIC Slider C 6x20

**Purpose** The DIC slider is used to set the DIC contrast technique.

**Position** The DIC slider is inserted into the 6x20 slot above the nosepiece.

The DIC slider is available in two versions:

- DIC Slider C 6x20 for objectives EC Epiplan 5x - 20x
- DIC Slider C 6x20 for objectives EC Epiplan 50x - 100x

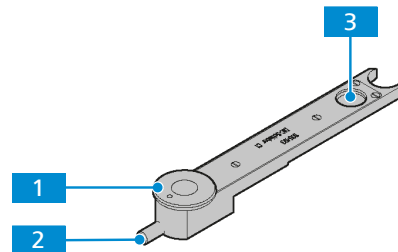


Fig. 51: DIC Slider C 6x20

- |                        |                        |
|------------------------|------------------------|
| <b>1</b> Setting wheel | <b>2</b> Setting screw |
| <b>3</b> DIC prism     |                        |

## 10.7 Condenser, achromatic-aplanatic 0.9 BF DF PhC DIC

**Purpose** Condensers are used to optimize the transmitted light illumination. The condenser is usable for brightfield, darkfield, phase contrast and DIC applications.

**Position** The condenser is mounted on the condenser carrier of the stand.


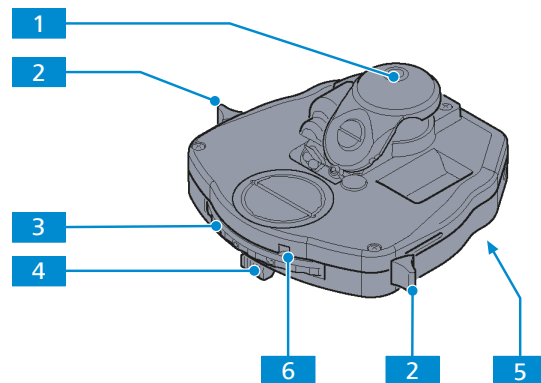


Fig. 52: Condenser, achromatic-aplanatic 0.9 BF DF PhC DIC


- |   |   |
|---|---|
| <b>1</b> Front lens   | <b>2</b> Lever for switching the front lens in/out (left/right) |
| <b>3</b> Knurled ring for adjusting the position of the 5-positions turret disk | <b>4</b> Sliding control for setting the aperture diaphragm     |
| <b>5</b> Dovetail ring mount  | <b>6</b> Display field of the adjusted turret disk position     |

### 10.7.1 Assembling the Condenser, achromatic-aplanatic 0.9 BF DF PhC DIC

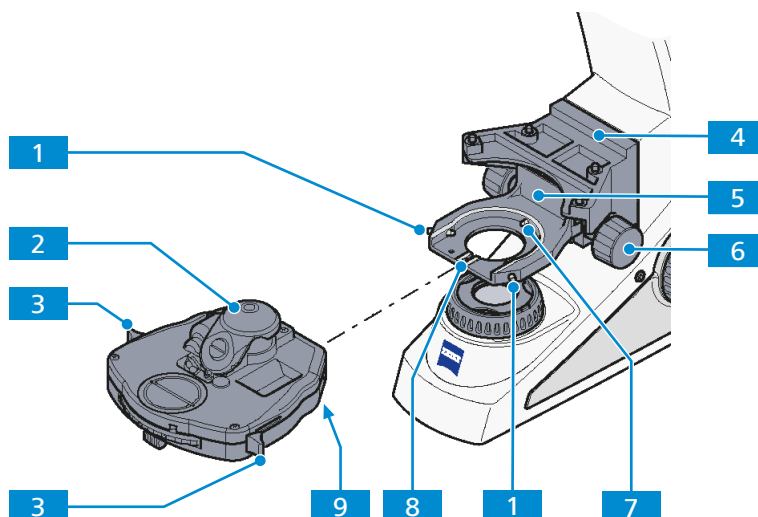

- Procedure**
1. Carefully move the stage carrier **4** with the focusing drive to the upper stop position.

**NOTICE** Make sure that the stage does not collide with the objective.

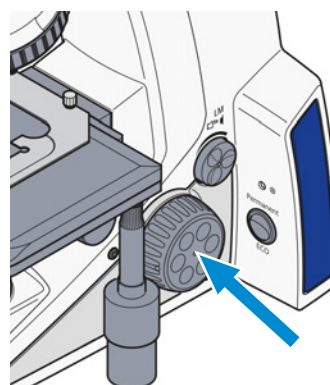


2. Swivel out the front lens **2** on the condenser using the lever **3**.
3. Unscrew both centering screws **1** on the condenser carrier **5** until their ends are no longer visible.
4. Using the knurled knob **6** for vertical adjustment, push the condenser carrier down as far as it will go.  
**NOTICE** If using a low-power system, make sure that this does not come to rest on the luminous-field diaphragm.
5. Insert the condenser between the condenser carrier and the stage carrier **4**. In doing so, align the screwed stud bolt **9** on the underside of the condenser with the groove **8** of the condenser carrier.
6. Press the condenser with the dovetail ring against the mainspring **7** of the condenser carrier until the condenser sits horizontally on the condenser carrier.
7. Screw in the centering screws **1** until they engage with the dovetail ring of the condenser.

### 10.7.2 Centering the Darkfield Diaphragm of the Condenser

**Parts and Tools**  2x Hex key, 1.5 mm

- Prerequisite**
- ✓ A suitable condenser with modulator disk is installed.
  - ✓ The illumination is adjusted for transmitted light brightfield microscopy.

- Procedure**
1. Set the modulator disks to position D (or DF = darkfield).
  2. Remove one eyepiece from the binocular tube or replace it with the auxiliary microscope.
  3. Observe the exit pupil of the objective.
  4. Turn the two centering screws, until the exit pupil of the objective appears homogeneously dark.

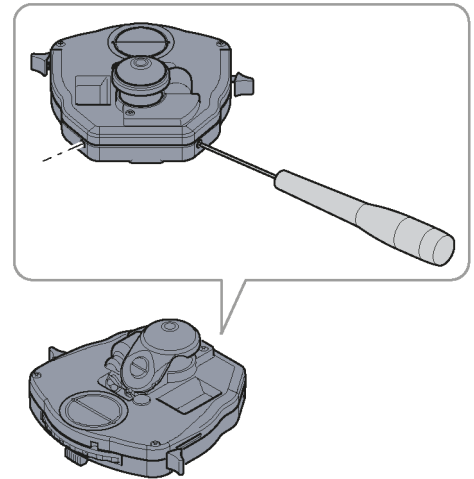


5. Insert the eyepiece.

### 10.7.3 Centering the Annular Phase Diaphragm of the Condenser


**Parts and Tools**  2x Hex key, 1.5 mm

- Prerequisite**
- ✓ A suitable condenser with modulator disk is installed.
  - ✓ The illumination is adjusted for transmitted light brightfield microscopy.

- Procedure**
1. Set the modulator disks to position **Ph** (phase contrast).
  2. Remove one eyepiece from the binocular tube or replace it with the auxiliary microscope.
  3. Observe the exit pupil of the objective.
  4. Check the centering and the overlap of the lighter annular phase diaphragm (in the condenser) with the darker phase ring (in the objective). Both rings must be centered and overlapping **B**.

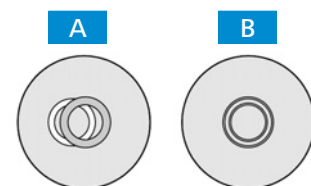

5. If the overlap is not exact **A**, recenter the lighter annular diaphragm.

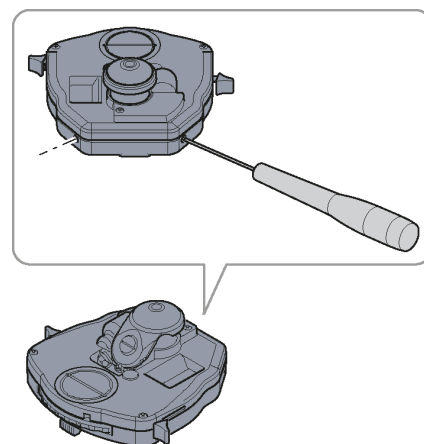

6. Remove the auxiliary microscope and replace the eyepiece.

## 10.8 Light Source for Transmitted Light

### **⚠ DANGER**

#### **Electric injury due to live parts**

When the microscope and its components are still switched on, coming in contact with live parts can lead to electric shock or burn.

- ▶ Switch off the microscope and its components prior to opening or cleaning.
- ▶ Disconnect live parts from the power supply.


### **⚠ CAUTION**

#### **Burning hazard due to hot light sources**

Light sources can become hot during processing.

- ▶ Allow it to cool down at least 15 min.

### 10.8.1 Removing the Cover (Transmitted Light Source)

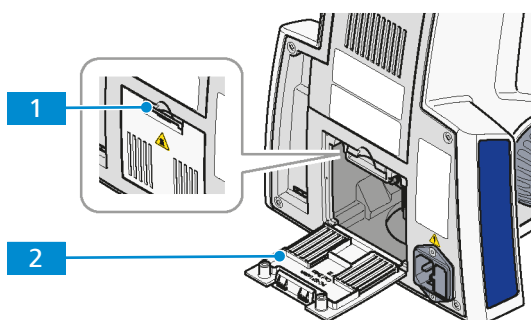


**Prerequisite** ✓ The microscope is switched off and disconnected from the mains.

- Procedure**
1. Press the clamping jaw **1** on the cover downwards.
  2. Swing the cover **2** down.
  3. Remove it from the retaining channels on the stand.
  4. Set the cover aside.

Proceed in the reverse order for installing.

### 10.8.2 Exchanging the Halogen Lamp

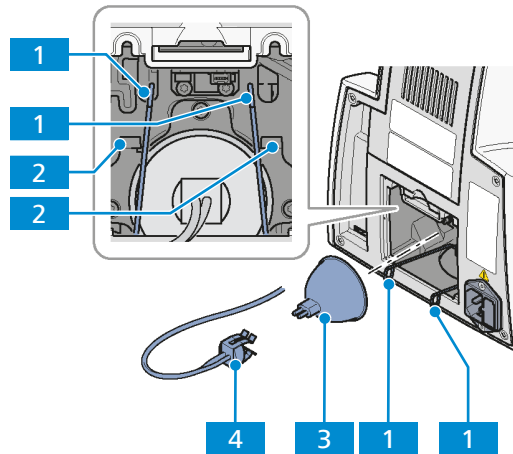

**Prerequisite** ✓ The cover is removed [▶ 132].

- Procedure**
1. Remove the lamp plug **4** from the halogen lamp **3**.
  2. Press the loops **1** on the securing clip of the lamp holder together and swing them out to the front.
  3. Remove the old halogen lamp.
  4. Position the new lamp with the lower front edge between the contact surface and securing clip.
  5. Lift the securing clip of the lamp holder with the lamp until they are completely enclosed in the lamp holder.
    - In doing so, slightly press the ends of the securing clip together and guide past the two upper retaining elements **2**.
    - Release the pressure until the clamping bracket opens and engages on both sides in the retaining elements.
  6. Check that the lamp is seated correctly.
  7. Push the lamp plug onto the pins of the lamp.
  8. Ensure that it engages properly to avoid bending the pins.
  9. Insert the cable of the lamp plug into the stand so that it is not damaged when the cover is attached.
  10. Insert the cover [▶ 132].



### 10.8.3 Exchanging the LED Light Source

Exchanging the LED light source is done in the following steps:

- removing the cover [▶ 132]
- removing the LED light source including adapter [▶ 135]
- exchanging the LED light source in the adapter [▶ 135]
- installing the LED light source including adapter [▶ 135]
- installing the cover [▶ 132]


#### 10.8.3.1 Removing the LED Light Source

##### CAUTION

##### Eye damage or skin irritation due to hazardous light emission

The light source belongs to Risk Group 2 as specified in IEC 62471 and emits LED radiation and UV radiation. Eye damage or skin irritation may result from exposure.


- ▶ Never look directly into the light-emitting aperture of the light source.
- ▶ Avoid exposure of skin to the radiation. Use suitable protective equipment/protective clothing if required.
- ▶ Before installing or removing the light source always make sure it is switched off.

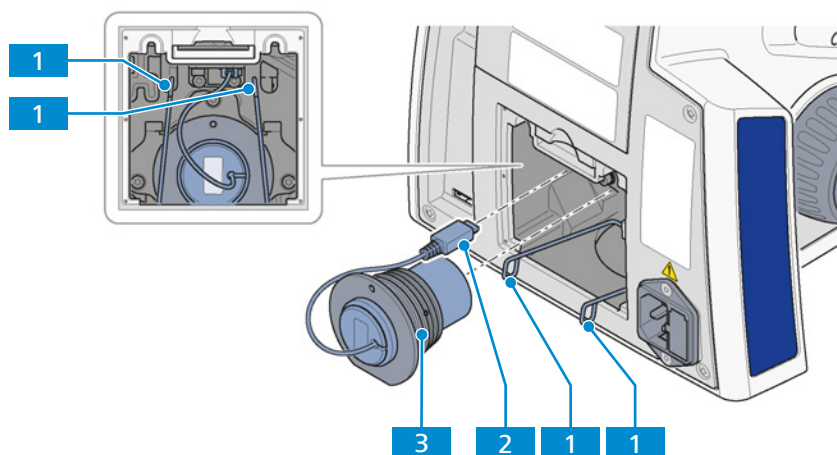

**Prerequisite** ✓ The cover is removed [▶ 132].

- Procedure**
1. Remove the LED light source plug **2** from the connector of the stand.
  2. Press the loops **1** on the securing clip of the lamp holder together and swing them out to the front.
  3. Pull the old LED light source with adapter **3** out of the holding tube.


### 10.8.3.2 Exchanging the LED Light Source in the Adapter

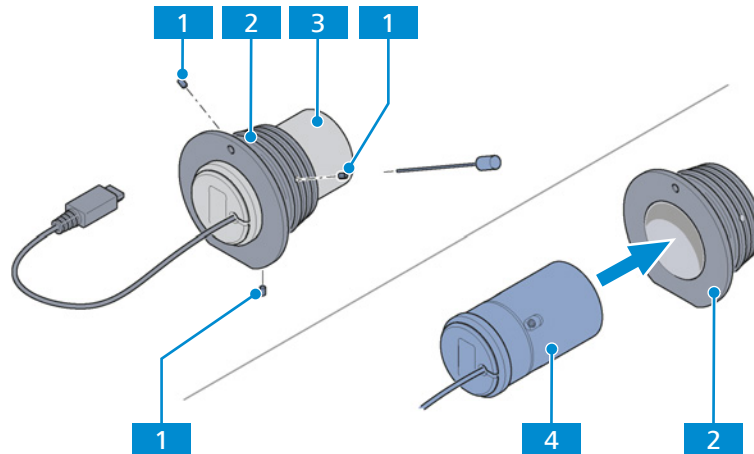

**Parts and Tools** Hex key, 2.5 mm

- Procedure**
1. Loosen the three side screws **1**.
  2. Remove the old LED light source **3** from the adapter **2**.
  3. Insert the new LED light source **4** into the adapter.
  4. Tighten the three side screws on the adapter.

### 10.8.3.3 Assembling the LED Light Source

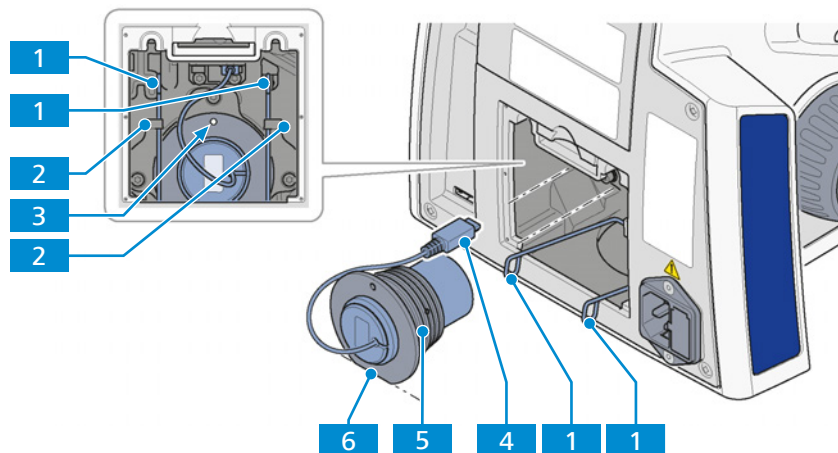

- Procedure**
1. Insert the new LED light source with the adapter **5** into the holding tube up to the stop.
  2. Position the new LED light source with the pinhole **3** at the top, or with the adapter edge **6** aligned between the contact surface and securing clips.
  3. Plug the light source plug **4** into the connector of the stand.
  4. Lift the securing clips of the lamp holder until they are completely enclosed in the lamp holder.
    - In doing so, slightly press the loops of the securing clips **1** together and guide past the two upper retaining elements **2**.
    - Release the pressure until the clamping bracket opens and engages on both sides in the retaining elements.
  5. *Insert the cover [▶ 132].*

## 10.9 Light Source for Reflected Light

### **DANGER**

#### **Electric injury due to live parts**

When the microscope and its components are still switched on, coming in contact with live parts can lead to electric shock or burn.

- Switch off the microscope and its components prior to opening or cleaning.
- Disconnect live parts from the power supply.


### **CAUTION**

#### **Burning hazard due to hot light sources**

Light sources can become hot during processing.

- Allow it to cool down at least 15 min.

### 10.9.1 Removing the Cover (Reflected Light Source)

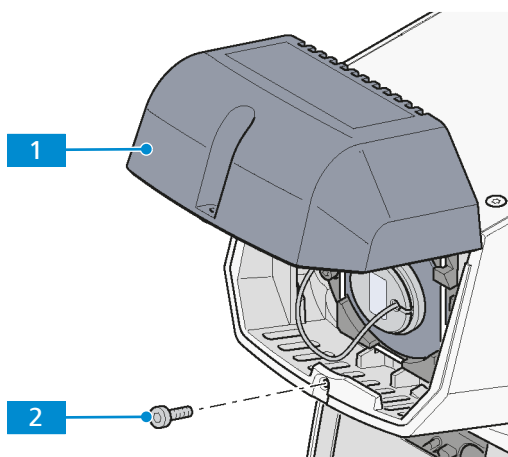


**Parts and Tools**  Hex key, 3.0 mm

**Prerequisite**  The microscope is switched off and disconnected from the mains.

- Procedure**
1. Unscrew the clamping screw **2** in the cover **1**.
  2. Swing the cover slightly upwards and apply pressure from below to remove it from the stand.

Proceed in the reverse order for installing.

### 10.9.2 Exchanging the Halogen Lamp

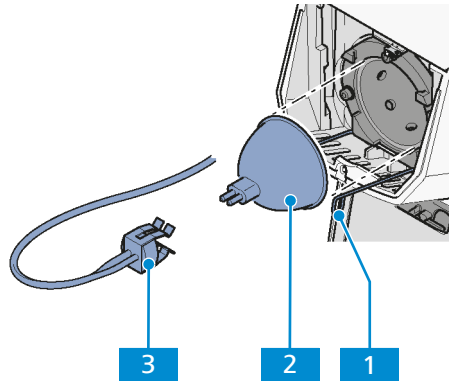

**Prerequisite** ✓ The cover is removed [▶ 136].

- Procedure**
1. Remove the lamp plug **3** from the halogen lamp **2**.
  2. Press the loops on the securing clip **1** of the lamp holder together and swing them out to the front.
  3. Remove the old halogen lamp.
  4. Position the new halogen lamp on the contact surface of the lamp holder (the lamp will be held securely by the groove).
  5. Press the securing clips on both sides of the lamp holder and swing them upwards until the securing clips are seated against the halogen lamp.
  6. Slowly release the securing clips so that they open and engage in the retaining elements to the right and left.
  7. Check that the halogen lamp is correctly seated.
  8. Push the lamp plug onto the pins of the lamp. Ensure that it engages properly to avoid bending the pins.
  9. Insert the cable of the lamp plug into the stand so that it is not damaged when the cover is attached.
  10. Insert the cover [▶ 136].

### 10.9.3 Exchanging the LED Light Source

exchanging the LED light source is done in the following steps:


- removing the cover [▶ 136]
- removing the LED light source including adapter [▶ 138]
- exchanging the LED light source in the adapter [▶ 139]
- assembling the LED light source including adapter [▶ 139]
- assembling the cover [▶ 136]

## 10.9.3.1 Removing the LED Light Source

**⚠ CAUTION****Eye damage or skin irritation due to hazardous light emission**

The light source belongs to Risk Group 2 as specified in IEC 62471 and emits LED radiation and UV radiation. Eye damage or skin irritation may result from exposure.


- ▶ Never look directly into the light-emitting aperture of the light source.
- ▶ Avoid exposure of skin to the radiation. Use suitable protective equipment/protective clothing if required.
- ▶ Before installing or removing the light source always make sure it is switched off.

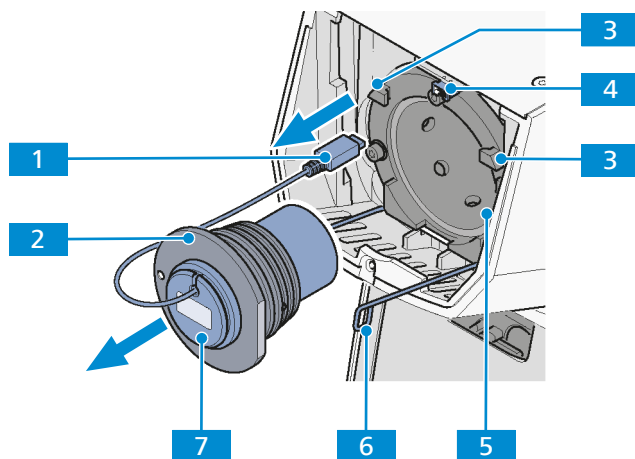

**Prerequisite** ✓ The cover is removed [▶ 136].

- Procedure**
1. Press the loops on the securing clip **6** of the lamp holder together and swing them out to the front.
  2. Remove the LED light source plug **1** from the connector **4** and the retaining elements **3** of the stand.
  3. Pull the old LED light source **7** with adapter **2** out of the holding tube **5**.


### 10.9.3.2 Exchanging the LED Light Source in the Adapter

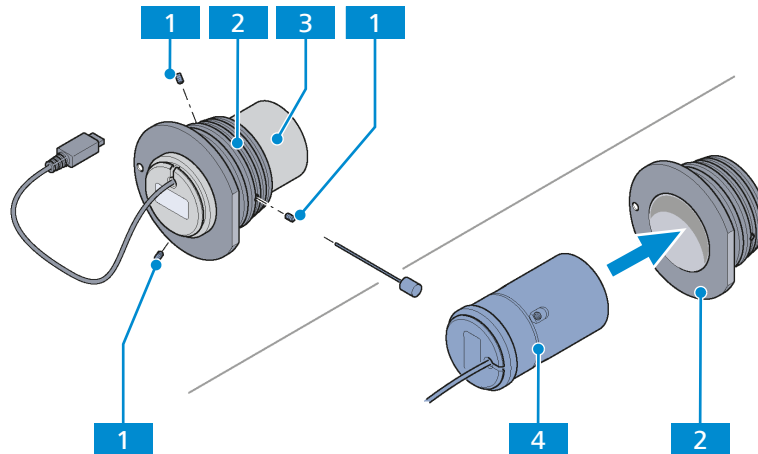

**Parts and Tools** Hex key, 2.5 mm

- Procedure**
1. Loosen the three side screws **1**.
  2. Remove the old LED light source **3** from the adapter **2**.
  3. Insert the new LED light source **4** into the adapter.
  4. Tighten the three side screws on the adapter.

### 10.9.3.3 Assembling the LED Light Source

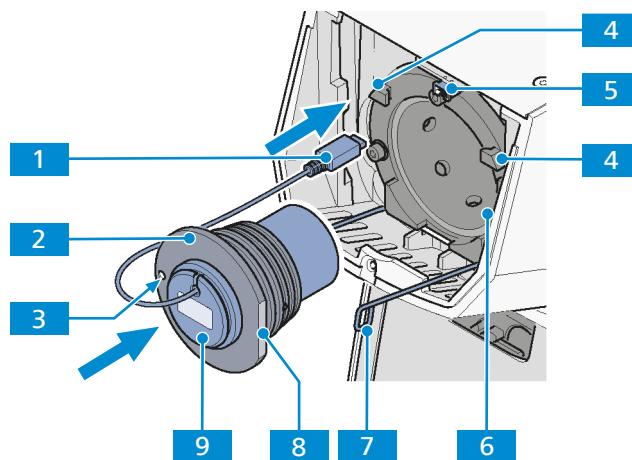

- Procedure**
1. Insert the new LED light source **9** with the adapter **2** into the holding tube **6** up to the stop.
  2. Position the new LED light source with the pinhole **3** at the left, or with the adapter edge **8** aligned to the right.
  3. Plug the light source plug **1** into the connector of the stand **5**.
  4. Lift the securing clips of the lamp holder until they are completely enclosed in the lamp holder.
    - In doing so, slightly press the loops of the securing clips **7** together and guide past the two retaining elements **4**.
    - Release the pressure until the clamping bracket opens and engages on both sides in the retaining elements.
  5. *Insert the cover [▶ 136].*

### 10.9.4 Exchanging the LED Modules of the Fluorescence Light Source


Exchanging the LED modules is done in the following steps:

- *removing the cover* [▶ 136]
- *removing the module holder* [▶ 140]
- *exchanging the LED modules* [▶ 140]
- *assembling the module holder* [▶ 141]
- *assembling the cover* [▶ 136]

#### 10.9.4.1 Removing the Module Holder

**Prerequisite** ✓ The cover is *removed* [▶ 136].

**Procedure** 1. Unscrew the module holder from the stand socket.

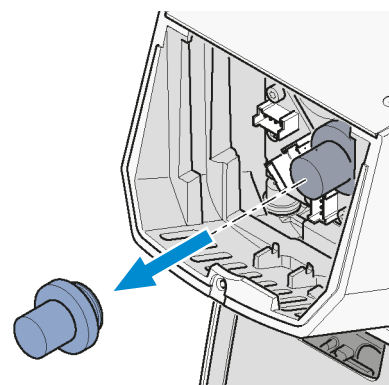


#### 10.9.4.2 Exchanging the LED Modules

##### **WARNING**

##### **Skin or eye injury due to hazardous light emission**

The light source belongs to Risk Group 3 as specified in IEC 62471 and emits LED radiation and UV radiation. Skin or eye injury can result from the exposure.

- ▶ Avoid any eye and skin exposure to the light-emitting aperture of the light source.
- ▶ Avoid exposure of skin to the radiation. Use suitable protective equipment/protective clothing if required.
- ▶ Before installing or removing the light source always make sure it is switched off.

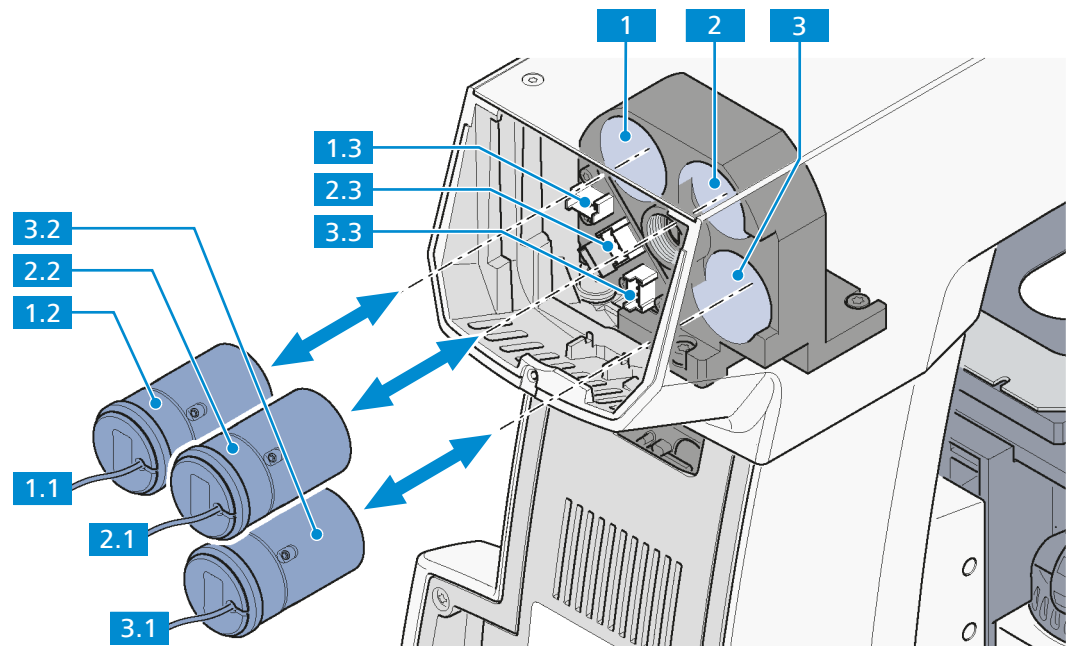

- Procedure**
1. Unplug the module plugs (1.1, 2.1, 3.1) from the connectors (1.3, 2.3, 3.3) of the stand socket.
  2. Remove the old LED modules (1.2, 2.2, 3.2) out of the LED tubes of the stand socket.
  3. Insert the new LED modules into the LED tubes (1, 2, 3) up to the stop.
  4. Ensure that the modules are inserted in the right holding positions.
  5. Plug the module plugs into the connectors of the stand socket.
  6. Ensure that the plugs are connected to the right connectors.

#### Info


The LED positions 1, 2 and 3 correspond to the UV, B and G labels on the right side of the Axiolab 5 Bio-TL/FL stand model.

#### Info

Due to the limited space in the LED module socket, the LED module in position 2 needs to be removed first when replacing the LED module in position 3.

### 10.9.4.3 Assembling the Module Holder

- Procedure**
1. Screw the module holder into the stand socket to fix the modules in place.

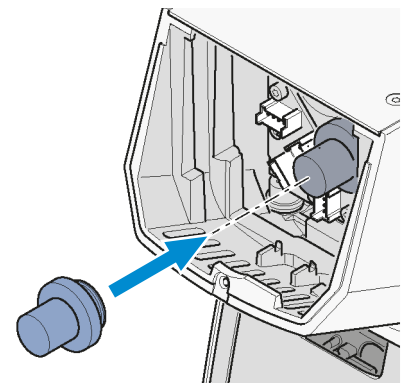

2. Insert the cover [▶ 136].



## 10.10 Axiocam 202 Mono/208 Color

**Purpose** The camera is used to snap photos or the microscopic image.

**Position** The Axiocam 202 mono or Axiocam 208 color is mounted on the camera port of the photo tube.


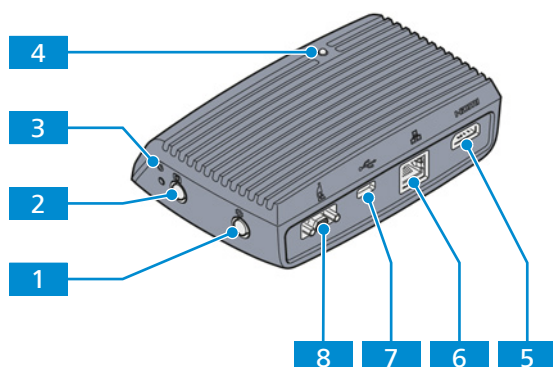


Fig. 53: Axiocam 202 mono/208 color

- |  |   |
|--|---|
| <b>1</b> OSD menu button   | <b>2</b> Image/video capture button   |
| <b>3</b> Camera factory reset button                                     | <b>4</b> Status LED   |
| <b>5</b> HDMI port for image data transfer to a monitor, TV or projector | <b>6</b> Gigabit Ethernet port (RJ45) for communication and image transfer                              |
| <b>7</b> Port for camera control and image transfer (USB 3.0)            | <b>8</b> Port for power supply and communication to the microscope stand (via Commercial Micro-D cable) |

### 10.10.1 Assembling the Axiocam 202 mono or Axiocam 208 color

- Parts and Tools**
- 🔧 C-mount camera adapter
  - 🔧 USB (Commercial Micro-D) cable (USB 2.0)
  - 🔧 HDMI cable

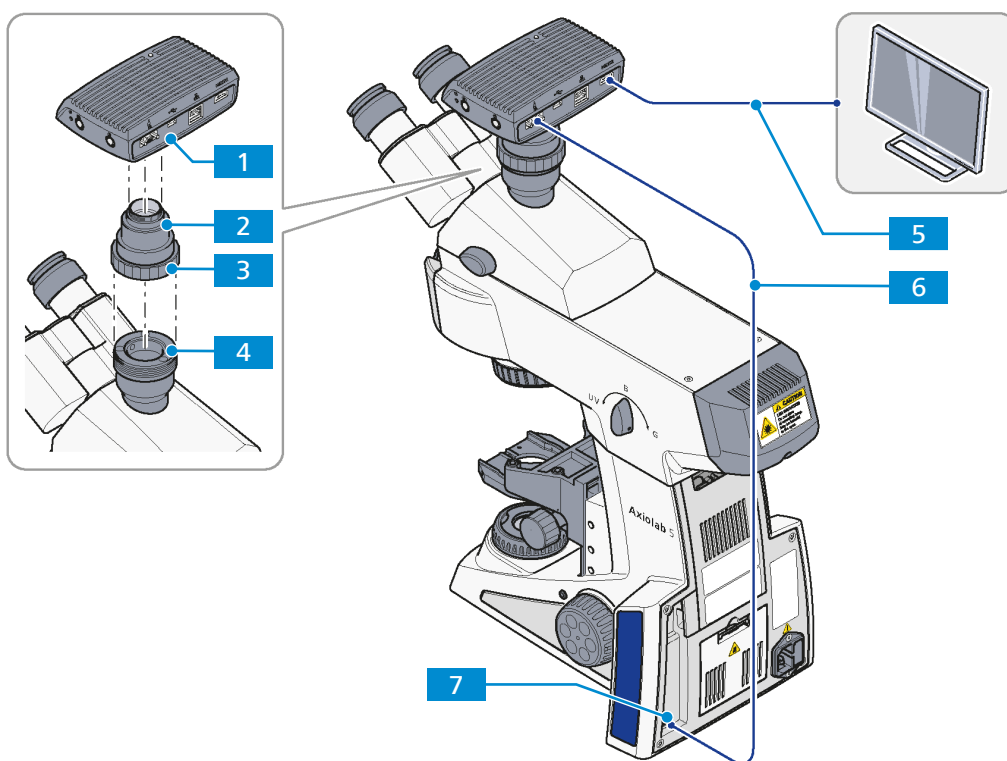

- Procedure**
1. Mount the C-mount camera adapter **2** on the Axiocam **1**.
  2. Attach the Axiocam with the adapter to the camera port **4** of the tube.
  3. Orient the camera to the stand and fix it in position by tightening the ring nut **3**.
  4. Connect the camera to the stand connector **7** via the USB (Commercial Micro-D) cable **6**.
  5. Connect the camera to an external monitor via an HDMI cable **5**.
  6. Alternatively, connect the camera to a WLAN router, USB Type-C drive or PC, see also *Operating Modes Using the Axiocam 202 mono/208 color* [▶ 143].

## 10.10.2 Operating Modes Using the Axiocam 202 mono/208 color

### 10.10.2.1 Axiocam as Standalone System

**Purpose** The camera is used to capture the microscopic image and store the data on the USB drive connected to the camera.

**Function** The camera acts as the control interface and is powered by the microscope via the USB (Commercial Micro-D power) cable.

A USB Type-C drive is included in the package and can be connected via the USB slot at the back of the camera for storing data. Then images are recorded and saved to the USB drive.

Functions of the microscope stand such as the Light Manager and encoding are automatically launched. The camera is equipped with image enhancement functions such as true color and noise reduction.

Functionality of the microscope:

- Light Manager
- Coded components
- Image enhancement (true color, noise reduction)
- Record and save images on the USB drive
- Record and save videos on the USB drive

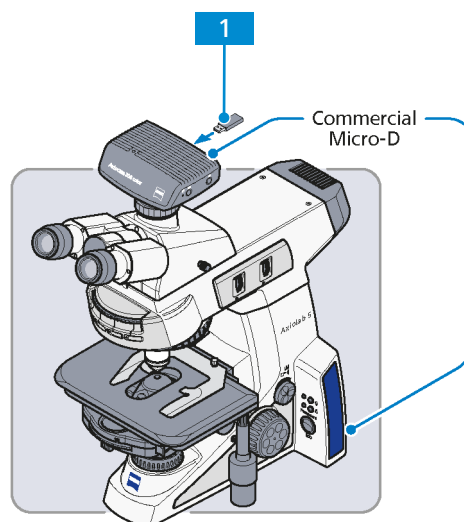


Fig. 54: Axiocam as standalone system

- 1** USB Type-C drive (included in package)

### 10.10.2.2 Axiocam connected to an HD monitor, TV or Projector

**Purpose** The camera is used to capture the microscopic image.

**Function** A monitor can be connected to the camera via an HDMI cable. The camera is powered by the microscope via the USB (Commercial Micro-D power) cable.

A monitor can be connected to the camera of the microscope via an HDMI cable. The camera is powered by the microscope via a USB (Commercial Micro-D power) cable. A USB hub can be connected via the USB port on the camera.

A wireless or wired mouse and keyboard can be connected to the camera via the USB hub, which together with the monitor, function as the control interface. Functions such as the Light Manager, encoding and image enhancement are automatically launched. Live images can be viewed on the monitor display and advanced features are available in the on-screen display (OSD).

When the microscope is operated together with the Colibri 3 light source, the one-key fluorescence function can be used. Images can be snapped and saved on the USB Type-C drive, which is connected via the USB hub.

Functionality of the microscope:

- Light Manager
- Coded components
- Image enhancement (true color, noise reduction)
- Observe live image on display
- Record and save image on the USB drive
- Record and save video on the USB drive
- One-key fluorescence (works only when the Axioscope is used with Colibri 3)
- Advanced features in OSD

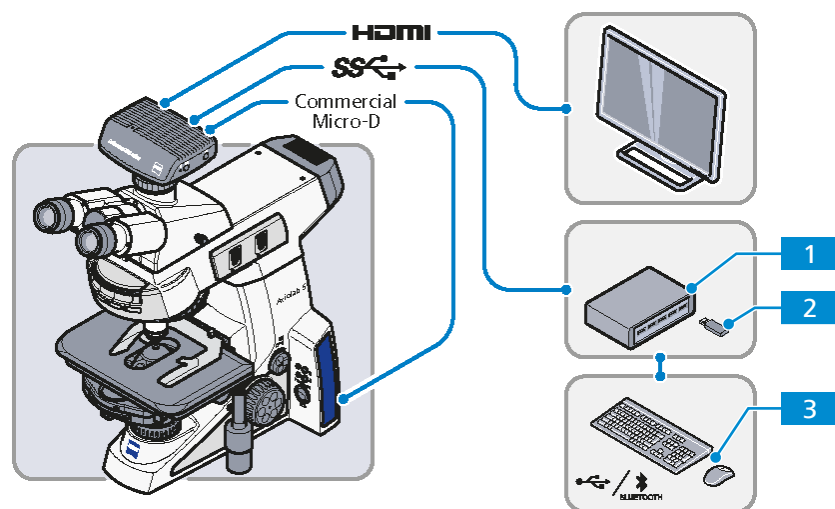


Fig. 55: Axiocam connected to an HD monitor, TV or projector

- 1** USB hub (input type C to output type A)    **2** USB Type-C drive provided in package  
**3** Mouse, keyboard

### 10.10.2.3 Axiocam connected with Labscope/Matscope via a Wi-Fi Dongle

**Purpose** The camera is used to capture the microscopic image.

**Function** The camera is powered by the microscope via the USB (Commercial Micro-D power) cable.

An optional monitor can be connected to the camera via an HDMI cable.

The recommended USB Wi-Fi dongle can be connected to the camera via the USB hub.

The control interface can be a PC or portable electronic device that uses Wi-Fi.

Functions such as the Light Manager, encoding, ECO mode and image enhancement are automatically launched.

When a monitor is connected, live images can be viewed on the monitor display. Live images can also be viewed on PC or portable devices and advanced features in Labscope/Matscope are available.

Functionality of the microscope:

- Light Manager
- Coded components
- ECO mode
- Image enhancement
- Observe live images
- Record and save images via software
- One-key fluorescence (this works only with Axiolab Bio-TL/FL)
- Advanced features in Labscope/Matscope

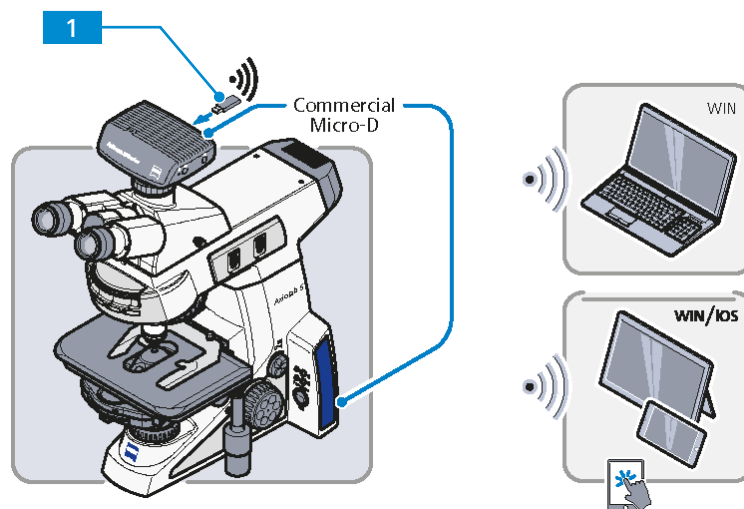


Fig. 56: Axiocam connected with Labscope/Matscope via a Wi-Fi dongle

- 1** USB Wi-Fi dongle (please see the ZEISS website for the recommended model)

#### 10.10.2.4 Axiocam connected with Labscope/Matscope via a WLAN Router

**Purpose** The camera is used to capture the microscopic image.

**Function** The camera is powered by the microscope via the USB (Commercial Micro-D power) cable.  
An optional monitor can be connected to the camera via an HDMI cable.

A router is connected to the camera via Ethernet.

The control interface can be a PC or portable electronic device controlled via Ethernet or Wi-Fi.

Functions such as the Light Manager, encoding, ECO mode and image enhancement are automatically launched.

When a monitor is connected, live images can be viewed on the monitor display. Live images can also be viewed on a PC or a portable device and advanced features in Labscope/Matscope are available.

Functionality of the microscope:

- Light Manager
- Coded components
- ECO mode
- Image enhancement
- Observe live images
- Record and save images via software
- One-key fluorescence (this works only with Axiolab Bio-TL/FL)
- Advanced features in Labscope/Matscope

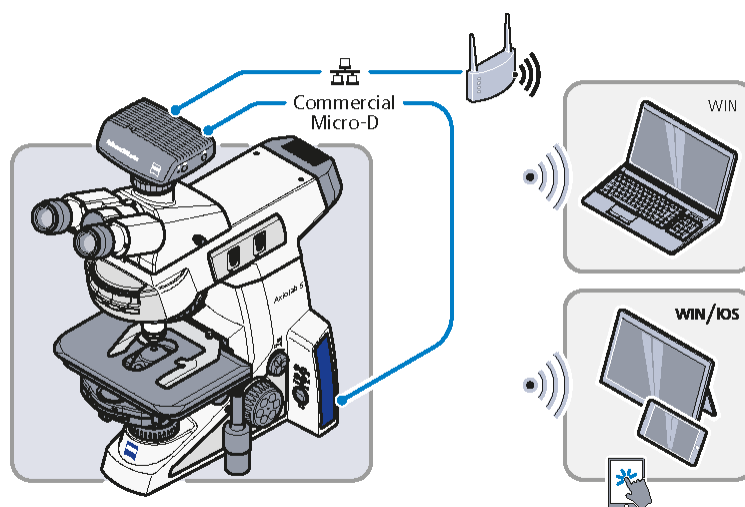


Fig. 57: Axiocam connected with Labscope/Matscope via a WLAN router

### 10.10.2.5 Axiocam connected with Labscope/Matscope via a USB

**Purpose** The camera is used to capture the microscopic image.

**Function** The camera is powered by the microscope via the USB (Commercial Micro-D power) cable.

An optional monitor can be connected to the camera via an HDMI cable.

A PC or Windows Surface can be connected to the camera via a USB cable.

Functions such as the Light Manager, encoding, ECO mode and image enhancement are automatically launched.

When a monitor is connected, live images can be viewed on the monitor display. Live images can also be viewed on a PC or Surface and advanced features in Labscope/Matscope are available.

Functionality of the microscope:

- Light Manager
- Coded components
- ECO mode
- Image enhancement
- Observe live images
- Record and save images via software
- One-key fluorescence (this works only with Axiolab Bio-TL/FL)
- Advanced features in Labscope/Matscope

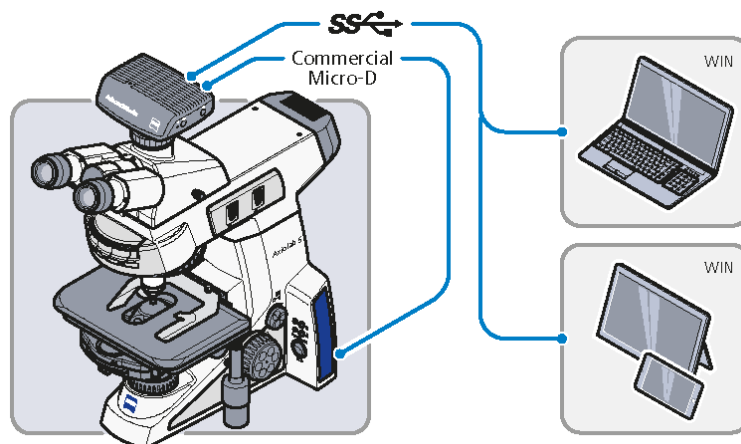


Fig. 58: Axiocam connected with Labscope/Matscope via a USB

### 10.10.2.6 Axiocam connected with ZEN software via a USB

**Purpose** The camera is used to capture the microscopic image.

**Function** The camera is powered by the microscope via the USB (Commercial Micro-D power) cable.

A workstation can be connected to the camera and the microscope stand via USB cables at the same time.

Functions such as the Light Manager, encoding and ECO mode are automatically launched.

Live images can also be viewed on the workstation and basic features in ZEN are available.

Functionality of the microscope:

- Light Manager
- Coded components
- ECO mode
- Image enhancement
- Observe live images
- Record and save images via software
- Basic features in ZEN


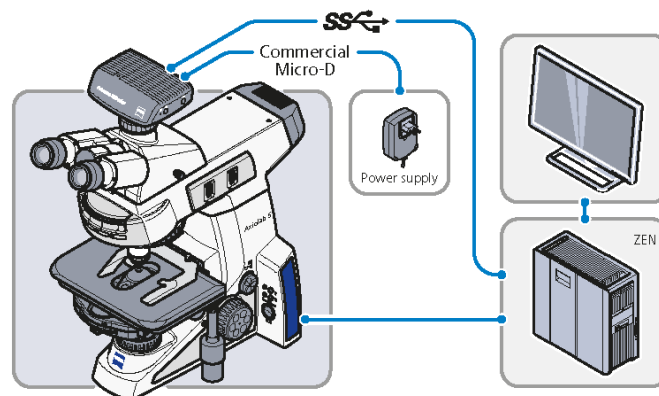


Fig. 59: Axiocam connected with ZEN software via a USB

### 10.11 Assembling the Filter 32x4 mm on the Luminous-field Diaphragm Operating Ring

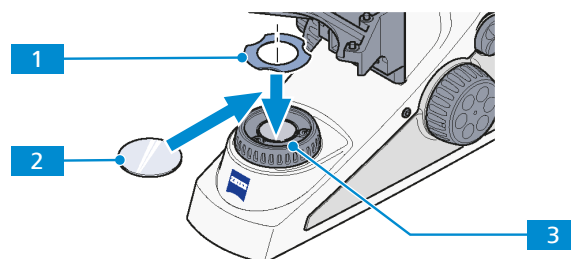

- Procedure**
1. If necessary, remove the filter clamp **1**.
  2. Place the filter **2** on the luminous-field diaphragm operating ring.
  3. To secure the filter, insert the filter clamp on the luminous-field diaphragm operating ring **3**.

Proceed in the reverse order for removal.

## 10.12 Dual Observation

**Purpose** The accessory unit allows two people to look at the samples through eyepieces at the same time. Both users will be able to have the simultaneous view of the sample.

**Position** The dual observation unit is mounted on the top of the stand.

**Function** The main observer (with the microscope stand) has a black stick to control the position of the arrow in the vision.

The following features and controls are available:

- Four-color LED (blue, green, red and white) arrow in the field of view to point at the region of interests
- Field of view 23 mm

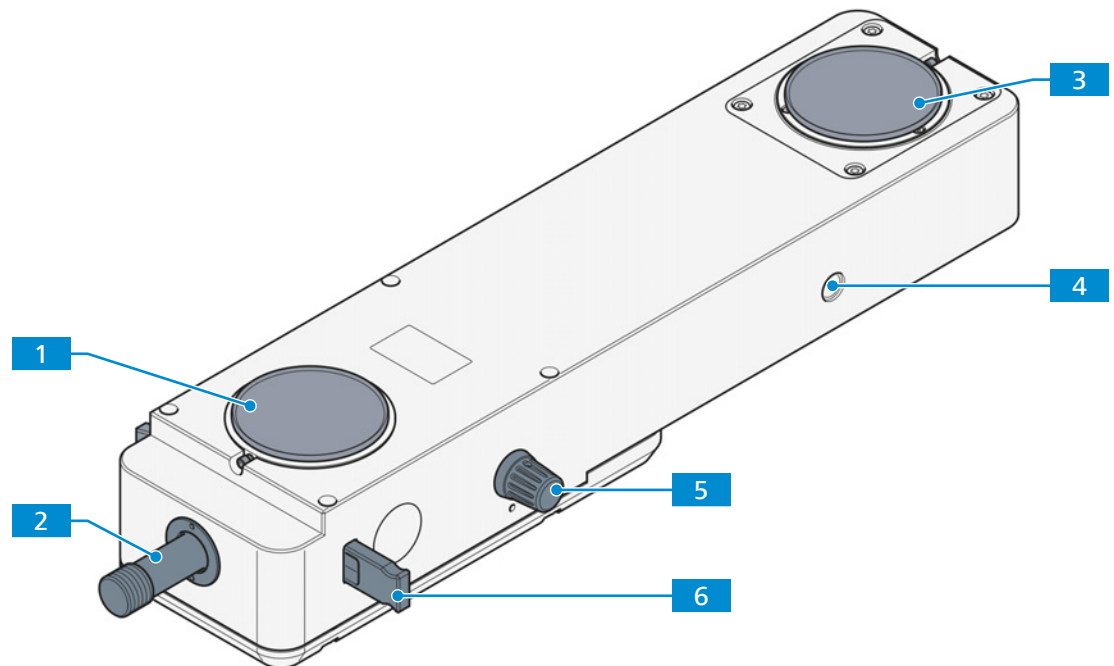


Fig. 60: Dual Observation

- |   |  |
|---|--|
| <b>1</b> Observation tube mount for main observer | <b>2</b> Joystick for LED arrow position control   |
| <b>3</b> Observation tube mount for co-observer   | <b>4</b> DC input socket (5V)                      |
| <b>5</b> ON/OFF switch                            | <b>6</b> Filter plate for LED arrow color changing |



### 10.12.1 Preparing the Binocular Tubes

1. Remove the tube lenses from all binocular tubes to be used [▶ 150].
2. Install the adaptation set to the binocular tube for the main observer [▶ 151].

#### 10.12.1.1 Removing the Tube Lens from the Binocular Tubes

##### NOTICE

##### Malfunction due to tube lens confusion


Binocular tubes and tube lenses are properly matched to one another.

- ▶ Mark binocular tubes and tube lenses in such a way that they can be assigned to each other again for later use.

**Parts and Tools**     Ring tool

**Prerequisite**     The binocular tube is removed from the microscope.

- Procedure**
1. Put the binocular tube **3** upside down on the table.

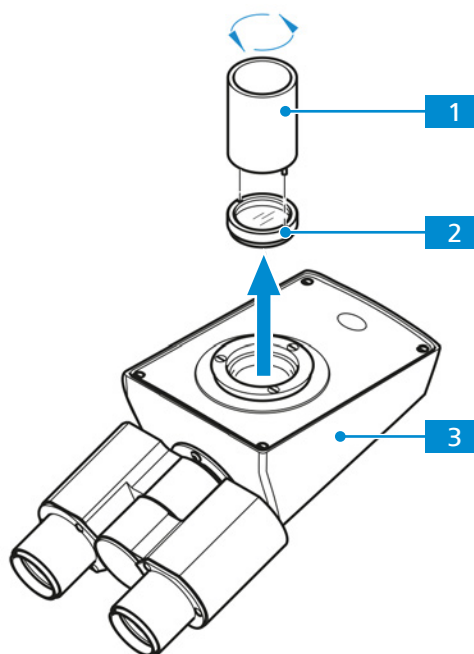

2. Remove the tube lens **2** using the ring tool **1**.
3. Store the tube lens in dust-free conditions.
4. Turn the binocular tube.

### 10.12.1.2 Inserting the Adaptation Set into the Binocular Tube for the Main Observer


The following glass filters for the adaptation set are available:

- neutral glass filter
- neutral density filter with 50% transmission
- neutral density filter with 25% transmission

**Parts and Tools**    Mounting tool

**Prerequisite**    The tube lens is *removed* [[▶ 150](#)].

- Procedure**
1. Put the adaptation set **4** with the retaining ring facing upwards on a clean, soft surface.
  2. Unscrew the retaining ring **2** using the mounting tool **1**.

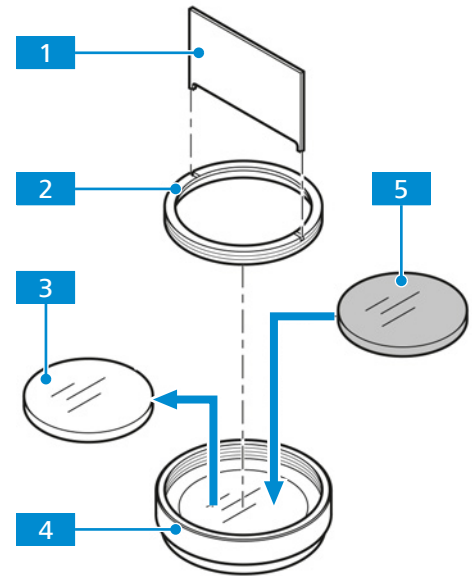

3. Carefully turn the adaption set upside down.  
→ The glass plate **3** slips out onto the soft surface.
4. Turn the adaption set.
5. Insert the required glass filter **5** into the adaptation set.
6. Tighten the retaining ring **2** using the mounting tool **1**.
7. Screw the adaptation set into the binocular tube.

## 10.12.2 Assembling the Dual Observation

**Info**

All tubes should be equipped with eyepieces of the same field diameter (preferably 23 mm) in order to obtain same-size fields of view for the main observer and the co-observer.

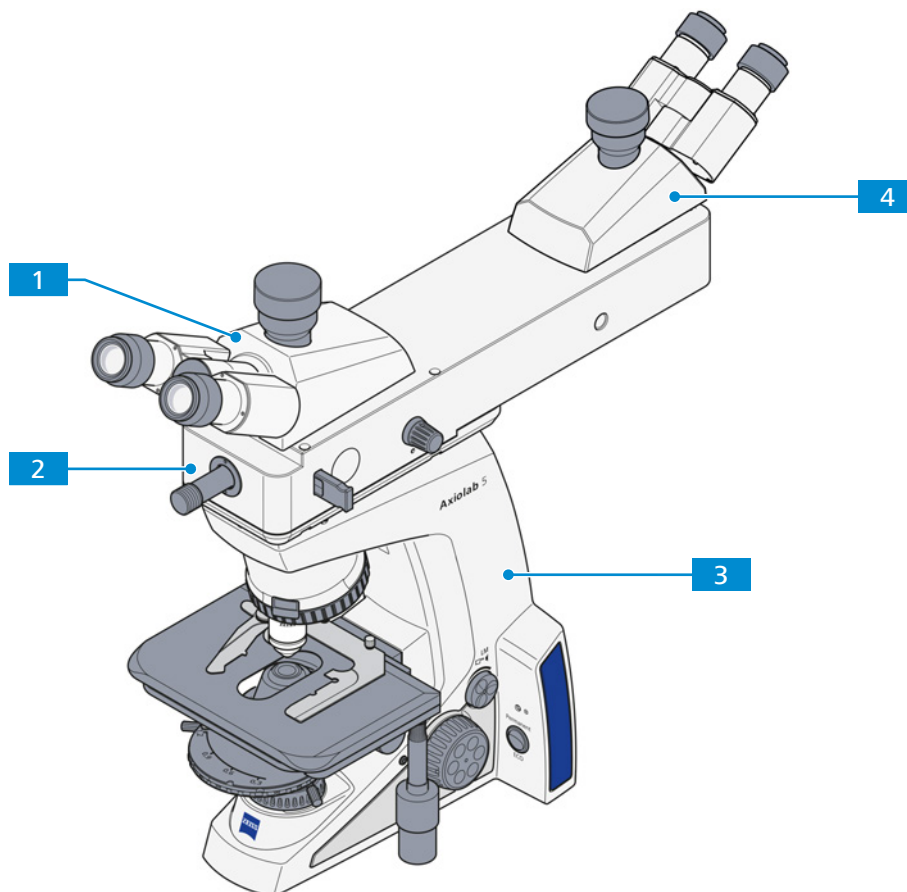


Fig. 61: Axiolab 5 with dual observation

- |   |   |
|---|---|
| <b>1</b> Binocular tube for the main observer | <b>2</b> Dual observation                   |
| <b>3</b> Microscope stand                     | <b>4</b> Binocular tube for the co-observer |


- Prerequisite**
- ✓ The tube lenses are *removed from both binocular tubes* [▶ 150].
  - ✓ The adaptation set is *installed to the binocular tube for main observer* [▶ 151].
  - ✓ The microscope is placed on a suitable table.

- Procedure**
1. Install the dual observation **2** on the microscope stand **3** with the joystick facing to the main observer.
  2. Install the *binocular tube* [▶ 53] for the main observer **1**.
  3. Install the *eyepieces* [▶ 55] for the main observer.
  4. Install the binocular tube for the co-observer **4**.
  5. Install the eyepieces for the co-observer.
  6. Connect the plug-in power cable to the microscope stand.

### 10.12.3 Operating the Dual Observation


- Prerequisite**
- ✓ The dual observation is assembled.
  - ✓ The microscope is switched on.

- Procedure**
1. Switch on the dual observation by turning the rotary button clockwise.

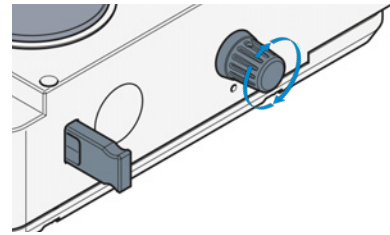


→ The blue LED below the knob lights up.


2. Adjust the light intensity by turning the rotary button.
3. Position the LED arrow in the field of view by means of the joystick.

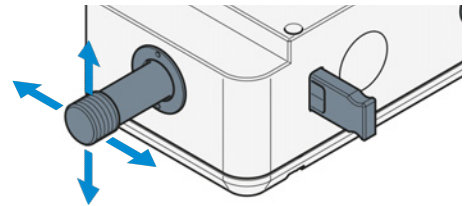

4. Set the color of the LED arrow by pulling or pushing the filter plate.

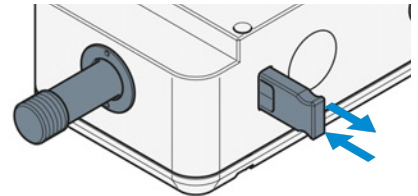

5. Re-focus the eyepieces of the co-observer by approx. 1/3 diopter to compensate the glass path of the removed tube lens.
6. Switch off the dual observation by turning the rotary button counterclockwise.

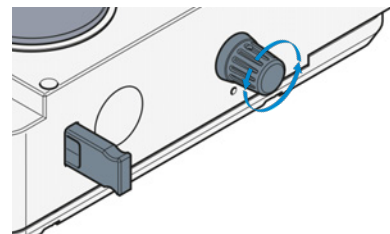

→ The blue LED below the knob goes out.

## Revision History

Revision	Date of Issue	Introduced Modifications
11	05/2024	<ul style="list-style-type: none"> <li>Editorial rework</li> <li>Correction of chapter(s):               <ul style="list-style-type: none"> <li>– <i>Performance Data and Specifications</i> [▶ 104]</li> <li>– <i>Accessories and System Expansions</i> [▶ 108]</li> <li>– <i>Binocular Photo Tube 30°/23 (100:0/30:70/0:100), Reversed Image</i> [▶ 114]</li> <li>– <i>Dual Observation</i> [▶ 149]</li> </ul> </li> </ul>
10	02/2024	<ul style="list-style-type: none"> <li>Editorial rework</li> <li>Changes on Light Manager</li> <li>Components and accessories added</li> </ul>
9	04/2023	<ul style="list-style-type: none"> <li>Editorial rework</li> <li>Implementation of UKCA marking</li> </ul>
8	05/2022	<ul style="list-style-type: none"> <li>Implementation of revision history</li> <li>Adaptation to Regulation (EU) 2017/746 (IVDR)</li> </ul>

Tab. 2: Revision History

# Glossary

---

**BF (Brightfield)**

Illumination and imaging system where direct light passes through the objective aperture and provides a bright background against which the image is viewed.

---

**C-DIC**

Differential Interference Contrast in circularly polarized light, a contrast method which employs the differential interference contrast technique with circularly polarized light, thus fully imaging sample structures which otherwise are only visible in a certain orientation

---

**DF (Darkfield)**

Illumination and imaging system that prevents direct light from entering the objective aperture.

---

**DIC (Differential Interference Contrast)**

An imaging light microscopy method that converts differences in the optical path length in the object into differences in the brightness of the image

---

**FL (Fluorescence)**

Phenomenon of a selective absorption of radiation with relatively short wavelength (i.e., relatively high energy) by matter with the result of the emission of radiation with longer wavelengths (i.e., lower energy), which persists only very briefly after the excitation has ceased.

---

**RL (Reflected Light)**

Designation for microscopy techniques to image light that was reflected by the object

---

**TIC (Total Interference Contrast)**

Total Interference Contrast in circularly polarized light is a technique for imaging and layer thickness measurement in materials microscopy. Contrary to traditional polarization interferometers, work in this technique is carried out in circularly polarized light.

---

**TL (Transmitted light)**

Light used for illuminating a object, where the light is transmitted through the object.

---

**ZEISS Sales & Service Partner**

The Sales & Service Partner is generally in the field for customer support in a regional area and / or a clearly defined customer group.

---

**ZEISS service representative**

Specially trained service expert, either ZEISS staff or authorized service partner of ZEISS.

---

**ZEN software**

Modular software, that is controlling all ZEISS light microscope systems and has a wide application field: acquiring images, processing images and analyzing images.

# Index

## A

Accessories	108
Adaptation set	151
Air Conditioning and Quality	104
Ametropia	64
Analyzer slider	123
Aperture diaphragm	40, 41, 129
Auxiliary microscope	55
Axiocam	142, 143
Axiolab 5 Bio-TL	21
Axiolab 5 Bio-TL/FL	22
Axiolab 5 Mat-TL/RL	24
Axiolab 5 Pol-TL	22
Axiolab 5 Pol-TL/conoscopy	23
Axiolab 5 Pol-TL/RL	23

## B

Binocular photo tube	34, 35, 36, 111, 112
Binocular tube	111
installing	53, 54
remove tube lens	150
Bireflectance	51
Birefringence	46, 75, 78
Brightfield	45, 50, 68, 83

## C

Camera	142
installing	143
Filter carrier	
installing	128
Circular polarization contrast	48, 77, 82
Circular polarizer	126
Clamping device	43
installing	120
removing	119
Cleaning	
Water-soluble contaminations	95
Color filter carrier	125, 127
Components of the stand	25, 26, 27, 28, 29, 31
Condenser	40, 129
darkfield	58
installing	57, 58, 130
mounting	59
Condenser carrier	39
Darkfield condenser	
installing	58
Conoscopic observation	50
Connecting to the mains	62
Contamination	103

## Cover

installing	132, 136
removing	132, 136
Crystals, determining the optical character	49, 81

## D

Darkfield	45, 51, 72, 86
Decontamination	103
Annular phase diaphragm	
centering	131
Darkfield diaphragm	
centering	131
DIC slider	129
Display elements	32
Disposal	103
Dual observation	
installing	152
operating	153
Dual observation	149

## E

ECO mode	68
Reversible eyecups	
installing	110
Eyepiece	38, 55
Eyepiece reticle	38
Eyepieces	38
Eyepiece reticle	
adjusting for ametropia	64

## F

Filter	125, 127
installing	148
Fluorescence	51, 89
Friction of coaxial knurled knobs of the stage	97
Fuse	
exchange	96

## G

General Safety Information	12
----------------------------	----

## H

Halogen lamp, replacing	133, 137
-------------------------	----------

Hazards	14
Biological	16
Electromagnetic interference	17
Ergonomic	15
Infection	16
Operating Environment	15
Optical radiation	17
Prevention	14
Skin irritation	16
Thermal	17
Transport	14

## I

Improper use	11
Intermediate plate	54
Interpupillary distance	64

## K

Key functions	32
---------------	----

## L

Lambda 4/plate	126
Lambda plate	125
Light Manager	
function	44
saving ratios	67
Switching Off	66
Switching On	66
Fluorescence light source	
exchanging LED	140
LED light source	
exchanging	134, 137
Low-power system	126
centering	127
installing	127

## M

Mains connection	104
Mains, connecting to	62
Maintenance	93
interval	94
schedule	94
Mechanical stage	42, 96, 116, 117
adjusting the friction	97
installing	115
removing	116
setting the drive length	96
Modulator disk	
installing	60

## N

Nosepiece	39
-----------	----

## O

Objective	39
centering	122
Operation	
Prerequisites	63
Optical character of crystals	49, 81
Optional system expansions	108
Installation	108

## P

Path differences	48, 76, 80
Performance data	104
Permanent mode	68
Phase contrast	45, 74
Photo tube	34, 35, 36, 111, 112
Pinhole diaphragm	55
Polarization	46, 50, 74, 87
Polarization direction	46, 76, 79



Polarization contrast	
circular	77, 82
Polarizer	124, 125, 126
installing	128
Prerequisites	
Operation	63

## R

Reflected light	
brightfield	50, 83
darkfield	51, 86
fluorescence	51, 89
polarization	51, 87
Reflector turret	43
Reflector module	
installing	61
removing	62
Reflector turret	61
Requirements	
for Operators	12
Reticle	38
Rotary stage	43, 118
centering	121
installing	120
removing	119

## S

Safe Operating Condition	12
Safety	11, 93
devices	19
interlocks	19
Sample guide	118
installing	120
removing	119
Sample holder	
installing	115
removing	116
Factory settings	101
Shutdown	102
Software	10
Spare parts	13
Specimen holder	42, 116, 117
Storage	102
Storage compartment	44
Switching off	92
Switching on	63

## T

Tool	44
Training	12
Transmitted light polarization	50
Transmitted light	
brightfield	45, 68
darkfield	45, 72
phase contrast	45, 74
polarization	46, 74
Transport	102

Troubleshooting	98
Tube	34, 35, 36, 111, 112
assembling components	55

## U

Unpacking	52
-----------	----

## V

Viewing height	64
----------------	----

## W

Warning	
labels	17, 18
lights	17
Weight and Sizes	104

## Z

ZEISS	
Portal	10
Service agreements	93



**Carl Zeiss Microscopy GmbH**

Carl-Zeiss-Promenade 10  
07745 Jena  
Germany

phone: +49 1803 33 63 34  
fax: +49 3641 64 3439

[info.microscopy.de@zeiss.com](mailto:info.microscopy.de@zeiss.com)  
[www.zeiss.com/microscopy](http://www.zeiss.com/microscopy)